

QuantStudio™ 6 Pro Real-Time PCR System and QuantStudio™ 7 Pro Real-Time PCR System

USER GUIDE

Publication Number MAN0018045

Revision H.0



For Research Use Only. Not for use in diagnostic procedures.

ThermoFisher
SCIENTIFIC



Life Technologies Holdings Pte Ltd | Block 33 | Marsiling Industrial Estate Road 3 | #07-06, Singapore 739256
For descriptions of symbols on product labels or product documents, go to thermofisher.com/symbols-definition.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. MAN0018045

Revision	Date	Description
H.0	28 July 2021	v1.5.0 <ul style="list-style-type: none">Updated the RNase P instrument verification after calibration to a recommendation only.Added instructions to select the language, to enable or disable the reminder to insert a plate, and to import a TaqMan™ Assay file.Updated the instructions to manage the naming convention for the data file.Removed the connection option of the instrument when connecting to Thermo Fisher™ Connect (connection options are mobile device and computer).Updated the Security, Auditing, and E-signature (SAE) Administrator Console to v2.1 or later.
G.0	17 February 2021	v1.4.0 Updated centrifuges for the TaqMan™ Array Card.
F.0	6 November 2020	v1.4.0 <ul style="list-style-type: none">Specified that the instrument and the desktop software must be connected to the same SAE server for an account password to be updated on both applications.Provided the option of a static IP address or a dynamic IP address to connect to the SAE server.Corrected standards on 384-well RNase P plate (Figure 14 on page 105).Added a TaqMan™ Array Card block, including system template file names, calibration workflow, instructions to edit the plate file, and instructions to load and unload from the instrument (QuantStudio™ 7 Pro System only).Added instructions to review the calibration status in the QuantStudio™ Design and Analysis Software v2.Added a peripheral self-test and updated the previous self-verification test to a hardware test.Added instructions to upload a certificate for a TaqMan™ Array Plate with RFID.Added the option of a voice recording for Smart Help.Added remote support.Updated the RFID labels (Figure 27 on page 216 and Figure 28 on page 217).
E.0	27 April 2020	v1.3.0 <ul style="list-style-type: none">Added protein thermal shift experiments to the instrument functions.Added a 96-well 0.1-mL block.Updated instructions to export data files when setting up a run and added instructions to import or delete export settings.Updated instructions to save an edited run method.Updated instructions to add or edit reagents and then to apply the reagents to the plate.Updated instructions to view the amplification plot after a run.Added instructions to view QC Check status after the run.Added information about a locked template file.Added functions for using the instrument with the SAE Administrator Console.

Revision	Date	Description
D.0	7 January 2020	<p>v1.2.1</p> <ul style="list-style-type: none"> Added information about position to set up facial authentication and to sign in with facial authentication. Added instructions to test the connection to the Connect cloud-based platform and to view the privacy policy for the Connect platform. Added instructions to repair the software after a software update. Changed port 7000 to 7443 for the instrument-to-computer connection. Corrected C_t to C_q in description of instrument verification.
C.0	3 December 2019	<p>v1.2.0</p> <ul style="list-style-type: none"> Added a 384-well plate block. Updated instructions to set up facial authentication. Added a new voice command for calibration. Added custom melt curve run calibration. Added Security, Auditing, and E-signature (SAE) v2.0 module. Added instructions to generate and reset an access key for external software to connect to the instrument. Added instructions to import an Assay Information File. Corrected that a Connect profile is required for the instrument schedule and Smart Help. Added new radio compliance standards.
B.0	22 August 2019	<p>v1.1.0</p> <ul style="list-style-type: none"> Corrected the text on the buttons for the following items: <ul style="list-style-type: none"> Add facial authentication to an existing instrument profile. View the license agreement. Reset the instrument (restore factory default). Added the following items: <ul style="list-style-type: none"> TET™ dye as a system dye. Information about sleep mode. Information about the automatic sign-out feature. The option to configure the instrument so that sign-in is required (guest access is not permitted). Instructions to set up email notification. Information about the speaker volume for each user. Instructions to back up the instrument and to restore a backup of the instrument. Information about multiple instruments and email notification in the comparison of a local instrument profile and a Connect profile. Instructions to unlink a Connect profile. Instructions to update a PIN for a Connect profile. Lighting guidelines for facial authentication. Radio compliance standards and RF transceiver specifications. Corrected the following items: <ul style="list-style-type: none"> Indicators for voice commands. Voice commands are only available in select countries. Instructions to remove VeriFlex™ Zones and the Auto Delta setting. Excitation and emission wavelengths of the filters in the calibration instructions. Add a four-digit PIN when signing into the instrument for the first time using a Connect profile. Add a four-digit PIN when linking a local instrument profile to a Connect profile.

Revision	Date	Description
B.0 (continued)	22 August 2019	<ul style="list-style-type: none"> – Select a region for the Connect cloud-based platform. – A run cannot be started if a block has never been calibrated on the instrument. – The options after a calibration run is complete. – The example files for the desktop software are embedded in the software and do not need to be downloaded. • Updated the following items: <ul style="list-style-type: none"> – C_q Export checkbox is selected as the default. – Instructions to sign in with facial authentication.
A.0	18 April 2019	New document (v1.0.1).

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

NOTICE TO PURCHASER: DISCLAIMER OF LICENSE: Purchase of this software product alone does not imply any license under any process, instrument or other apparatus, system, composition, reagent or kit rights under patent claims owned or otherwise controlled by Thermo Fisher Scientific, either expressly, or by estoppel.

Trademarks: All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. TaqMan is a registered trademark of Roche Molecular Systems, Inc., used under permission and license. Cy is a registered trademark of Global Life Sciences Solutions Germany GmbH. Amazon, Alexa, and all related logos are trademarks of Amazon.com, Inc. or its affiliates. Windows is a trademark of Microsoft Corporation. Chrome is a trademark of Google LLC. Pentium is a trademark of Intel Corporation or its subsidiaries in the U.S. and/or other countries. Macintosh is a trademark of Apple Inc., registered in the U.S. and other countries.

©2021 Thermo Fisher Scientific Inc. All rights reserved.

Contents

■ CHAPTER 1 Product information	14
Network and password security requirements	14
Network configuration and security	14
Password security	14
Instrument hardware description	15
Overview of the instrument	15
Features of each instrument	16
Instrument filters and supported dyes	16
Overview of data collection	18
VeriFlex™ Zones	18
AutoDelta settings	19
Parts of the instrument	20
Instrument status indicator	22
Hands-free features	22
Facial authentication	22
Voice commands	22
Indicators for the hands-free features	23
Software description	24
Parts of the home screen	24
Thermo Fisher™ Connect definitions	26
Instrument, desktop software, and Connect platform features	26
Overview of file locations and files	27
Components of the files	30
Third-party software	31
Network connection options	31
Types of runs	31
Standard curve run	31
Relative standard curve run	32
Comparative C_t ($\Delta\Delta C_t$) run	32
Genotyping run	33
Presence/absence run	33
Melt curve run	33

System template file names	34
System template file names for the 96-well 0.2-mL block	34
System template file names for the 96-well 0.1-mL block	35
System template file names for the 384-well block	37
System template file names for the TaqMan™ Array Card block	38
Workflow	40
CHAPTER 2 Get started	41
Precautions for use	41
Installation and instrument verification	42
Overview of guest access	42
Use the instrument as a guest	42
Power on the instrument	43
Sign in	43
Sign in with a PIN	43
Sign in with facial authentication	44
Enable or disable the microphone	44
Use voice commands	45
Sign out	46
CHAPTER 3 Create and run plates on the instrument	47
Options to run a plate	47
Create and run a plate from a system template or a saved plate file	47
Repeat the last run	49
Overview of customized data export	50
(Optional) Scan a barcode using a barcode scanner	51
Load a plate into the instrument	51
Load a TaqMan™ Array Card into the instrument	52
Unload a plate or a TaqMan™ Array Card from the instrument	52
CHAPTER 4 Edit a plate before starting a run	54
Overview of a template file with restricted editing	54
Options to edit the properties of a plate file	55
Import a sample setup	55
Import an auto export setting	55
Delete an auto export setting	55
Options to edit the method of a plate file	56
Method elements	57
Edit the method	57
Add or remove steps	58
Add or remove stages	58
Add or remove data collection locations	58

Set up or remove VeriFlex™ Zones	59
Set up or remove the Auto Delta feature	61
Configure ramp rates	62
Add or remove a pause	62
Options to edit the well details of a plate file	62
Options to view a plate layout	63
Assign or edit a well for a plate	65
Assign or edit a well for a TaqMan™ Array Card	66
Add or edit the reagent details for a plate	66
Apply the reagents to a plate	67
Delete the reagent details	67
Add or edit the reagent details for a TaqMan™ Array Card	68
Apply the reagents to a TaqMan™ Array Card	68
Edit a passive reference dye	68
Import an Assay Information File (AIF)	69
Import a TaqMan™ Assay file	69
■ CHAPTER 5 Options for a run	70
Screens available during a run	70
Edit a plate file during a run	71
Pause an instrument run	72
Stop an instrument run	72
View the method during a run	73
View real-time data and plots during a run	73
Adjust the display of real-time plots during a run	74
View the time remaining during a run	74
Lock the touchscreen during a run	74
Unlock the touchscreen during a run	75
Instrument Schedule	75
Overview of the instrument schedule	75
Schedule a run	75
Cancel a scheduled run	75
■ CHAPTER 6 View and manage files	77
Manage plate files	77
Copy plate files	77
Delete plate files	78
Manage data files	78
View QC Check status after a run	78
View an amplification plot after a run	80
Transfer a data file immediately after a run	81
Transfer data files at a later time	81

View the list of data files on the instrument	82
Delete data files from the instrument	82
■ CHAPTER 7 Calibrate and verify instrument performance	84
Instrument and block calibration	84
Calibration and verification schedule	85
Calibration descriptions	86
View the calibration status and set reminders	87
View the calibration status on the instrument	87
Set calibration frequency and reminders on the instrument	88
View calibration status and set reminders on the Connect platform	88
Review calibration status	89
Export a calibration report	89
Perform ROI/uniformity, background, and dye calibrations for plate blocks	89
Calibration workflow	90
Prepare a calibration plate	90
Perform calibrations	91
View the calibration images	93
Transfer calibration results	94
Transfer calibration data files	94
Troubleshoot calibration failure	95
Identify contamination	95
Create a background plate (<i>optional</i>)	96
Perform ROI/uniformity, background, dye, and normalization calibrations for the TaqMan® Array Card block	97
Calibration workflow	97
Prepare a calibration TaqMan® Array Card	98
Perform calibrations	100
View the calibration images	101
Troubleshoot normalization calibration failure	102
Perform verification using RNase P plates	103
Instrument verification description	104
RNase P instrument verification plate	104
Performance specifications pass criteria	106
Prepare an RNase P plate	106
Perform RNase P verification	107
Troubleshoot verification failure	108
Perform verification using RNase P with a TaqMan® Array Card	109
RNase P instrument verification TaqMan® Array Card	109
Prepare an RNase P TaqMan® Array Card	110
Perform RNase P verification	112

Calibrate custom dyes	113
Overview of custom dyes	113
Use a dilution series to determine an optimal custom dye concentration	114
Calibrate the custom dye	117
Calibrate for a custom melt curve run	120
CHAPTER 8 Manage profiles	122
Require sign-in	122
Recommended order to set up profiles	123
Overview of local instrument profiles and Connect profiles	123
Connect instrument profile roles and functions	124
Test the connection to the Connect platform	124
Link the instrument to your Connect account	126
Create a new local instrument profile	127
Link a local profile to a Connect profile	127
Unlink a Connect account	129
If you link when you <i>are</i> signed in to the instrument	129
If you link when you <i>are not</i> signed in to the instrument	130
Manage individual instrument profiles	131
Create an administrator instrument profile during initial start-up	131
Set up facial authentication at the initial creation of an instrument profile	131
Add facial authentication to an existing instrument profile	132
Change the PIN for an instrument profile	133
Change the PIN for a Thermo Fisher™ Connect profile	133
Manage all instrument profiles as an administrator	133
View all instrument profiles	133
Reset the PIN for a profile	134
Add a profile	134
Enable or disable administrator privileges	135
Delete a profile	135
CHAPTER 9 Use the instrument with the Security, Auditing, and E-signature (SAE) Administrator Console	136
Overview of the SAE Administrator Console components	137
Overview of the instrument features when the SAE module is enabled	137
SAE Administrator Console functions	138
Functions that are controlled	138
Functions that can be audited	139
Functions that can be signed	139
SAE Administrator Console roles	139
Dialog boxes in the SAE module	140
Profiles when the SAE module is enabled	141

SAE-enabled system components	141
Compatibility between SAE-enabled and SAE-disabled components	142
Enable SAE functions	143
Workflow: Enable SAE functions	143
Configure SAE Administrator Console application profiles	143
Enable SAE on the instrument and specify the SAE server (administrator only)	143
Connect to the SAE server	144
Enable SAE functions in QuantStudio™ Design and Analysis Software v2	145
Disable SAE on the instrument (administrator only)	145
Sign in as a local administrator with SAE enabled	146
Sign in with SAE enabled	147
Sign out with SAE enabled	148
Change your SAE account password on the instrument	149
Sign a plate setup	150
Enter an audit reason	151
View signing records for a plate file	152
View audit records	153
Use the instrument when the SAE server is offline	154
E-signature requirements to start a run	154
Signing in after automatic screen locking	155
SAE error messages and actions	155
■ CHAPTER 10 Maintain the instrument	156
Install a block	156
Change the block	157
Remove the block	158
Decontaminate the sample block	159
Materials required	159
Clean the sample block	159
Detailed procedures for cleaning the sample block	160
Replace the instrument fuses	161
Materials required	161
Replace the fuses	161
Power on or off, store, and move	163
Enable and edit sleep mode	163
Power on the instrument	164
Power off the instrument	164
Prepare the instrument to store, move, or ship	165
Move the instrument	165

■ CHAPTER 11 Configure the instrument settings	166
View the instrument details	166
View the license agreement	167
View the privacy policy for the Connect cloud-based platform	167
Configure the instrument settings	168
Enable or disable demonstration mode	168
Edit the instrument name (administrator only)	168
Set the date and the time	168
Configure the heated cover settings	169
Adjust the speaker volume	169
Configure the network	169
Select the region for the Connect platform	170
Select a disk storage management mode	171
Enable and edit automatic sign-out	171
Set an access key for external software to connect to the instrument	171
Reset the access key for external software to connect to the instrument	172
Configure maintenance and service settings	172
Request technical support with Smart Help	172
Request instrument service with Smart Help	173
Perform and view self-verification tests	173
View the instrument log	178
Export the instrument log	179
View the instrument statistics	179
Enable or disable remote monitoring	180
Update the software	180
Repair the software	181
Back up the instrument	181
Restore a backup of the instrument	182
Upload a certificate for a TaqMan™ Array Plate with RFID	182
Overview of an instrument reset (restore factory default)	183
Reset the instrument (restore factory default)	183
Request remote support	183
Configure the instrument for the hands-free features	185
Manage the run history	185
Manage the file name convention of the data file	185
Enable or disable the reminder to insert a plate	186
Select the language	186
View the notifications	187
Enable and disable home screen notifications	187
Set up email notifications	187
Clear the notifications	188

■ APPENDIX A Instrument specifications and layout	189
Instrument dimensions	189
Instrument and computer clearances	190
Configured system dimensions	191
Light guidelines	192
Electrical requirements	192
Environmental requirements	193
■ APPENDIX B Connect the instrument to a network	194
Connect the computer to the instrument directly or to a LAN	194
Connect the instrument to a wired network	195
Connect the instrument to a wireless network	196
Instrument and computer connections	197
Download and install the desktop software	198
Computer requirements for the desktop software	198
Download the desktop software	198
Install the desktop software	199
Networking	199
Supported options for instrument and computer connections	199
Control and monitor networked instruments	201
Networking guidelines and best practices	201
Firewall ports that must be open	201
Ethernet port overview	202
Third-party software	202
■ APPENDIX C Detailed procedures for preparation of a TaqMan™ Array Card	203
Guidelines for preparation of a card	203
TaqMan™ Array Card diagram	204
Load the PCR reaction mix	205
Centrifuge the card	205
Seal the card	207
■ APPENDIX D Parts and materials	210
Kits, consumables, accessories, and reagents	210
96-well 0.2 mL consumables	210
96-well 0.1 mL consumables	211
384-well consumables	211
TaqMan™ Array Card consumables	212
Accessories	213

■ APPENDIX E Safety	214
Symbols on this instrument	214
Standard safety symbols	215
Additional safety symbols	216
Location of safety labels	216
Control and connection symbols	220
Conformity symbols	220
Safety information for instruments not manufactured by Thermo Fisher Scientific	221
Instrument safety	221
General	221
Physical injury	222
Electrical safety	223
Laser safety	224
Cleaning and decontamination	225
Instrument component and accessory disposal	225
Safety and electromagnetic compatibility (EMC) standards	225
Safety standards	226
EMC standards	226
Environmental design standards	227
Radio compliance standards	227
RF transceiver specifications	229
Chemical safety	230
Biological hazard safety	231
■ APPENDIX F Documentation and support	232
Related documentation	232
Customer and technical support	233
Limited product warranty	233
Index	234

■ Network and password security requirements	14
■ Instrument hardware description	15
■ Hands-free features	22
■ Software description	24
■ Types of runs	31
■ System template file names	34
■ Workflow	40

Network and password security requirements

Network configuration and security

The network configuration and security settings of your laboratory or facility (such as firewalls, anti-virus software, network passwords) are the sole responsibility of your facility administrator, IT, and security personnel. This product does not provide any network or security configuration files, utilities, or instructions.

If external or network drives are connected to the software, it is the responsibility of your IT personnel to ensure that such drives are configured and secured correctly to prevent data corruption or loss. It is the responsibility of your facility administrator, IT, and security personnel to prevent the use of any unsecured ports (such as USB, Ethernet) and ensure that the system security is maintained.

Password security

Thermo Fisher Scientific strongly recommends that you maintain unique passwords for all accounts in use on this product. All passwords should be reset upon first sign in to the product. Change passwords according to your organization's password policy.

It is the sole responsibility of your IT personnel to develop and enforce secure use of passwords.

Instrument hardware description

Overview of the instrument

The Applied Biosystems™ QuantStudio™ 6 Pro and 7 Pro Real-Time PCR Systems use fluorescent-based polymerase chain reaction (PCR) reagents to perform the following functions:

- Quantitative detection of target nucleic acid sequences (targets)
- Qualitative detection of targets (endpoint analysis)
- Qualitative analysis of the PCR product (post-PCR target nucleic acid sequences melt curve analysis)
- Protein thermal shift experiments



Each instrument is available with an interchangeable block. The following blocks are available:

- 96-well 0.2-mL block
- 96-well 0.1-mL block
- 384-well block
- TaqMan™ Array Card block (QuantStudio™ 7 Pro System only)

The instrument can be used directly from the touchscreen to create and start runs. To design runs or to analyze data, the instrument can be integrated with the desktop software or the Connect platform.

Each instrument has the following features:

- Sign-in with facial authentication
- An interchangeable block
- Microphone for voice commands
- Speakers for feedback from voice commands and for instructional videos

- Smart Help to request technical support or instrument service directly from the instrument
- Barcode scanner for tracking plates or TaqMan™ Array Cards

Features of each instrument

Feature	QuantStudio™ 6 Pro System	QuantStudio™ 7 Pro System
Filter set ^[1]	Coupled five-color filter set	Decoupled six-by-six color filter set
VeriFlex™ Zones ^[2]	Three VeriFlex™ Zones	Six VeriFlex™ Zones
TaqMan™ Array Card block	—	Available
Security, Auditing, and E-signature (SAE) v2.0 module	—	Available ^[3]

[1] See “System dyes” on page 16.

[2] See “VeriFlex™ Zones” on page 18.

[3] See Chapter 9, “Use the instrument with the Security, Auditing, and E-signature (SAE) Administrator Console”.

Instrument filters and supported dyes

System dyes

The QuantStudio™ 6 Pro Real-Time PCR System has a coupled five-color filter set.

The QuantStudio™ 7 Pro Real-Time PCR System has a decoupled six-by-six color filter set.


For more information about available spectral dye calibration kits, contact Support.

Peak filter	Color	Filter wavelength (nm) ^[1]		System dyes	Example custom dyes
		Excitation	Emission		
x1-m1	Blue	470 ± 15	520 ± 15	FAM™ dye, SYBR GREEN™ dye	SYT09
x2-m2	Green	520 ± 10	558 ± 11	VIC™ dye	JOE™ dye, HEX™ dye, TET™ dye ^[2]
x3-m3	Yellow	550 ± 11	586 ± 10	NED™ dye, TAMRA™ dye, ABY™ dye	Cy®3 dye
x4-m4	Orange	580 ± 10	623 ± 14	ROX™ dye, JUN™ dye	Texas Red™ dye
x5-m5	Red	640 ± 10	682 ± 14	MUSTANG PURPLE™ dye, Cy®5	LIZ™ dye
x6-m6	Deep-Red	662 ± 10	711 ± 12	None ^[3]	Cy®5.5 dye

[1] The central wavelengths are the optimized wavelengths.

[2] The HEX™ dye and TET™ dye from Thermo Fisher Scientific fall within the emission wavelength range of the system, therefore they can be added and adapted for use on the instrument.

[3] This filter set currently does not support any dyes supplied by Thermo Fisher Scientific.

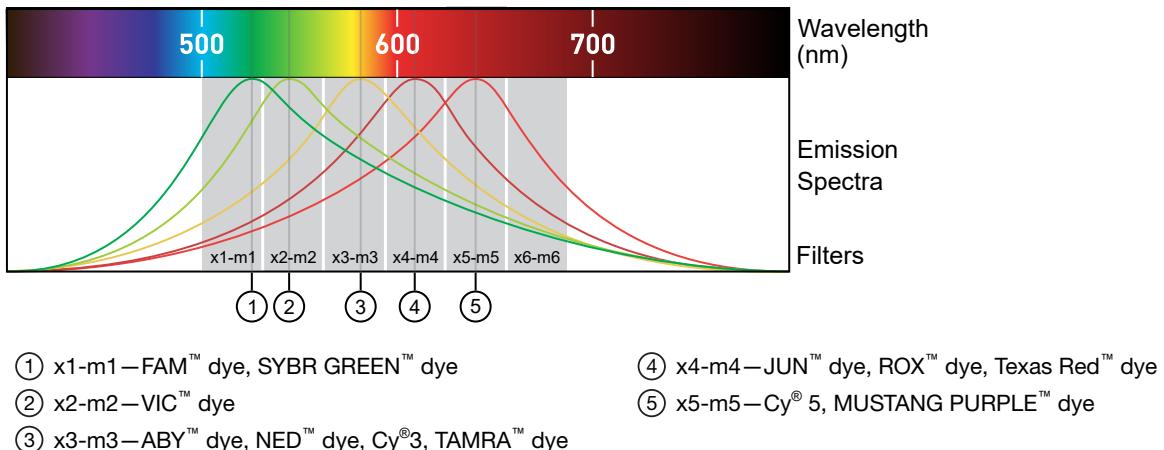

Custom dyes

The instrument can run assays designed with custom dyes. Custom dyes include the following options:

- Dyes that are not supplied by Thermo Fisher Scientific.
- Dyes or formulations of dyes that are not system dyes for the instrument.

To use a custom dye on the instrument, review the following requirements.

- Calibrate the instrument for the custom dye (see “Calibrate custom dyes” on page 113).
- Ensure that the custom dye excites between 455–672 nm and emits between 505–723 nm.
- Select a custom dye that does not overlap with other dyes used in the run (see “System dyes” on page 16).
- The custom dye for use in calibration must be attached to the 5' end of a short DNA oligonucleotide consisting of the first two bases of the probe sequence without a quencher at the 3' end.

Overview of data collection

The instrument collects raw fluorescence data at different points during the PCR cycle, depending on the type of run performed.

When you create a plate file, you can customize the optical filter channels through which the instrument collects data. You can specify a filter channel set for all PCR thermal protocols and for melt curve stages.

Run type	Run type	Data collection point
Real-time	<ul style="list-style-type: none"> Standard curve Relative standard curve Comparative C_t ($\Delta\Delta C_t$) Melt curve 	<p>During the thermal cycling protocol.</p> <p>Typical timing is to collect data at each cycle of a PCR stage or continuously during a melt stage.</p>
Post-PCR (endpoint)	<ul style="list-style-type: none"> Genotyping Presence/absence 	<ul style="list-style-type: none"> After thermal cycling is completed. For Presence/Absence and Genotyping analysis, data collection before the PCR cycle is optional, but recommended. (Optional) During the thermal cycling protocol. Collecting data during the run can confirm genotyping results by viewing traces in allelic discrimination plots or viewing genotyping calls at earlier cycles.

VeriFlex™ Zones

VeriFlex™ Zones provide independent temperature zones that offer enhanced functionality and precise control over your real-time PCR runs.

The independent zones are ideal for real-time PCR optimization. They also provide the ability to perform multiple runs at the same time. Unlike standard gradients which give a sigmoidal temperature curve across the columns, blocks with VeriFlex™ Zones help deliver accurate temperatures across every zone.

The QuantStudio™ 6 Pro System has three VeriFlex™ Zones (Figure 1 on page 19). The QuantStudio™ 7 Pro System has six VeriFlex™ Zones (Figure 2 on page 19).

Note: VeriFlex™ Zones are only applicable for the 96-well 0.2-mL block and the 96-well 0.1-mL block.

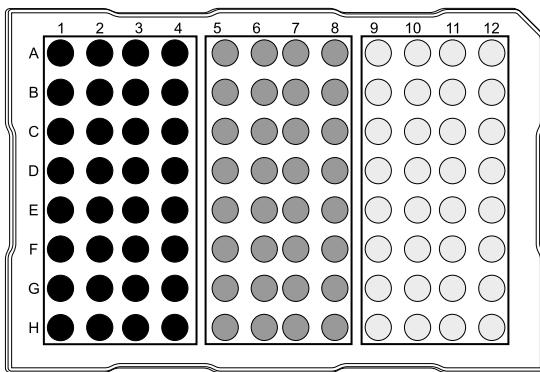


Figure 1 Three VeriFlex™ Zones

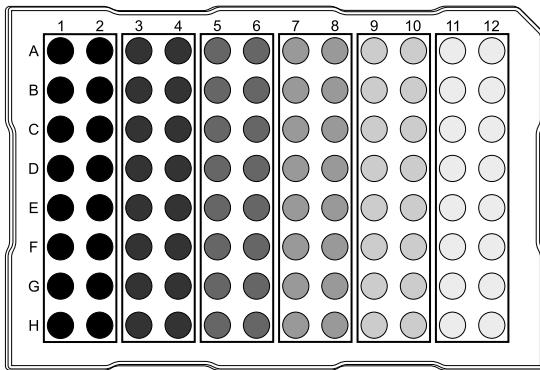


Figure 2 Six VeriFlex™ Zones

Note: Plate files are not compatible between the QuantStudio™ 6 Pro System and the QuantStudio™ 7 Pro System. For more information, see “Compatibility of plate files” on page 29.

AutoDelta settings

AutoDelta enables an incremental increase or decrease in the temperature or time in a PCR cycle.

To use AutoDelta, specify a difference in the time and the temperature. Specify the PCR cycle for this change to be applied.

Parts of the instrument

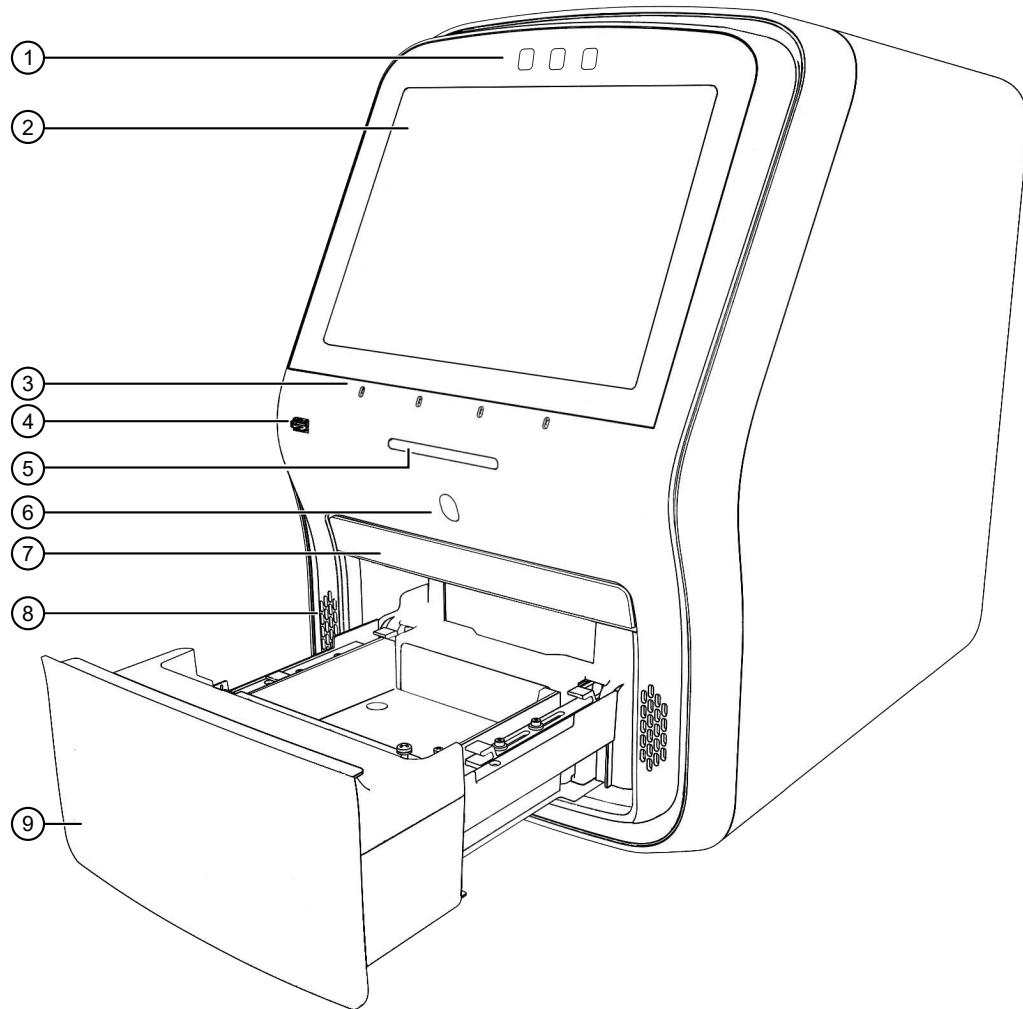


Figure 3 Front of the instrument

① Cameras	⑥ Proximity sensor
② Touchscreen	⑦ Access door for block change
③ Microphones	⑧ Speakers
④ USB port	⑨ Drawer
⑤ Indicator light	

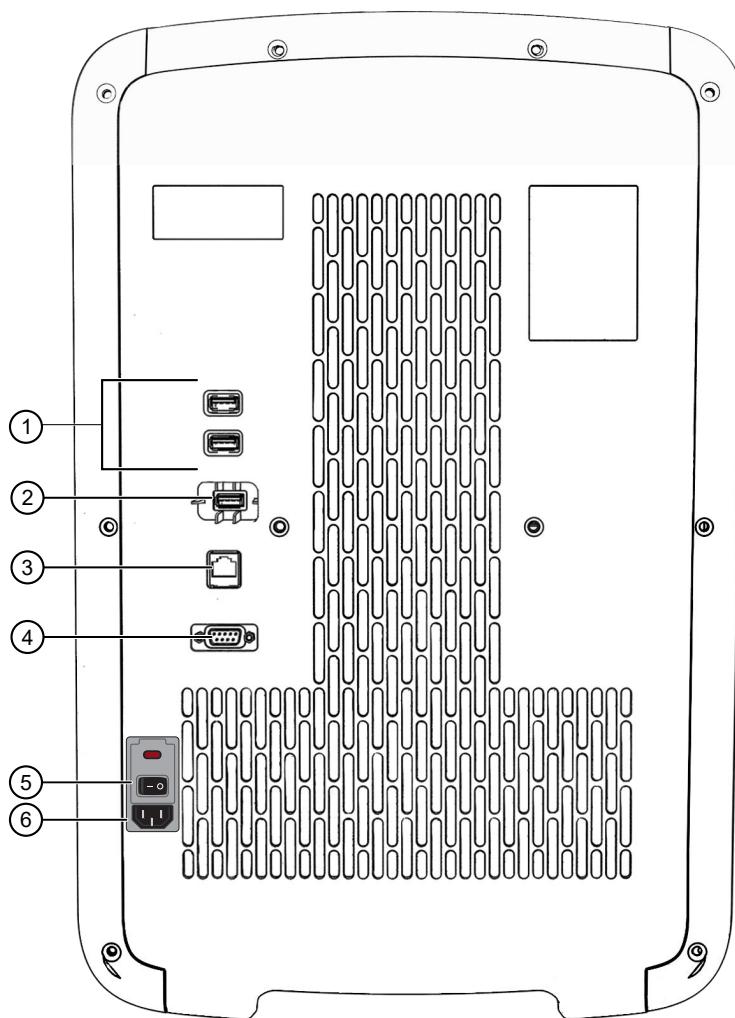


Figure 4 Back of the instrument

① USB ports	④ RS-232 port (service use only)
② Wi-Fi dongle port	⑤ ON/OFF switch
③ Ethernet port	⑥ Power inlet receptacle

Note: The instrument recognizes only one external storage device at a time for data transfer.

The USB ports at the rear of the instrument can be used for external devices, for example, a mouse or a keyboard.

Instrument status indicator

Indicator	Instrument status
Off	Powered off Sleep mode
Blue light changing brightness slowly	Powering on
Blue light changing brightness	Standby
Blue light moving inwards towards middle of instrument	Drawer opening
Blue light moving outwards towards side of instrument	Drawer closing
Blue light moving back and forth	Block is being changed
Blue light on	Instrument is ready to start a run or is performing a run
Amber blinking	Error

Hands-free features

Facial authentication

The instrument allows hands-free facial authentication to sign in a user. The instrument camera detects a face. The instrument will sign in the user after the face is matched to the photo associated with a profile.

Voice commands

The instrument recognizes a set of commands for hands-free operation. For list of supported voice commands, see “[Use voice commands](#)” on page 45.

Voice commands are only available when a user is signed in with a Connect account. Voice commands are not available with a local instrument profile.

Voice commands are only available in select countries.

Voice commands use Amazon™ Alexa™ for Business.

Indicators for the hands-free features

Indicators for facial authentication

Display	Status of facial authentication
The Sign In screen displays  .	Facial authentication for the instrument is enabled.
The Sign In screen does not display  .	<ul style="list-style-type: none">Facial authentication for the instrument is disabled.All users must sign in with a PIN. <p>Note: To enable facial authentication, see “Configure the instrument for the hands-free features” on page 185.</p>

Indicators for voice commands

For the location of the proximity sensor, see Figure 3 on page 20.

Display	Status of voice commands
 (Gray)	The voice command function is not available because the proximity sensor does not detect a person.
 (Gray)	The voice command function is not available if any of these conditions is true: <ul style="list-style-type: none">The voice command function for the instrument is disabled by an administrator.The user is signed in with a local profile.The microphone was disabled by the user and the proximity sensor does not detect a person. The icon will return to a red state when the proximity sensor detects a person.
 (Amber)	The instrument is preparing to accept voice commands. <p>Note: The amber display is an intermediate state and will only appear for several seconds.</p>
 (Blue)	The voice command function is available if all of these conditions are true: <ul style="list-style-type: none">The voice command function for the instrument is enabled by an administrator.The user is signed in with a Connect instrument profile.The proximity sensor detects a person.
 (Red)	The microphone was disabled by the user (see “Enable or disable the microphone” on page 44).

To enable voice commands, see “Configure the instrument for the hands-free features” on page 185.

For more information about linking a local instrument profile to a Connect profile, see “Link the instrument to your Connect account” on page 126 and “If you link when you are signed in to the instrument” on page 129.

Software description

Parts of the home screen



Figure 5 Home screen

① Instrument name	⑦ Block installed
② Instrument calendar	⑧ Current user
③ Microphone	⑨ Connectivity icons
④ Help	⑩ Buttons to access plate files
⑤ Eject	⑪ Instrument settings
⑥ Status dial	

Table 1 Parts of the home screen

Element of the home screen	Function
Instrument name	Set by the administrator to identify the instrument.
Calendar	Schedule time for a run or see runs that are scheduled.
Microphone	Shows the status of the microphone (see “Indicators for voice commands” on page 23 and “Enable or disable the microphone” on page 44).
Help	Access the Help.
Eject	Open or close the drawer to insert or remove a plate or a TaqMan™ Array Card.

Table 1 Parts of the home screen (continued)

Element of the home screen	Function
Status dial	<ul style="list-style-type: none">When the instrument is not in use—Displays Set up run.When the instrument is in use—Displays the sample block temperature, the elapsed run time, and the run status.When the instrument is locked while in use—Displays Locked within the status dial.When a run is complete—Displays Run complete.
Block type	<p>The block that is installed on the instrument.</p> <p>If applicable, the following items are noted:</p> <ul style="list-style-type: none">If there is no block and heated cover set installed.If the block and heated cover are mismatched.
Current user	<p>Displays the current signed-in user.</p> <p>Guest is displayed if no user is signed in.</p>
Connectivity	<p>Only the icons that apply to the connectivity status of the instrument at the time are displayed.</p> <ul style="list-style-type: none">—The instrument is wirelessly connected.—The instrument is connected to a wired network.—A USB drive is plugged into the instrument.—The instrument is connected to a network drive.—The instrument is connected to a Connect account.
Buttons to access plate files ^[1]	<p>Access a plate file to edit it or start a run.</p> <ul style="list-style-type: none">Set up run—Open a system template or a saved plate file.(Load plate file)—Open a system template or a saved plate file.(Run last)—Open the last plate file that was used to start a run. The last plate file is specific to the profile.
Instrument settings	<ul style="list-style-type: none">Edit the file name convention.View the run history, and transfer or delete data files.Select the language.Install, change, or remove the block and the heated cover.Configure the instrument settings.Configure the maintenance and service settings.View notifications.Configure hands-free operation.

^[1] Plate files are applicable to TaqMan™ Array Cards.

Thermo Fisher™ Connect definitions

The terms Connect, Cloud, Thermo Fisher Connect, and Thermo Fisher Cloud are used interchangeably on the instrument touchscreen. The terms Thermo Fisher™ Connect and Connect are used in this document.

The following icons are used interchangeably on the instrument touchscreen:

- 
- 

The  icon is used in this document.

Instrument, desktop software, and Connect platform features

The instrument and the software include the features described below.

Note: System templates cannot be edited. They can be edited, then saved as a separate plate file.

Feature	Instrument	Desktop	Connect
Use as guest (no sign in)	✓	✓	—
Schedule a run	✓	—	✓
Open system templates	✓	✓	✓
Open system templates, edit them, and save them as plate files	✓	✓	✓
Open plate files created on the desktop software	✓	✓	✓
Open plate files created on the Connect platform	✓	✓	✓
Open plate file created on the instrument	✓	✓	✓
Send plate files to the instrument run queue	—	✓	—
Access the plate file associated with last run on the instrument	✓	—	—
Load plate in instrument	✓	—	—
Start run	✓	—	—
View real-time data during a run	✓	—	✓
View results after a run is complete	✓ (Limited)	✓	✓
See instrument status (run in progress)	✓	—	✓
Send email notifications	✓	—	—
Check calibration status	✓	✓	✓
Analyze and export results	✓	✓	✓

(continued)

Feature	Instrument	Desktop	Connect
View analysis of results	—	✓	✓
Set calibration reminders	✓	—	✓
Review exported calibration or RNase P verification results	—	✓	✓ (RNase P only)

Overview of file locations and files

Note: The term plate file is applicable to both plates and TaqMan™ Array Cards.

File locations

File location	Description and function
Run queue	Plate files sent from the desktop software.
Public	<ul style="list-style-type: none"> Plate files saved by guest users are stored in this folder. Data files from runs started by guest users are stored in this folder.
Templates	<ul style="list-style-type: none"> System templates are stored in this folder. System templates cannot be edited. They must be saved as a separate plate file and stored in a different folder. Data files cannot be saved in this folder.
Connect	<ul style="list-style-type: none"> Plate files can be retrieved from a Connect profile. Data files can be saved to a Connect profile. You must have a Connect profile.
Network drive	<ul style="list-style-type: none"> Plate files can be retrieved from a network drive. Data files can be saved to a network drive. The instrument must be connected to a network or a computer.
USB drive	<ul style="list-style-type: none"> Plate files can be retrieved from a USB drive. Data files can be saved to a USB drive. A USB drive must be inserted into the instrument.
My instrument	<ul style="list-style-type: none"> This location is only available to a user who is signed in. Guest users cannot access this location to retrieve plate files or save data files. This location is specific to a user. Files are not accessible to any other users, except for administrators. Plate files can be retrieved from this location. Data files are always saved to this location.

(continued)

File location	Description and function
Post-read	<ul style="list-style-type: none"> In endpoint PCR, the pre-read data is saved in this folder for post-read analysis.

Files

IMPORTANT! Plate files and data files are not included in a back-up of the instrument and are not restored when a backup of the instrument is restored.

Plate files and data files must be transferred separately. See Chapter 6, “View and manage files”. System templates are included in the back-up and are restored when a backup of the instrument is restored.

All file types are applicable to both plates and TaqMan™ Array Cards.

Item	Description and function	Supported in
System template	<ul style="list-style-type: none"> A non-editable template. Accessed from  (Load plate file) and the Templates folder (only system templates are stored in this folder). Accessed from the Set up run button. Is saved as a separate plate file if it is edited. Contains the thermal protocol for the real-time PCR. Does not contain the sample information. Does not contain the target information. Sample information and target information must be added (saved as a separate plate file). The sample information and target information can be added on the desktop software or on the Connect platform after the run is complete. 	Instrument Desktop Connect
Plate File From  (Load plate file)	<ul style="list-style-type: none"> A file that has been saved from a system template. Can be edited and saved. Can be edited and saved as a separate plate file. Contains the thermal protocol for the real-time PCR. Can contain the sample information. Can contain the target information. Sample information and target information must be added if not present. The sample information and target information can be added on the desktop software or on the Connect platform after the run is complete. 	Instrument Desktop Connect

(continued)

Item	Description and function	Supported in
Plate File From  (Run last)	<ul style="list-style-type: none">The plate file associated with the last run for the profile.Can be edited and saved.Can be edited and saved as a separate plate file.Contains the thermal protocol for the real-time PCR.Contains the sample information. The sample information can be edited on the desktop software or on the Connect platform after the run is complete.Contains the target information. The target information can be edited on the desktop software or on the Connect platform after the run is complete.Sample information and target information can be edited on the instrument.If the sample information and the target information for the last run were added on the desktop software or on Connect after the run was complete, they will not be included on the plate file that is on the instrument.	Instrument
Data file	<ul style="list-style-type: none">Contains the real-time PCR data.Can be viewed on the instrument only immediately after the run is complete.Transfer to Connect or to the desktop software to view and analyze data.	Instrument (limited) Desktop Connect

Compatibility of plate files

The plate files are specific to the instrument.

A plate file for a QuantStudio™ 6 Pro System can only be opened on this instrument. It cannot be opened on a QuantStudio™ 7 Pro System.

A plate file for a QuantStudio™ 7 Pro System can only be opened on this instrument. It cannot be opened on a QuantStudio™ 6 Pro System.

Components of the files

Component	Instrument	Desktop	Connect
Properties	<ul style="list-style-type: none"> • Data file name • Reagent information <ul style="list-style-type: none"> – Name – Type – Lot number – Barcode – Catalog number – Expiry date • Plate barcode or TaqMan™ Array Card barcode • Location to send data after the run is complete 	<ul style="list-style-type: none"> • Experiment file name • Plate barcode or TaqMan™ Array Card barcode • User name • Instrument type • Block type • Analysis module • Chemistry (reagent information) • Run mode • Comments • (<i>Connect platform only</i>) Notifications 	
Method	Thermal cycling conditions <ul style="list-style-type: none"> • Temperature • Time • Cover temperature • Reaction volume • Data collection points • Ramp rates • VeriFlex™ Zones • Auto Delta settings 	Thermal cycling conditions	
Plate	Sample names Targets SNP assays	Define and assign samples, targets or SNP assays, and tasks in the Plate setup screen.	
Run	Monitor a run in progress View time remaining, temperature, method, and amplification plots Pause, resume, or stop a run	Monitor a run in progress View time remaining	Monitor a run in progress (link to Connect Instrument Details) View time remaining, temperature, method, and amplification plots
Results	Not applicable	Review plots	
Export	Export results Data files are transferred to the desktop software or the Connect platform for analysis	Export results	




Third-party software

Before installing third-party software on the computer running the desktop software, confirm that the third-party software will not do the following:

- Restrict Ethernet communication.
- Interfere with instrument or computer operation.

Network connection options

You can connect an instrument to a network or computer in the configurations listed below. For more information, see Appendix B, “Connect the instrument to a network”.

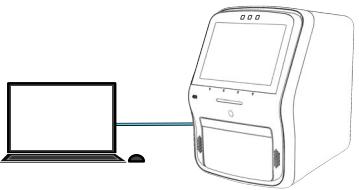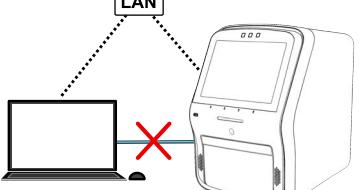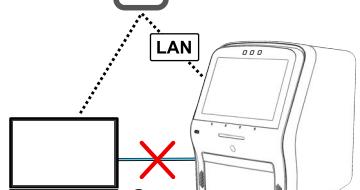Direct connection	Local area network (LAN) connection	Thermo Fisher™ Connect connection
wired	wired or wireless	wired or wireless
		

Types of runs

Standard curve run

A standard curve run determines the absolute target quantity in samples.

1. The software measures amplification of the target in a standard dilution series and in test samples.
2. The software generates a standard curve using data from the standard dilution series.
3. The software uses the standard curve to interpolate the absolute quantity of target in the test samples.

Relative standard curve run

A relative standard curve run determines the relative target quantity in samples.

1. The software measures amplification of the target of interest and of an endogenous control target in a standard dilution series, in a reference (calibrator) sample, and in test samples.

The endogenous control is a target that is expressed equally in all samples; examples of endogenous controls are β -actin, GAPDH, and 18S ribosomal RNA. The software can algorithmically incorporate multiple endogenous control targets in relative quantification calculations.

The reference sample is used as the basis for relative quantification results (or 1 \times sample). For example, in a study of drug effects on gene expression, an untreated control is an appropriate reference sample.

2. The software generates standard curves for the target of interest and the endogenous control using data from the corresponding standard dilution series.
3. The software uses the standard curves to interpolate the quantities of the target of interest and the endogenous control in each sample. The target quantity in each sample is then normalized to the sample's endogenous control quantity.
4. To determine the relative quantity of the target in test samples, the software divides the normalized target quantity in the sample by the normalized target quantity in the reference sample.

Comparative C_t ($\Delta\Delta C_t$) run

A comparative C_t ($\Delta\Delta C_t$) run determines the relative target quantity in samples.

1. The software measures amplification of the target of interest and of an endogenous control target in a reference (calibrator) sample and in test samples.

The endogenous control is a target that is expressed equally in all samples; examples of endogenous controls are β -actin, GAPDH, and 18S ribosomal RNA. The software can algorithmically incorporate multiple endogenous control targets in relative quantification calculations.

The reference sample is used as the basis for relative quantification results (or 1 \times sample). For example, in a study of drug effects on gene expression, an untreated control is an appropriate reference sample.

2. The measurements for the target of interest are normalized to the endogenous control.
3. To determine the relative quantity of the target in test samples, the software compares the normalized ΔC_q (ΔC_t or ΔC_{rt}) for the sample to the normalized ΔC_q (ΔC_t or ΔC_{rt}) for the reference sample.

Genotyping run

A genotyping run detects the single nucleotide polymorphism (SNP) variants of a target nucleic acid sequence.

Genotyping runs use preformulated TaqMan™ SNP Genotyping Assays that include the following components:

- Two sequence-specific primers for amplification of sequences containing the SNP of interest
- Two allele-specific TaqMan™ probes for Allele 1 and Allele 2

1. The software normalizes the fluorescence of the reporter dyes to the fluorescence of the passive reference dye in each well.
2. The software plots the normalized reporter dye signal of each sample well on an Allelic Discrimination Plot, which contrasts the reporter dye intensities of the allele-specific probes.
3. The software algorithmically clusters the sample data, and assigns a genotype call to the samples of each cluster according to its position on the plot.

Presence/absence run

Presence/absence run determines the presence or absence of a target nucleic acid sequence in a sample.

The software calls the target present or absent based on an algorithmically determined call threshold. (The call threshold is different from the C_t threshold; the C_t threshold is not used to make calls.)

Melt curve run

Melt curve run determines the melting temperature (T_m) of the amplification products of a PCR that used intercalating dyes.

In the software, melt curve analysis is included in the default run method for any run type that uses intercalating dyes.

1. The software plots a melt curve based on the fluorescence of the dye with respect to change in temperature.
2. Using the melt curve, the software calculates the melting temperature (T_m).

System template file names

The first part of the file name varies, based on the instrument (QuantStudio™ 6 Pro System or QuantStudio™ 7 Pro System).

The instrument will display the system templates applicable to the block that is installed.

System template file names for the 96-well 0.2-mL block

System template name	Run type
QS7Pro-96-Well-0-2ml_GT_Fast.edt	<ul style="list-style-type: none"> • Genotyping run • 96-well, 0.2-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-2ml_GT_Std.edt	<ul style="list-style-type: none"> • Genotyping run • 96-well, 0.2-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-2ml_Melt_Fast.edt	<ul style="list-style-type: none"> • Melt curve run (no PCR stage) • 96-well, 0.2-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-2ml_Melt_Std.edt	<ul style="list-style-type: none"> • Melt curve run (no PCR stage) • 96-well, 0.2-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-2ml_PA_Fast.edt	<ul style="list-style-type: none"> • Presence/absence run • 96-well, 0.2-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-2ml_PA_Std.edt	<ul style="list-style-type: none"> • Presence/absence run • 96-well, 0.2-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-2ml_PCR_Melt_Fast.edt	<ul style="list-style-type: none"> • Melt curve run (with PCR stage) • 96-well, 0.2-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-2ml_PCR_Melt_Std.edt	<ul style="list-style-type: none"> • Melt curve run (with PCR stage) • 96-well, 0.2-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-2ml_RQ_Fast.edt	<ul style="list-style-type: none"> • Comparative C_t run • 96-well, 0.2-mL Standard plate • Fast cycling mode

(continued)

System template name	Run type
QS7Pro-96-Well-0-2ml_RQ_Std.edt	<ul style="list-style-type: none"> Comparative C_t run 96-well, 0.2-mL Standard plate Standard cycling mode
QS7Pro-96-Well-0-2ml_RSC_Fast.edt	<ul style="list-style-type: none"> Relative standard curve run 96-well, 0.2-mL Standard plate Fast cycling mode
QS7Pro-96-Well-0-2ml_RSC_Std.edt	<ul style="list-style-type: none"> Relative standard curve run 96-well, 0.2-mL Standard plate Standard cycling mode
QS7Pro-96-Well-0-2ml_SC_Fast.edt	<ul style="list-style-type: none"> Standard curve run 96-well, 0.2-mL Standard plate Fast cycling mode
QS7Pro-96-Well-0-2ml_SC_Std.edt	<ul style="list-style-type: none"> Standard curve run 96-well, 0.2-mL Standard plate Standard cycling mode

System template file names for the 96-well 0.1-mL block

System template name	Run type
QS7Pro-96-Well-0-1ml_GT_Fast.edt	<ul style="list-style-type: none"> Genotyping run 96-well, 0.1-mL Standard plate Fast cycling mode
QS7Pro-96-Well-0-1ml_GT_Std.edt	<ul style="list-style-type: none"> Genotyping run 96-well, 0.1-mL Standard plate Standard cycling mode
QS7Pro-96-Well-0-1ml_Melt_Fast.edt	<ul style="list-style-type: none"> Melt curve run (no PCR stage) 96-well, 0.1-mL Standard plate Fast cycling mode
QS7Pro-96-Well-0-1ml_Melt_Std.edt	<ul style="list-style-type: none"> Melt curve run (no PCR stage) 96-well, 0.1-mL Standard plate Standard cycling mode

(continued)

System template name	Run type
QS7Pro-96-Well-0-1ml_PA_Fast.edt	<ul style="list-style-type: none"> • Presence/absence run • 96-well, 0.1-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-1ml_PA_Std.edt	<ul style="list-style-type: none"> • Presence/absence run • 96-well, 0.1-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-1ml_PCR_Melt_Fast.edt	<ul style="list-style-type: none"> • Melt curve run (with PCR stage) • 96-well, 0.1-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-1ml_PCR_Melt_Std.edt	<ul style="list-style-type: none"> • Melt curve run (with PCR stage) • 96-well, 0.1-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-1ml_RQ_Fast.edt	<ul style="list-style-type: none"> • Comparative C_t run • 96-well, 0.1-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-1ml_RQ_Std.edt	<ul style="list-style-type: none"> • Comparative C_t run • 96-well, 0.1-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-1ml_RSC_Fast.edt	<ul style="list-style-type: none"> • Relative standard curve run • 96-well, 0.1-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-1ml_RSC_Std.edt	<ul style="list-style-type: none"> • Relative standard curve run • 96-well, 0.1-mL Standard plate • Standard cycling mode
QS7Pro-96-Well-0-1ml_SC_Fast.edt	<ul style="list-style-type: none"> • Standard curve run • 96-well, 0.1-mL Standard plate • Fast cycling mode
QS7Pro-96-Well-0-1ml_SC_Std.edt	<ul style="list-style-type: none"> • Standard curve run • 96-well, 0.1-mL Standard plate • Standard cycling mode

System template file names for the 384-well block

System template name	Run type
QS7Pro-384-Well-GT-Fast.edt	<ul style="list-style-type: none">• Genotyping run• 384-well plate• Fast cycling mode
QS7Pro-384-Well-GT-Std.edt	<ul style="list-style-type: none">• Genotyping run• 384-well plate• Standard cycling mode
QS7Pro-384-Well-Melt-Fast.edt	<ul style="list-style-type: none">• Melt curve run (no PCR stage)• 384-well plate• Fast cycling mode
QS7Pro-384-Well-Melt-Std.edt	<ul style="list-style-type: none">• Melt curve run (no PCR stage)• 384-well plate• Standard cycling mode
QS7Pro-384-Well-PA-Fast.edt	<ul style="list-style-type: none">• Presence/absence run• 384-well plate• Fast cycling mode
QS7Pro-384-Well-PA-Std.edt	<ul style="list-style-type: none">• Presence/absence run• 384-well plate• Standard cycling mode
QS7Pro-384-Well-PCR-Melt-Fast.edt	<ul style="list-style-type: none">• Melt curve run (with PCR stage)• 384-well plate• Fast cycling mode
QS7Pro-384-Well-PCR-Melt-Std.edt	<ul style="list-style-type: none">• Melt curve run (with PCR stage)• 384-well plate• Standard cycling mode
QS7Pro-384-Well-RQ-Fast.edt	<ul style="list-style-type: none">• Comparative C_t run• 384-well plate• Fast cycling mode
QS7Pro-384-Well-RQ-Std.edt	<ul style="list-style-type: none">• Comparative C_t run• 384-well plate• Standard cycling mode

(continued)

System template name	Run type
QS7Pro-384-Well-RSC-Fast.edt	<ul style="list-style-type: none"> Relative standard curve run 384-well plate Fast cycling mode
QS7Pro-384-Well-RSC-Std.edt	<ul style="list-style-type: none"> Relative standard curve run 384-well plate Standard cycling mode
QS7Pro-384-Well-SC-Fast.edt	<ul style="list-style-type: none"> Standard curve run 384-well plate Fast cycling mode
QS7Pro-384-Well-SC-Std.edt	<ul style="list-style-type: none"> Standard curve run 384-well plate Standard cycling mode

System template file names for the TaqMan™ Array Card block

Note: The TaqMan™ Array Card is only available for the QuantStudio™ 7 Pro System.

System template name	Run type
QS7Pro-TAC-PA-Fast.edt	<ul style="list-style-type: none"> Presence/absence run TaqMan™ Array Card Fast cycling mode
QS7Pro-TAC-PA-Std.edt	<ul style="list-style-type: none"> Presence/absence run TaqMan™ Array Card Standard cycling mode
QS7Pro-TAC-RQ-Fast.edt	<ul style="list-style-type: none"> Comparative C_t run TaqMan™ Array Card Fast cycling mode
QS7Pro-TAC-RQ-Std.edt	<ul style="list-style-type: none"> Comparative C_t run TaqMan™ Array Card Standard cycling mode
QS7Pro-TAC-SC-Fast.edt	<ul style="list-style-type: none"> Standard curve run TaqMan™ Array Card Fast cycling mode

(continued)

System template name	Run type
QS7Pro-TAC-SC-Std.edt	<ul style="list-style-type: none">• Standard curve run• TaqMan™ Array Card• Standard cycling mode
QS7Pro-TAC-RQ-Melt-Fast.edt	<ul style="list-style-type: none">• Comparative C_t run with melt curve• TaqMan™ Array Card• Fast cycling mode
QS7Pro-TAC-RQ-Melt-Std.edt	<ul style="list-style-type: none">• Comparative C_t run with melt curve• TaqMan™ Array Card• Standard cycling mode
QS7Pro-TAC-SC-Melt-Fast.edt	<ul style="list-style-type: none">• Standard curve run with melt curve• TaqMan™ Array Card• Fast cycling mode
QS7Pro-TAC-SC-Melt-Std.edt	<ul style="list-style-type: none">• Standard curve run with melt curve• TaqMan™ Array Card• Standard cycling mode

Workflow

Sign in to your instrument profile or your Connect account
(PIN or facial authentication)



Open a plate file
(System template or saved plate file)



(Optional) Edit the plate file



Load the plate into the instrument



Start the run



Unload the plate from the instrument



(Optional) Transfer the data file

(A location to transfer the data file might have been selected before the run. The data file only needs to be transferred manually after the run in order to transfer to a new location.)

■ Precautions for use	41
■ Installation and instrument verification	42
■ Overview of guest access	42
■ Power on the instrument	43
■ Sign in	43
■ Enable or disable the microphone	44
■ Use voice commands	45
■ Sign out	46

Precautions for use



CAUTION! PHYSICAL INJURY HAZARD. Do not remove the instrument cover. There are no components inside the instrument that you can safely service yourself. If you suspect a problem, contact technical support.



CAUTION! FIRE HAZARD. For continued protection against the risk of fire, replace fuses only with listed and certified fuses of the same type and rating as those currently in the instrument.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the sample block temperature can reach 100°C. Allow it to cool to room temperature before handling.



CAUTION! Before using a cleaning or decontamination method other than those recommended by Thermo Fisher Scientific, confirm with Thermo Fisher Scientific that the proposed method will not damage the instrument.



CAUTION! Use flat caps for tubes. Rounded caps can damage the heated cover.

Installation and instrument verification

The instrument is installed by a representative from Thermo Fisher Scientific. The following tasks will be performed during the installation process:

- Unpack the instrument and the computer
- Install the instrument and the computer
- Install QuantStudio™ Design and Analysis Software v2 on the computer
- Perform calibrations and RNase P verification

There is a recommended order to create instrument profiles and link Connect accounts to the instrument (see “Recommended order to set up profiles” on page 123).

Overview of guest access


The instrument can be used without signing in. The following limits are in place if you do not sign in to the instrument:

- All actions are recorded to the guest profile.
- You have access only to the public folder for selecting and storing experiments.
- You cannot transfer data files to Connect.
- Data files can only be transferred to a USB drive or a network drive.

The instrument can be configured so that guest access is not permitted (see “Require sign-in” on page 122).

Use the instrument as a guest

If the **Guest** button is not visible on the **Sign In** screen, guest access has been disabled. An administrator must enable guest access (see “Require sign-in” on page 122).

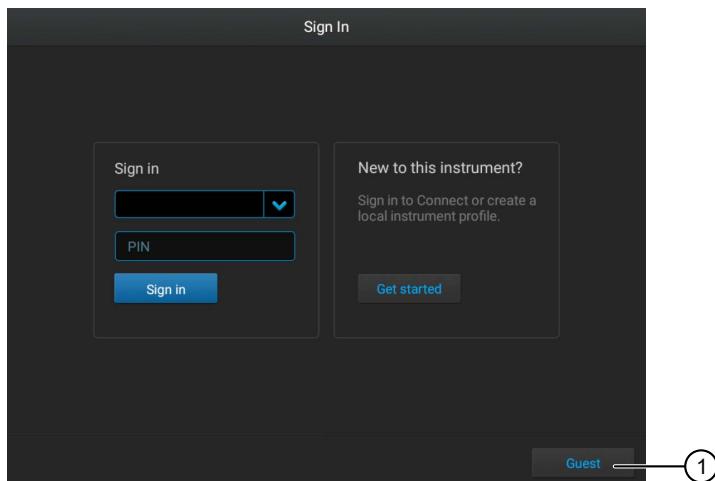

① Guest button

In the **Sign-In** screen, tap **Guest**.

Power on the instrument

If left unattended, the instrument automatically enters sleep mode to conserve power. The default time that an instrument is inactive before going into sleep mode is 15 minutes.

1. Tap anywhere on the touchscreen to determine if the instrument is in sleep mode.
If the screen becomes active, the instrument is already powered on.
2. If the home screen does not display, power on the instrument by pressing the ON/OFF switch on the rear panel.

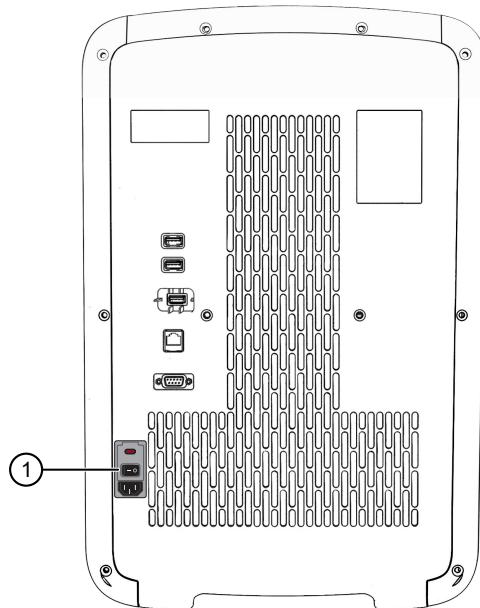


Figure 6 Location of the ON/OFF switch

① ON/OFF switch

Note: The length of time that the instrument is inactive before going into sleep mode can be customized (see “Enable and edit sleep mode” on page 163).

Sign in

Sign in with a PIN

Create an instrument profile before signing into the instrument (see “Create a new local instrument profile” on page 127 and “Link the instrument to your Connect account” on page 126).

1. In the **Sign In** screen, tap the **PIN** field, then enter your PIN.
If your instrument profile is not the default that is displayed, tap **▼**, then select your instrument profile.
2. Enter your PIN, then tap **Enter**.

You are signed in and the home screen is displayed.

See “Parts of the home screen” on page 24.

Sign in with facial authentication

For information on indicators for facial authentication, see “Indicators for facial authentication” on page 23.

For lighting guidelines, see “Light guidelines” on page 192.

Create an instrument profile before signing into the instrument (see “Create a new local instrument profile” on page 127 or “Link the instrument to your Connect account” on page 126). The instrument profile must be set up with facial authentication (see “Add facial authentication to an existing instrument profile” on page 132).

The instrument must accept facial authentication (see “Configure the instrument for the hands-free features” on page 185).

The instrument must have the sign-in screen displayed. Sign out the current user to return to the sign-in screen (see “Sign out” on page 46).

1. Stand in front of the instrument.

Center yourself in front of the instrument. The ideal position is approximately 27 inches (69 cm) from the camera.

Note: You do not need to position yourself as closely to the instrument to sign in as when you took the photos to set up facial authentication.

2. Wait while the instrument searches for a match with the photos associated with profiles.

3. *(Optional)* Reposition yourself if the instrument indicates that the face is out of position.

Ensure that you are centered in front of the instrument. Move away from the instrument and towards the instrument to find the correct position.

The user is signed into the instrument and the home screen is displayed.

Enable or disable the microphone

The microphone can be disabled if the voice command function is available. A blue microphone icon indicates that the voice command function is available.

The microphone cannot be manually disabled if the voice command function is already not available. A gray microphone icon indicates that the voice command function is unavailable.

To enable or disable the voice command function for the instrument, see “Configure the instrument for the hands-free features” on page 185.

For a description of the indicators, see “Indicators for voice commands” on page 23.

Tap  **(Microphone)**.

The microphone icon will change to red with a bar to indicate that it is disabled ().

If the proximity sensor does not detect a person, the microphone icon will change to gray with a bar (ⓘ). It will change back to red with a bar when the proximity sensor detects a person.

Use voice commands

The instrument must accept voice commands (see “Configure the instrument for the hands-free features” on page 185).

A Connect profile must be used. For information about linking a local instrument profile to a Connect profile, see “Link a local profile to a Connect profile” on page 127.

Voice commands are not available when the instrument is used as a guest. Voice commands are not available with a local instrument profile.

The instrument must display a screen where a voice command is supported.

For information on indicators for voice commands, see “Indicators for voice commands” on page 23.

Say “Alexa™, open pro instrument”, wait for the instrument audio prompt and, then issue the voice command.

The instrument audio prompt depends on the voice command that was issued.

Voice command	Instrument audio output	Instrument action
Home	—	The home screen is displayed.
Open door	—	Instrument drawer opens.
Shut door	—	Instrument drawer closes.
Begin Run	“Are you sure you want to start the run? Yes or no.”	If yes, the instrument starts the run. If no, the Plate properties screen is displayed.
End run	“Are you sure you want to stop the run? Yes or no.”	If yes, the instrument stops the run and displays the Plate properties screen. If no, the run continues.
Lock screen	—	The instrument locks the screen. The status dial displays Locked .
Change Block ^[1]	—	If the instrument has a block installed, the Change Block screen is displayed. If there is no block in the instrument, the Install Block screen is displayed.
Instrument scheduler ^[1]	—	The instrument schedule is displayed.
Help	—	The help topic associated with the current page on the screen is displayed.
Smart support ^[1]	—	The Smart Help screen is displayed.
Calibrations	—	The Calibration screen is displayed.

^[1] Unsaved changes are discarded after the instrument action is completed.

Sign out

1. In the home screen, tap  **(Profile)**.
2. Tap **Sign out**, then tap **Yes** to confirm.

You are signed out and the **Sign In** screen is displayed.

■ Options to run a plate	47
■ Overview of customized data export	50
■ (Optional) Scan a barcode using a barcode scanner	51
■ Load a plate into the instrument	51
■ Load a TaqMan™ Array Card into the instrument	52
■ Unload a plate or a TaqMan™ Array Card from the instrument	52

Options to run a plate

Create and run a plate from a system template or a saved plate file

A system template is a plate file that is pre-loaded on the instrument and not editable. It can be edited and saved as a new custom plate file.

Plate files can be edited in Connect or the desktop software.

The **Run Queue** folder contains plate files sent to the instrument from the desktop software.

1. In the home screen, tap one of the following options:

-  **(Load plate file)**
- **Set up run**

2. Tap the location of the system template or the plate file in the left column.

-  **Run Queue**
-  **Public**
-  **My Instrument**

Note: **My instrument** is not available if you are using the instrument as a guest. You must be signed in.

-  **Templates**
-  **Connect**

Note: You must be signed in with a Connect profile.

-  **Network Drive**

-  **USB Drive**

3. Tap the template file name in the right column.

The **Plate Properties** screen is displayed.

4. *(Optional)* In the **Plate Properties** screen, edit the **Data file name**.

The default file name is the template name appended with the date and time.

To edit the default file name, see “Manage the file name convention of the data file” on page 185.

5. *(Optional)* Use the barcode scanner.

See “*(Optional)* Scan a barcode using a barcode scanner” on page 51.

6. Tap the location to send the data file.

-  **Connect platform**

Note: You must be signed in with a Connect profile.

-  **USB drive**
-  **Network drive**
-  **This instrument**

Note: This option is not available if you are using the instrument as a guest. You must be signed in.

The connection status of each location is displayed.

If the connection between the instrument and the location that was selected to save the data file is interrupted, the instrument completes the run. The data file is saved on the instrument. The data file must be transferred to another location manually (see “Transfer data files at a later time” on page 81).

7. *(Optional)* Select the **Export run results with** checkbox.

For more information about this setting, see “Overview of customized data export” on page 50.

To import a different export setting, see “Import an auto export setting” on page 55.

8. *(Optional)* Edit the run method (see page 57).

9. *(Optional)* Edit the plate (see page 65).

10. Load a plate into the instrument (see page 51).

11. Tap **Start run**, then confirm that a plate was loaded.

The confirmation page can be disabled (see “Enable or disable the reminder to insert a plate” on page 186).

The status dial is displayed. The status dial contains the following information:

- The block temperature
- The elapsed time of the run
- The status of the run



Figure 7 Home screen during a run

For options during a run, see Chapter 5, “Options for a run”.

Repeat the last run

The last run is associated with the profile. It is not the last run for the instrument.

1. In the home screen, tap  **(Run last)**.

The **Plate Properties** screen is displayed.

2. *(Optional)* In the **Plate Properties** screen, edit the **Data file name**.

The default file name is the template name appended with the current date and time.

To edit the default file name, see “Manage the file name convention of the data file” on page 185.

3. *(Optional)* Use the barcode scanner.

See “*(Optional)* Scan a barcode using a barcode scanner” on page 51.

4. Tap the location to send the data.

-  **Connect platform**

Note: You must be signed in with a Connect profile.

-  **USB drive**

-  **Network drive**

-  **This instrument**

Note: This option is not available if you are using the instrument as a guest. You must be signed in.

The connection status of each location is displayed.

If the connection between the instrument and the location that was selected to save the data file is interrupted, the instrument completes the run. The data file is saved on the instrument. The data file must be transferred to another location manually (see “Transfer data files at a later time” on page 81).

5. *(Optional)* Select the **Export run results with** checkbox.

For more information about this setting, see “Overview of customized data export” on page 50.

To import a different export setting, see “Import an auto export setting” on page 55.

6. *(Optional)* Edit the run method (see page 57).

7. *(Optional)* Edit the plate (see page 65).

8. Load a plate into the instrument (see page 51).

9. Tap **Start run**, then confirm that a plate was loaded.

The confirmation page can be disabled (see “Enable or disable the reminder to insert a plate” on page 186).

The status dial is displayed. The status dial contains the following information:

- The block temperature
- The elapsed time of the run
- The status of the run

See Figure 7 on page 49.

For options during a run, see Chapter 5, “Options for a run”.

Overview of customized data export

The instrument can perform the primary analysis and the secondary analysis of the data, then export the results.

The instrument has a default export setting.

The export settings can be customized. The instrument will perform the primary and secondary analysis according to the export setting file.

The export settings are customized in the QuantStudio™ Design and Analysis Software v2, then transferred to the instrument. The export settings are in a JSON file format.

The export settings cannot be created or edited on the instrument.

(Optional) Scan a barcode using a barcode scanner

The instrument is compatible with an optional Handheld Barcode Scanner (Cat. No. [4488442](#), purchased separately). The barcode scanner reads Code 128 (alphanumeric), which supports 128 ASCII character barcodes.

For safety information, see “Laser safety” on page 224.

1. Tap the **Barcode** field.
2. Hold the scanner 20–30 cm away from a plate or container label and aim at the center of the barcode, then press the trigger.
3. Slowly move the scanning beam across the barcode until the scanner emits a high-pitched tone.

When the scanner scans a barcode, it automatically transmits the following information:

- Transmits the alphanumeric equivalent of the barcode to the **Barcode** field.
- Transmits other reagent information.

For more information about the hand-held barcode scanner, see the user documentation provided with the barcode scanner.

Load a plate into the instrument

IMPORTANT! The instrument should be used by trained operators who have been warned of the moving parts hazard.



CAUTION! Use flat caps for 0.2-mL tubes and 0.1-mL tubes. Rounded caps can damage the heated cover.


1. Tap  **(Eject)** to open the drawer.
2. Load the plate onto the block.
 - Well A1 of the plate is in the top-left corner of the block.
 - If the plate has a barcode, the barcode faces the front of the instrument.
3. Tap  **(Eject)** to close the drawer.

Load a TaqMan™ Array Card into the instrument

IMPORTANT! The instrument should be used by trained operators who have been warned of the moving parts hazard.

1. Tap  **(Eject)** to open the drawer.
2. Load the plate onto the block.
 - Well A1 of the array card is in the top-left corner of the block.
 - If the array card has a barcode, the barcode faces the front of the instrument.

The block has guides for the position of the array card.

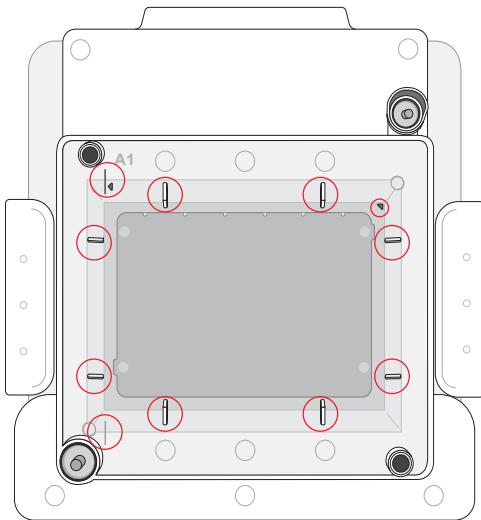

The circles indicate the guides for the position of the array card.

3. Tap  **(Eject)** to close the drawer.

Unload a plate or a TaqMan™ Array Card from the instrument



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the plate temperature can reach 100°C. Allow it to cool to room temperature before handling.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the TaqMan™ Array Card temperature can reach 100°C. Allow it to cool to room temperature before handling.

The status dial displays **Run complete** when the run is finished. If the automatic sign-out feature is enabled and the set time has elapsed, the instrument will display the **Sign In** screen.

1. Tap  **(Eject)** to open the instrument drawer, then remove the plate.
2. Tap  **(Eject)** to close the instrument drawer.

For information on plate disposal, see “Chemical safety” on page 230 and “Biological hazard safety” on page 231.

Edit a plate before starting a run

■ Overview of a template file with restricted editing	54
■ Options to edit the properties of a plate file	55
■ Options to edit the method of a plate file	56
■ Options to edit the well details of a plate file	62


Overview of a template file with restricted editing

A template file can be set up with restricted editing in the QuantStudio™ Design and Analysis Software v2. If a template file is set up with restricted editing, only certain parts of the template file can be edited. The parts of the template file that cannot be edited are defined by the user in QuantStudio™ Design and Analysis Software v2.

A user cannot set up the restricted editing properties on the instrument.

If a template file has restricted editing, it is noted in the following locations:

- The top of the screen
- Under the **Data file properties** heading
- In the **Data file name** field

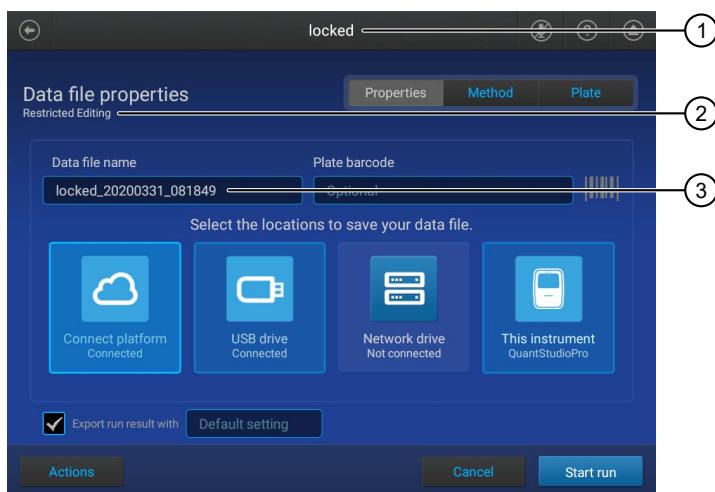

- ① The top of the screen
- ② Under the **Data file properties** heading
- ③ In the **Data file name** field

Options to edit the properties of a plate file

The following files can be edited:

- A system template (see “Create and run a plate from a system template or a saved plate file” on page 47)
A system template that is edited must be saved as a plate file. A system template cannot be overwritten.
- A saved plate file (see “Create and run a plate from a system template or a saved plate file” on page 47)
- The plate file associated with the last run (see “Repeat the last run” on page 49)

The following items can be edited in the **Properties** tab:

- Import the sample setup (see page 55)

Import a sample setup

This feature allows you to import a sample setup file. A sample setup file shows the position of each sample on the plate. The sample setup file is a TXT file or a CSV file.

1. In the **Properties** tab, tap **Actions**, then tap **Import sample setup**.
2. In the **Import Sample Setup** screen, tap the appropriate folder, then tap the appropriate file.

Import an auto export setting

The export settings can be customized. For more information about the export settings, see “Overview of customized data export” on page 50.

The export settings file is in a JSON file format.

The export settings must be customized in the QuantStudio™ Design and Analysis Software v2, then they are imported onto the instrument.

1. In the **Properties** tab, tap the **Export run results with** field.
The **Auto Export Settings** screen is displayed.
2. Tap **Import**.
3. Tap the location of the file in the left pane, then tap the file in the right pane.

Delete an auto export setting

1. In the **Properties** tab, tap the **Export run results with** field.
The **Auto Export Settings** screen is displayed.
2. Tap **Delete**.

Options to edit the method of a plate file

The following files can be edited:

- A system template (see “Create and run a plate from a system template or a saved plate file” on page 47)
A system template that is edited must be saved as a plate file. A system template cannot be overwritten.
- A saved plate file (see “Create and run a plate from a system template or a saved plate file” on page 47)
- The plate file associated with the last run (see “Repeat the last run” on page 49)

The following items can be edited in the **Methods** tab:

- Time and temperature for each PCR step (see page 57)
- Cover temperature and volume (see page 57)
- Number of cycles (see page 57)
- Steps (see page 58)
- Stages (see page 58)
- Data collection points (see page 58)
- VeriFlex™ Zones (see page 59)
- Auto Delta (see page 61)
- Ramp rates (see page 62)
- Pauses (see page 62)

Method elements



① Method tab	⑥ Number of cycles for the stage
② Heated cover temperature	⑦ Temperature for the step
③ Reaction volume	⑧ Hold time for the step
④ Stage of thermal protocol	⑨ Data collection point
⑤ Step within a stage	

Edit the method

1. In the **Method** tab, tap **Edit**.
2. Tap a field, then enter the changes.
 - Cover temperature
 - Volume
 - Temperature for any steps
 - Time for any steps
 - Number of PCR cycles
3. Tap **Save**.
4. Enter a name for the plate file, then tap **Done**.
5. In the **Save Template** screen, select a storage location for the plate file, then tap **Done**.
A confirmation is displayed.
6. Tap **Close**.

Add or remove steps

Add a PCR step or remove a PCR step.

1. In the **Method** tab, tap **Edit** ▶ **Advanced** ▶ **Manage steps**.

Note: Use < or > to navigate through the method if it is not all displayed.

2. Tap **Add step**, then tap + (Add) in the location of the step to be added.
3. Enter the temperature and time, then tap **Enter**.
4. Tap **Remove step**, then tap — (Remove) in the appropriate step.
5. Tap **Done**.

The **Method** tab is displayed with the updated method.

Add or remove stages

Add a PCR stage or remove a PCR stage.

1. In the **Method** tab, tap **Edit** ▶ **Advanced** ▶ **Manage stages**.

Note: Use < or > to navigate through the method if it is not all displayed.

2. Tap **Add Stage**, then tap + (Add) in the location of the stage to be added.
3. Tap the type of stage.
 - **Hold**
 - **Pre-Read**
 - **Post-Read**
 - **Infinite**
 - **PCR**
 - **Melt Curve**

The new stage is added. The time and temperature must be edited as part of the method (see “Edit the method” on page 57).

4. Tap **Remove stage**, then tap — (Remove).
5. Tap **Done**.


Add or remove data collection locations

1. In the **Method** tab, tap **Edit** ▶ **Advanced** ▶ **Data collection location**.

Note: Use < or > to navigate through the method if it is not all displayed.

2. Tap  **(Data collection location)** to designate a step as a data collection location or to remove data collection from the step.

The  **(Data collection location)** button functions as an on-off switch.

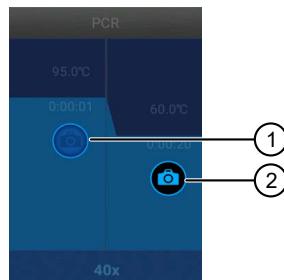

① Data collection is off
② Data collection is on

3. Tap **Done**.

Set up or remove VeriFlex™ Zones

VeriFlex™ Zones are only available for 96-well blocks.

- The QuantStudio™ 6 Pro System has three VeriFlex™ Zones.
- The QuantStudio™ 7 Pro System has six VeriFlex™ Zones.

Note: For more information about compatibility of plate files with VeriFlex™ Zones between the two instruments, see “VeriFlex™ Zones” on page 18.

1. In the **Method** tab, tap **Edit** ▶ **Advanced** ▶ **VeriFlex™ Zones / Auto Delta**.

Note: Use < or > to navigate through the method if it is not all displayed.

Note: The PCR stage provides the option of VeriFlex™ Zones or the Auto Delta feature. Ensure that the **VeriFlex™ Zones** option is selected.



Figure 8 VeriFlex™ Zones and Auto Delta settings

① **VeriFlex™ Zones** is selected

2. Tap the temperature field in the appropriate VeriFlex™ Zone.
3. Enter a temperature, then tap **Enter**.
The temperatures in adjacent zones must have a difference of $\leq 5^{\circ}\text{C}$.
4. Tap **Done**.
A stage with VeriFlex™ Zones is denoted with a  icon.
5. In the **Method** tab, tap  **Remove** **Done** to remove the VeriFlex™ Zones.

Note: The VeriFlex™ Zones setting is removed from the **Method** tab.



① Indicates that this is the **Method** tab where the VeriFlex™ Zones settings can be removed

② Indicates that VeriFlex™ Zones are assigned. Tap to remove the VeriFlex™ Zones.

Set up or remove the Auto Delta feature

Note: The Auto Delta feature is only available in the PCR stage.

1. In the **Method** tab, tap **Edit** ▶ **Advanced** ▶ **VeriFlex™ Zones / Auto Delta**.

Note: Use < or > to navigate through the method if it is not all displayed.

2. Tap  (**Edit**) in one of the PCR stages.

3. Tap **Auto Delta settings**.

See Figure 8 on page 60.

4. Enter the starting cycle to apply Auto Delta settings.

5. Set the **Delta temperature** slider to the + position or the — position, then enter the numerical value for the difference in the temperature.

6. Set the **Delta time** slider to the + position or the — position, then enter the numerical value for the difference in the time.

7. Tap **Done**.

A PCR step with an Auto Delta setting applied to it is denoted with an .

8. In the **Method** tab, tap  ▶ **Remove** ▶ **Done** to remove the Auto Delta setting.

Note: The Auto Delta setting is removed from the **Method** tab.



① Indicates that this is the **Method** tab where Auto Delta settings can be removed.

② Indicates that Auto Delta is assigned. Tap to remove the Auto Delta settings.

Configure ramp rates

The ramp rate is how quickly the temperature changes from one stage or step to another stage or step. The maximum ramp rate for the instrument is 6.5°C/second.

1. In the **Method** tab, tap **Edit** ▶ **Advanced** ▶ **Ramp rates**.

Note: Use < or > to navigate through the method if it is not all displayed.

2. Tap a ramp rate, then edit the value.
3. Tap **Done**.

Add or remove a pause

A pause allows you to open the drawer to add additional reagents to the plate.

1. In the **Method** tab, tap **Edit** ▶ **Advanced** ▶ **Add pause**.

Note: Use < or > to navigate through the method if it is not all displayed.

2. Tap **II (Pause)**.
3. Enter a temperature and the cycle.
A pause in the thermal cycling protocol is denoted with a . The cycle of the pause and the temperature of the pause are also denoted.
4. Tap **II (Pause)** to remove the pause.
5. Tap **Done**.

Options to edit the well details of a plate file

The following files can be edited:

- A system template (see page 47)
A system template that is edited must be saved as a plate file. A system template cannot be overwritten.
- A saved plate file (see “Create and run a plate from a system template or a saved plate file” on page 47)
- The plate file associated with the last run (see “Repeat the last run” on page 49)

The following items can be edited in the **Plate** tab:

- Sample name and sample type (see “Assign or edit a well for a plate” on page 65 and “Assign or edit a well for a TaqMan™ Array Card” on page 66)
- Target name, reporter dye, and quencher dye (see page 66)
- Reagents (see “Add or edit the reagent details for a plate” on page 66 and “Add or edit the reagent details for a TaqMan™ Array Card” on page 68)
- Passive reference dye (see page 68)
- Import an assay information file (AIF; see page 69)


A plate layout file for preplated assays will already be populated with the target name, reporter dye, and quencher dye.

Options to view a plate layout

The following views are available for a plate layout:

- Plate view
 - 
 - Shows each well on the plate
 - Tap to provide information about the sample name, sample type, target name, reporter dye, or quencher dye
 - An individual well
 - Multiple wells (tap each individual well)
 - A row (tap the letter)
 - A column (tap the number)
 - Tap **(Select all)** to select all of the wells in the plate
- Well table view
 - 
 - Shows the well ID, the sample, the target, the reporter, the quencher, the type, and the reagents


The plate view and the well table view can show the sample color or the target color with the **Color By** dropdown list.

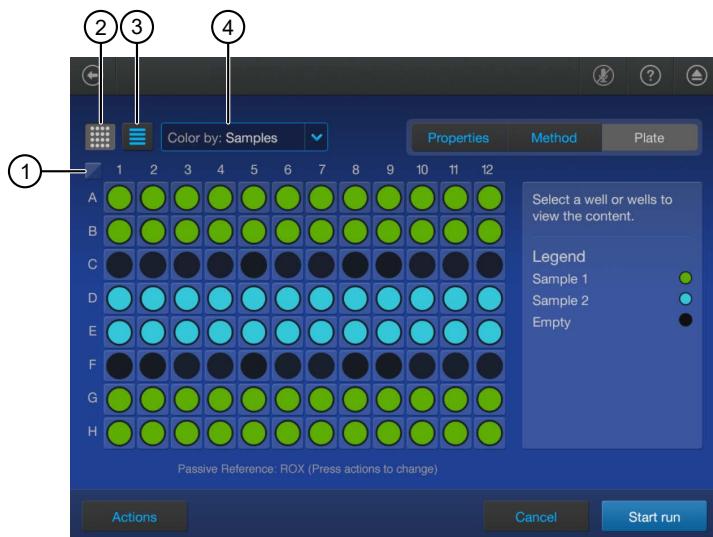

- ① Select all button
- ② Plate view button
- ③ Well table view button
- ④ Color by dropdown list

View a 384-well plate

The **Plate** tab displays a 384-well plate if the 384-well block is installed. The plate can be edited in the 384-well view. You can zoom in to a 96-well section of the 384-well plate in order to edit the plate.

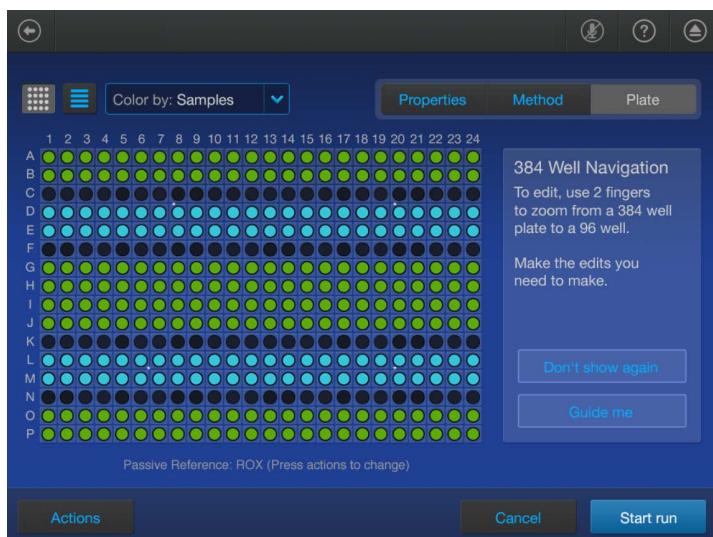

1. In the region of the 384-well plate that you want to zoom in on, pinch outwards with two fingers.
2. (Optional) To change the region that you are zoomed in on, use two fingers to drag the plate in the appropriate direction.

The 384-well plate can zoom in on any 96-well section. It is not limited to specific 96-well quadrants.

If you selected a full row or column in the 96-well view, only the set of wells that was visible in the 96-well view is selected. The full row or column of the 384-well plate is not selected.

3. To return to the full 384-well view from the 96-well section, pinch inwards with two fingers.

View a TaqMan™ Array Card

The **Plate** tab displays a TaqMan™ Array Card if the TaqMan™ Array Card block is installed.

The samples and reagents are edited by port. The samples and reagents are not edited by well.

The targets are added by an assay information file. The targets cannot be edited on the instrument touchscreen.



Figure 9 Plate tab for a TaqMan™ Array Card

① Port

Assign or edit a well for a plate

A target or an assay must be defined (see “Add a target or an assay” on page 66).

A sample setup file can be imported instead of assigning or editing a well manually (see “Import a sample setup” on page 55).

1. In the **Plate** tab, tap a well or multiple wells.
2. Tap the **Sample Information** pane, then tap **Samples**.

Note: If no wells are selected, the pane is labeled **Legend**.

3. Enter a sample name.
4. Select a sample type from the dropdown list.

5. Tap **Target/SNP**, then select **Target** or **Assay** from the dropdown list.
6. Tap a row in the table, then tap **Done**.

Add a target or an assay

1. In the **Plate** tab, tap a well or multiple wells.
2. Tap the **Sample Information** pane, then tap **Target/SNP**.
3. Tap a row in the table, then tap **Add New**.
4. Enter the target, then select the reporter dye and the quencher dye from the respective dropdown lists.
5. Tap **Add Target**, then tap **Close**.

Assign the target or the assay to a well (see “Assign or edit a well for a plate” on page 65).

Assign or edit a well for a TaqMan™ Array Card

An assay information file must be imported to assign targets (see “Import an Assay Information File (AIF)” on page 69). Targets cannot be assigned on the instrument touchscreen.

A sample can only be edited by port. It cannot be edited by individual well.

A sample setup file can be imported instead of assigning or editing a well manually (see “Import a sample setup” on page 55).

1. In the **Plate** tab, tap a port.
See Figure 9 on page 65.
2. Tap the **Legend** pane, then tap **Samples**.
3. Enter a sample name.
4. Select a sample type from the dropdown list.
5. Tap **Done**.

Add or edit the reagent details for a plate

1. In the **Plate** tab, tap one of the following items.
 - A single well
 - Multiple wells
 - **(Select all)**
2. Tap the **Sample Information** pane, then tap **Reagents**.

Note: If no wells are selected, the pane is labeled **Legend**.

3. Tap a row in the table, then enter the information.
 - Reagent name
 - Reagent type
 - Lot number
 - SKU (Part number)
 - Expiration date
4. Tap **Save reagent**.
5. Tap **Done**.

Apply the reagents to a plate

Add the reagents before they are applied to a plate (see “Add or edit the reagent details for a plate” on page 66).

1. In the **Plate** tab, tap one of the following items.
 - A single well
 - Multiple wells
 - **(Select all)**
2. Tap the **Sample Information** pane, then tap **Reagents**.

Note: If no wells are selected, the pane is labeled **Legend**.

3. Tap a row in the table to apply it to the plate.
4. Tap **Done**.

Delete the reagent details

1. In the **Plate** tab, tap one of the following items.
 - A single well
 - Multiple wells
 - **(Select all)**
2. Tap the **Sample Information** pane, then tap **Reagents**.

Note: If no wells are selected, the pane is labeled **Legend**.

3. *(Optional)* Tap **Clear all**.

4. Tap a row in the table, then delete the information in each field.
 - Reagent name
 - Reagent type
 - Lot number
 - SKU (Part number)
 - Expiration date
 - Quencher dye
5. Tap **Save reagent**.
6. Tap **Done**.

Add or edit the reagent details for a TaqMan™ Array Card

1. In the **Plate** tab, tap a port.
See Figure 9 on page 65.
2. Tap the **Legend** pane, then tap **Reagents**.
3. Tap a row in the table.
4. Tap **Add new**, then enter the information.
 - Reagent name
 - Reagent type
 - Lot number
 - SKU (Part number)
 - Expiration date
5. Tap **Save reagent**, then tap **Done**.

Apply the reagents to a TaqMan™ Array Card

Add the reagents before they are applied to a plate (see “Add or edit the reagent details for a TaqMan™ Array Card” on page 68).

A reagent can only be applied by port. It cannot be applied by individual well.

1. In the **Plate** tab, tap a port.
See Figure 9 on page 65.
2. Tap the **Legend** pane, then tap **Reagents**.
3. Tap a row in the table to apply it to the port.
4. Tap **Done**.

Edit a passive reference dye

A passive reference dye normalizes the fluorescent reporter signal in real-time PCR.

1. In the **Plate** tab, tap **Actions** ▶ **Edit passive reference dye**.
The **Passive Reference Dye** screen is displayed.
2. Select the passive reference dye.

The passive reference dye is displayed under the plate diagram.

Import an Assay Information File (AIF)

The Assay Information File is provided in TXT format and/or XML and HTML format, depending on the product line and order date. The HTML format is a reference for users. The XML or the TXT format is imported.

For a detailed description of the AIF, see *Understanding Your Shipment* (Pub. No. MAN0017153)

1. In the **Plate** tab, tap **Actions** ▶ **Import AIF file**.
2. In the **Import AIF file** screen, tap the appropriate folder, then tap the appropriate file.

Import a TaqMan™ Assay file

The TaqMan™ Assay file can be imported to the instrument directly from the Thermo Fisher Scientific website. The instrument must be connected to the internet.

1. In the **Plate** tab, tap **Actions** ▶ **Import TaqMan™ Assay files**.
2. On the **Select A Product** screen, select one of the following options.
 - **Fixed TaqMan™ Array Plates**
 - **Custom TaqMan™ Array Plates**
 - **TaqMan™ Assays**
3. Enter the information.

Option	Information
Fixed TaqMan™ Array Plates	<i>Part number</i> <i>Batch number</i>
Custom TaqMan™ Array Plates	<i>Sales order</i> <i>Lot number</i>
TaqMan™ Assays	<i>Sales order</i> <i>Rack ID</i>

4. Click **Import**.

The plate layout contains the targets and samples.

■ Screens available during a run	70
■ Edit a plate file during a run	71
■ Pause an instrument run	72
■ Stop an instrument run	72
■ View the method during a run	73
■ View real-time data and plots during a run	73
■ View the time remaining during a run	74
■ Lock the touchscreen during a run	74
■ Unlock the touchscreen during a run	75
■ Instrument Schedule	75

Screens available during a run

From the home screen, you can swipe or tap the arrows to navigate to different screens. The following screens are available:

- A screen displaying the method (swipe left once or tap the right arrow (>) once)
- The real-time amplification plot (swipe left twice or tap the arrow right (>) twice)
- The time remaining in the run (swipe left three times or tap the right arrow (>) three times)

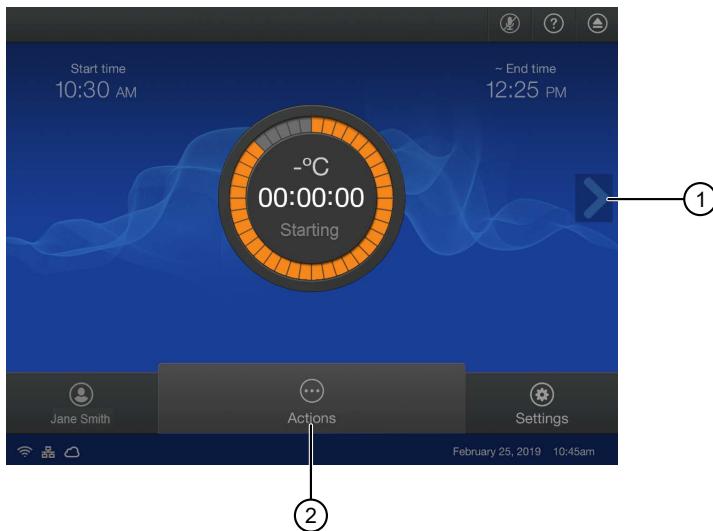


Figure 10 Home screen during a run

- ① Arrow
- ② Actions menu

Edit a plate file during a run

The targets and samples can be added or edited after a run has started. The method cannot be edited after a run has started.

1. In the home screen during a run, tap  **(Actions)**.

See Figure 10 on page 71.

2. Tap **Set up**.

3. In the **Plate** tab, tap a well or multiple wells.

Note: Use the plate view to edit or assign a well (see “Options to view a plate layout” on page 63).

4. Tap the **Sample Information** pane, then tap **Samples**.

Note: If no wells are selected, the pane is labeled **Legend**.

5. Enter a sample name, then select a sample type from the dropdown list.

- Unknown
- Standard
- No template control (NTC)

6. Tap **Target/SNP**, then select **Target** or **Assay** from the dropdown list.

7. Tap a row in the table, then tap **Close ▶ Done**.

Pause an instrument run

Pause a run to add reagents to the plate or to edit the number of cycles.

1. In the home screen during a run, tap  **(Actions)**.
See Figure 10 on page 71.
2. Tap **Pause run**, then enter a pause temperature.
3. *(Optional)* Tap **Edit** to change the number of cycles.
4. *(Optional)* Tap  **(Eject)** to open the drawer, then tap  **(Eject)** to close the drawer.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the plate temperature can reach 100°C. If you want to access the plate during a run pause, enter room temperature as the pause temperature and allow the plate to cool to room temperature before handling.

Stop an instrument run




CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the plate temperature can reach 100°C. If you want to access the plate during a run pause, enter room temperature as the pause temperature and allow the plate to cool to room temperature before handling.

1. In the home screen during a run, tap  **(Actions)**.
See Figure 10 on page 71.
2. Tap **Stop run**.
3. Unload the plate (see “Unload a plate or a TaqMan™ Array Card from the instrument” on page 52).

View the method during a run

1. In the home screen during a run, swipe left once or tap the right arrow (>) once.

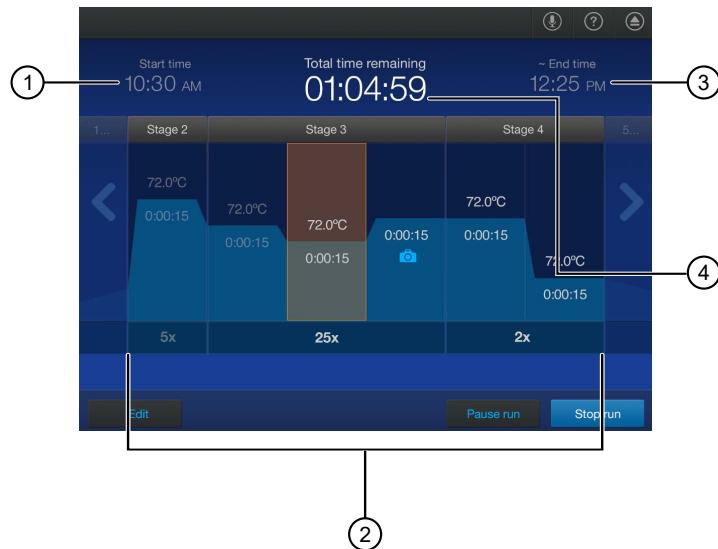

- ① Start time
- ② Method, including the current stage
- ③ End time
- ④ Time remaining

2. Swipe right once or tap the left arrow (<) once to return to the home screen.

View real-time data and plots during a run

1. In the home screen during a run, swipe left twice or tap the right arrow (>) twice.



2. *(Optional)* Tap the dropdown list then select the view for the plot.
 - **Linear**
 - **Log 10**
3. *(Optional)* Tap **Actions** ▶ **Analyze amp plots**.
4. *(Optional)* Tap **Actions** ▶ **View well details** to see the amplification for a specific well.
5. Swipe right twice or tap the left arrow (⟨) twice to return to the home screen.

Adjust the display of real-time plots during a run

1. In the real-time amplification plot, tap  **(Zoom)** or pinch.
2. Tap the arrows to pan left or right, or swipe left or right within the amplification plot.

View the time remaining during a run

1. In the home screen during a run, swipe left three times or tap the right arrow (⟩) three times. The time remaining in the run is displayed.
2. Swipe right three times or tap the left arrow (⟨) three times to return to the home screen.

Lock the touchscreen during a run

After you have started a run, you can lock the touchscreen so that other users cannot interfere with instrument operation.

Note: You must be signed in.

1. In the home screen during a run, tap  **(Profile)**.
2. Tap **Lock screen** ▶ **Lock**.
The status dial displays **Locked**.

The touchscreen automatically unlocks when the run is complete.

The touchscreen can be unlocked during the run (see “Unlock the touchscreen during a run” on page 75).

Unlock the touchscreen during a run

The touchscreen can be unlocked by the user who locked it or an administrator. The touchscreen automatically unlocks when the run is complete.

1. Tap anywhere on the touchscreen.
2. Tap the **PIN Code** field, then enter your PIN.
The status dial is displayed without **Locked**.

Instrument Schedule

Overview of the instrument schedule

The instrument schedule allows a user to reserve the instrument at a particular date and time.

A Connect profile is required.

The instrument is not locked to prevent any other users at the time.

Schedule a run

1. In the home screen, tap  **(Calendar)**.
2. Select a view.
 - **Day**
 - **Week**
 - **Month**
3. Tap a time.
The day on the calendar can be selected in the month view.
The day view and the week view show the schedule in 30-minute increments.
4. Edit the start date, end date, start time, and end time.
5. *(Optional)* Add notes.
6. Tap **Save**, then tap **Done** to exit the schedule.

Cancel a scheduled run

1. In the home screen, tap  **(Calendar)**.
2. Select a view.
 - **Day**
 - **Week**
 - **Month**

3. Tap the reservation.

The **Edit Reservation** screen is displayed.

4. Tap **Cancel Reservation**, then tap **Yes**.

6

View and manage files

■ Manage plate files	77
■ Manage data files	78

Manage plate files

Copy plate files

1. In the home screen, tap  **Load plate file** or **Set up run**.
2. Tap the appropriate folder.
 -  **Run Queue**
 -  **Public**
 -  **Connect**
 -  **Network Drive**
 -  **USB Drive**
3. Tap **Manage**.
4. Tap the file, then tap **Copy files**.
Multiple files can be selected.
5. Select the destination folder.
 -  **Public**
 -  **Connect**
 -  **Network Drive**
 -  **USB Drive**
6. Tap **Paste**.

Tap  **(Back)** to return to the home screen.

Delete plate files

System templates cannot be deleted. A plate file that was saved from a system template can be deleted.

For descriptions of files, see “Files” on page 28.

1. In the home screen, tap  **Load plate file** or **Set up run**.
2. Tap the appropriate folder.
 -  **Run Queue**
 -  **Public**
 -  **Connect**
 -  **Network Drive**
 -  **USB Drive**
3. Tap **Manage**.
4. Tap the file, then tap **Delete files**.
Multiple files can be selected.
5. Tap **Yes** to confirm.

Tap  **(Back)** to return to the home screen.

Manage data files

View QC Check status after a run

The status dial displays **Run complete** when a run is complete. After **Done** is tapped, the amplification plot can no longer be viewed on the instrument.

1. In the home screen after a run is complete, see the status of the QC Check.
The status of the QC Check is displayed under the status dial.

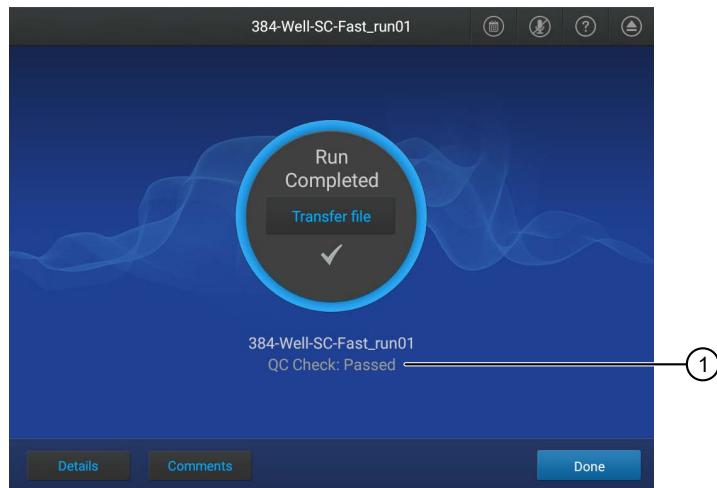


Figure 11 QC Check passed

① QC Check

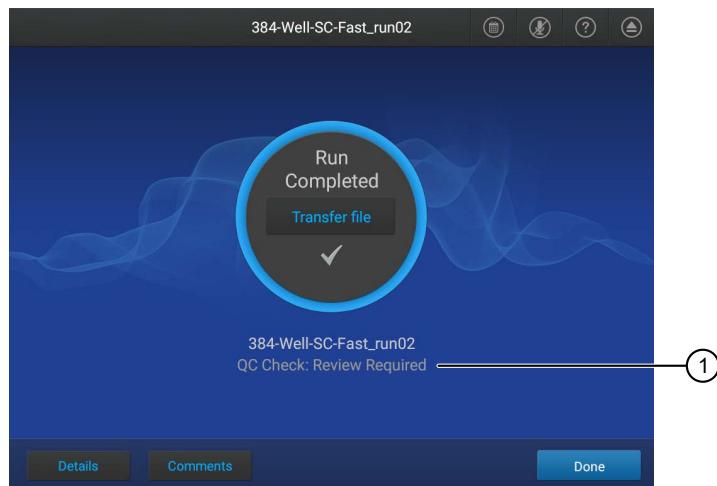



Figure 12 Review required for QC Check

① QC Check

2. See the QC data in the QuantStudio™ Design and Analysis Software v2 or the QuantStudio™ Design and Analysis v2.

View an amplification plot after a run

The status dial displays **Run complete** when a run is complete. The amplification plot can only be viewed immediately after a run. After **Done** is tapped, the amplification plot can no longer be viewed on the instrument.

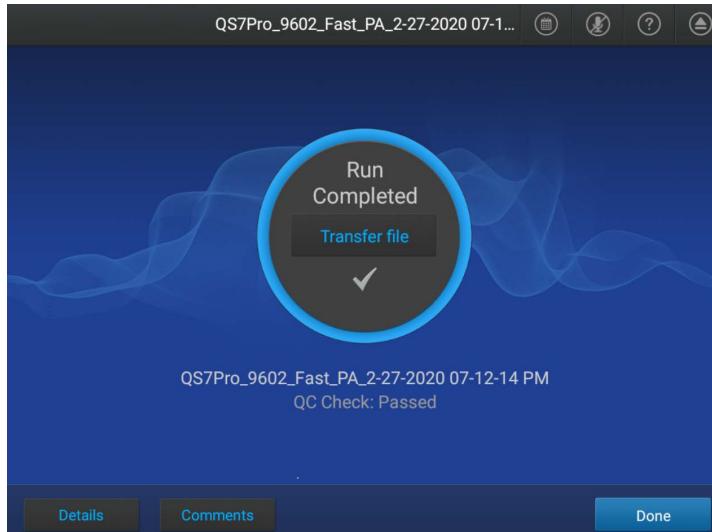

If the automatic sign-out feature is enabled and the set time has elapsed, the instrument will display the **Sign In** screen. The amplification plot cannot be viewed on the instrument after the run.

1. In the home screen after a run is complete, tap **Details**.
The amplification plot is displayed. The display can be adjusted (see “Adjust the display of real-time plots during a run” on page 74).
2. *(Optional)* On the **Details** screen, tap the **Plot Types** dropdown list, then select a plot type.
 - **Rn vs. Cycle**
 - **dRn vs. Cycle**
3. *(Optional)* On the **Details** screen, tap the **Y-Scales** dropdown list, then select the view for the plot.
 - **Linear**
 - **Log 10**
4. *(Optional)* Tap **Well details**, then tap **Samples** or **Targets**.
5. Select an amplification plot preview, then tap **View amp plot**.
6. *(Optional)* Tap a specific well on the right side of the amplification plot to see the details for the single well.
7. Tap **Close** to return to the home screen.

Transfer a data file immediately after a run

This section describes the transfer of a data file immediately after a run. To transfer a data file or a set of data files at a later time, see “Transfer data files at a later time” on page 81.

In the home screen, the status dial displays **Run complete** when a run is complete. The home screen also shows the location that the data file was transferred to. The location for transfer was selected before the run was started (see “Create and run a plate from a system template or a saved plate file” on page 47 and “Repeat the last run” on page 49).

If the automatic sign-out feature is enabled and the set time has elapsed, the instrument will display the **Sign In** screen. Data files cannot be transferred from the home screen. They can be transferred at a later time (see “Transfer data files at a later time” on page 81).

Data files will be transferred to the location that was selected before the run was started, even if the user was automatically signed out.

1. In the home screen after a run is complete, tap **Transfer file**.
2. Select the destination for the data file.

-  **Public**
-  **Connect**

Note: You must be signed in with a Connect profile.

-  **Network Drive**
-  **USB Drive**
-  **My Instrument**

Note: This option is not available if you are using the instrument as a guest. You must be signed in.

3. Tap **OK**, then tap **Transfer**.

Transfer data files at a later time

This section describes the transfer of a data file or a set of data files at any time. To transfer a data file immediately after a run, see “Transfer a data file immediately after a run” on page 81.

- Guests can only transfer guest data files.
- A user that is signed in can transfer their own data files.
- Administrators can transfer all data files.

1. Tap  **(Settings)** ▶ **Run history**.

The **Run History** screen is displayed. It contains a list of the runs performed on the instrument associated with the profile. Administrators can view all of the runs performed.

2. *(Optional)* Tap the table header to sort by that parameter.

- **Run name**
- **Date**
- **Transferred**
- **User** (Administrator only)

3. Tap the data files to be transferred or tap **Select all**.

Multiple data files can be selected.

4. Tap **Transfer**, then tap the storage location.

5. Tap **Done**.

View the list of data files on the instrument

- Guests can only view the list of guest data files.
- A user that is signed in can view the list of their own data files.
- Administrators can view all data files.

1. In the home screen, tap  **(Settings)** ▶ **Run history**.

The **Run History** screen is displayed.

2. Tap an individual run record, then tap **View**.

The following details are displayed:

• The file name	• The duration of the run
• The run status	• The reaction volume
• The user name	• The cover temperature
• The start time and the end time	

3. *(Optional)* Tap **Transfer** to transfer the data file.

4. *(Optional)* Tap **Delete** to delete the data file from the instrument.

5. Tap **Cancel** to return to the **Run History** screen.

Delete data files from the instrument

- Guests cannot delete any data files.
- A user that is signed in can delete their own data files.
- Administrators can delete all the data files.

1. In the home screen, tap  **(Settings)** ▶ **Run history**.

The **Run History** screen is displayed. It contains a list of the runs performed on the instrument associated with the profile. Administrators can view all of the runs performed.

2. *(Optional)* Tap the table header to sort by that parameter.

- **Run name**
- **Date**
- **Transferred**
- **User** (Administrator only)

3. Tap the data files to be deleted or tap **Select all**.

Multiple data files can be selected.

4. Tap **Delete**, then tap **Yes** to confirm.

Calibrate and verify instrument performance

■ Instrument and block calibration	84
■ Calibration and verification schedule	85
■ Calibration descriptions	86
■ View the calibration status and set reminders	87
■ Export a calibration report	89
■ Perform ROI/uniformity, background, and dye calibrations for plate blocks	89
■ Perform ROI/uniformity, background, dye, and normalization calibrations for the TaqMan™ Array Card block	97
■ Perform verification using RNase P plates	103
■ Perform verification using RNase P with a TaqMan™ Array Card	109
■ Calibrate custom dyes	113
■ Calibrate for a custom melt curve run	120

Instrument and block calibration

The QuantStudio™ 6 Pro System and the QuantStudio™ 7 Pro System have interchangeable blocks. The calibration status applies to the combination of the instrument and the block.

If a block is moved to an instrument that it has not been calibrated on, that combination of block and instrument needs to be calibrated. This applies even if the block has previously been calibrated on a different instrument.

If a block is moved back to an instrument that has previously been calibrated on, it does not need to be calibrated again. It needs to be calibrated again if the calibration is expired.

A run can still be started if a calibration of the block on the instrument is expired. A run cannot be started if a block has never been calibrated on the instrument.

Calibration and verification schedule

The instrument and block will be calibrated during the installation. To ensure optimal performance, perform calibrations at the recommended frequency.

Note: After instrument calibration, we recommend performing instrument verification using the provided RNase P plate.

IMPORTANT! Perform calibrations and instrument runs under the specified environmental conditions (see “Environmental requirements” on page 193). Exposure to extreme temperatures can adversely affect the instrument performance and shorten the life span of the instrument components.

To set calibration frequencies and reminders, see one of the following sections:

- View the calibration status on the instrument (see page 87)
- View calibration status and set reminders on the Connect platform (see page 88)

Calibration	Recommended frequency
ROI/Uniformity	<ul style="list-style-type: none">• Every two years (recommended).• Always perform new Background and Dye calibrations after an ROI/Uniformity calibration. You will be prompted to run a Background calibration after the ROI/Uniformity calibration is complete. The ROI/Uniformity calibration and the Background calibration must pass in order to accept the results. It is recommended to transfer the files if the calibration fails in order to troubleshoot. <p>Note: Performing an ROI/Uniformity calibration invalidates all other calibrations.</p>
Background	<ul style="list-style-type: none">• Every month (recommended).• Background calibration can also be performed, as needed:<ul style="list-style-type: none">– To check for contamination (depends on usage and laboratory conditions).– To obtain the most accurate data for the removal of background fluorescence. <p>Note: Performing a Background calibration does <i>not</i> invalidate any other calibration.</p>
Dye	<ul style="list-style-type: none">• Every two years (recommended).• During a Dye calibration, only the dyes on the given spectral calibration plate are calibrated. <p>Note: Performing a Dye calibration for a given dye plate does <i>not</i> invalidate any other calibration.</p>
Normalization (<i>TaqMan™ Array Card block only</i>)	Every two years (recommended).

(continued)

Calibration	Recommended frequency
RNase P instrument verification	<ul style="list-style-type: none"> After installing or moving the instrument. (Recommended) After performing any of the calibrations. As needed to confirm instrument performance.

Note: To prepare custom dye plates and to perform custom calibrations, see “Calibrate custom dyes” on page 113.

Calibration descriptions

The combination of the instrument and the block is calibrated (see “Instrument and block calibration” on page 84).

Calibration description and purpose	Pass criteria
ROI/Uniformity	
<ul style="list-style-type: none"> The software captures images for each optical filter. The software uses calibration data to map the increase in fluorescence to the plate wells during subsequent runs and to evaluate well-to-well consistency of the signals. 	The image for each filter distinguishes all wells of the plate. Each well in the image is distinct.
Background	
<ul style="list-style-type: none"> The software captures background images for each optical filter in the absence of sample and reagent, and it checks that the fluorescence from each well is below a fluorescence threshold. The software uses calibration data to remove background fluorescence during a run. <p>Note: You can also run this calibration to determine if contamination is related to the sample block or the plate.</p>	The plate images for all filters are free of abnormal fluorescence.
Dye	
<ul style="list-style-type: none"> The software extracts a spectral profile for each dye standard, then produces a set of spectral profiles plotted as fluorescence vs filter. The software uses calibration data to characterize and distinguish the individual contribution of each dye in the total fluorescence signals collected by the instrument. 	Dye spectra peak within the same filter as their group.

(continued)

Calibration description and purpose	Pass criteria
Normalization (TaqMan™ Array Card block only)	
<ul style="list-style-type: none"> The software extracts a spectral profile for each normalization standard. The software uses the normalization standards to obtain a correction factor for the specific dyes, for example, FAM™ dye or VIC™ dye. The correction factor is used to generate comparable component signal levels across instruments. 	Normalization standards have sufficient signals to generate a correction factor.

View the calibration status and set reminders

View the calibration status on the instrument

The **Status and Reminders** screen also allows you to perform the following functions:

- Transfer the calibration results (see page 94)
- Transfer the calibration data files (see page 94)
- View the calibration images (see page 93)

- In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Status and reminders**.

The **Status and Reminders** screen is displayed. It shows the following information:

- Type of calibration

Note:

- For dye calibration, the type of calibration is listed as the dye that was calibrated, for example, **FAM**.
- Normalization calibration is required for a TaqMan™ Array Card block. The type of calibration is listed as **FAM/ROX** or **VIC/ROX**.

- Status of the calibration
- Last calibration
- Date that the calibration expires
- If a reminder is set

- Tap a calibration row, then tap **View** to view the history of that specific calibration type.
 - Date performed
 - Result of the calibration
 - Whether the calibration is current
 - Operator
- Tap **Cancel** to return to the **Status and Reminders** screen.

Set calibration frequency and reminders on the instrument

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Status and reminders**.
The **Status and Reminders** screen is displayed.
2. Tap a calibration row, then tap **Edit**.
3. Tap on the **Expiration Interval** field, then enter the number of days.
The default is 730 days (2 years).
4. Tap on the **Reminder** field, then enter the number of days
The **Next Reminder** field will be populated.
5. Set the **On/Off** slider to the **On** position to enable a reminder about calibration.
6. Tap **Save**.

The calibration reminder can be set as a home screen notification (see “Enable and disable home screen notifications” on page 187).

View calibration status and set reminders on the Connect platform

Note: The calibration reminders feature requires a connection between the instrument and a computer network.

1. Sign in to Connect on a browser.
2. Tap  **(Instrument)** to open InstrumentConnect.
3. Select any of your registered instruments.
4. In the **Summary** tab:
 - Tap  **Calibrations** to view the status of each calibration type.
 - *(Optional)* Tap  **Calibration Reminders** to set the calibration reminder time table and enter the notification email address or addresses.

Note: More than one email address can receive the calibration reminders.

Note: The settings are automatically saved.

- *(Optional)* In the **Downloads** section, tap **Maintenance Summary.pdf** to download the calibration status report.

5. In the **Calibrations History** tab:
 - View the history of each calibration type.
 - *(Optional)* Tap  to download the calibration history report.

Review calibration status

The calibration status can be reviewed in the QuantStudio™ Design and Analysis Software v2. The instrument must be added to the **Manage Instruments** list.

1. In the home screen, click **Manage Instruments**.

To access instruments from a different screen, click  **System** ▶ **Instruments**.

2. In the **Summary** tab for the instrument, the calibration status is displayed in the **Calibration Status** column.

Export a calibration report

The calibration report is exported as a TXT file.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Status and reminders**.
2. Tap **Export**, then select a destination.
 -  **Connect**
 -  **Network Drive**
 -  **USB Drive**

The connectivity status of each location is displayed.

3. Tap **Done**.

Perform ROI/uniformity, background, and dye calibrations for plate blocks

To calibrate the TaqMan™ Array Card block, see “Perform ROI/uniformity, background, dye, and normalization calibrations for the TaqMan™ Array Card block” on page 97.

Calibration workflow

Perform an ROI/uniformity calibration

You are automatically prompted to perform background calibration.



Perform a background calibration

Perform any time that ROI/uniformity calibrations are current.



Perform system dye calibrations

Perform any time that ROI/uniformity and background calibrations are current.



(Optional) Perform custom dye calibrations (see “Calibrate custom dyes” on page 113)

Perform any time that ROI/uniformity and background calibrations are current.

Prepare a calibration plate

Required materials to prepare calibration plates

Note: We recommend calibrating with all Spectral Dye Calibrations Plates available for your block configuration even if you are not using all the dyes in the plates. This is recommended for optimal system performance.

Unless otherwise indicated, all materials are available through thermofisher.com. "MLS" indicates that the material is available from fisherscientific.com or another major laboratory supplier.

Item	Source
Calibration plates	
ROI/Uniformity plate	
Background calibration plate	
Dye calibration plates	
Equipment and consumables	
Centrifuge with plate adapter; buckets cleaned before use	MLS
Powder-free gloves	MLS
Safety glasses	MLS

Thaw, vortex, and centrifuge a calibration plate

IMPORTANT! Keep calibration plates protected from light until you perform the calibration. Do not remove the plate from its packaging until you are ready to use it. Prolonged exposure to light can diminish the fluorescence of the dyes in the wells of calibration plates.

1. Remove the calibration plate from the freezer, then thaw the plate in its packaging for 30 minutes.

IMPORTANT! Use each plate within 2 hours of thawing.


2. While wearing powder-free gloves, remove the calibration plate from its packaging. Do not remove the optical film.

Note: Do not discard the packaging for the calibration plate. Each calibration plate can be used up to three time if the following conditions are met:

- The plate is stored in its packing sleeve at -25°C to -15°C .
- The plate is used within 6 months after opening.
- The plate is used before the plate expiry date.

3. Vortex the plate for 5 seconds, then centrifuge at 750–1,000 $\times g$ for 2 minutes.

4. Confirm that the liquid in each well is at the bottom of the well and free of bubbles. If it is not, centrifuge the plate again.

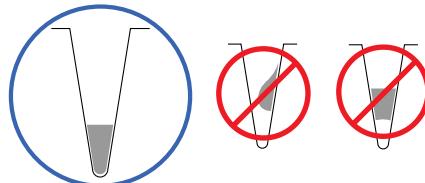

IMPORTANT! Keep the bottom of the plate clean. Fluids and other contaminants on the bottom of the plate can contaminate the sample block and cause an abnormally high background signal.

Perform calibrations

1. In the home screen tap **(Settings)** **Maintenance and service** **Calibration**.
The **Calibration** screen is displayed.

2. In the **Calibration** screen, select a calibration type.

Calibration type	Tap
ROI/uniformity	ROI and uniformity
Background	Background
Dye	System dye

Note: You will be prompted to perform a background calibration after a ROI/Uniformity calibration.

3. *(System dye calibration only)* Select the calibration plate.
4. Follow the instructions on the screen to start the calibration.
5. Load the plate into the instrument.



CAUTION! The instrument should be used by trained operators who have been warned of the moving parts hazard.

6. Tap **Start**.

When the ROI/Uniformity run is complete, you will be prompted to perform a background calibration.

When the Background calibration is complete, the **Calibration Status** will display. It will indicate whether the calibration passed or failed.

7. *(For calibrations that pass)* Select an option.

Option	Description
Details	View the calibration images.
Transfer EDS	Transfer the calibration file to Connect, a USB drive, or a network drive.
Reject results	The calibration file is deleted.
Accept results	The calibration file is saved.

If the results are accepted and the calibration file is saved, previous calibration files associated with the combination of the instrument and the block will no longer be used.

8. *(For calibrations that fail)* Select an option.

Option	Description
Details	View the calibration images.
Transfer EDS	Transfer the calibration file to Connect, a USB drive, or a network drive (recommended for troubleshooting).
Close	The calibration file is deleted.

See “Troubleshoot calibration failure” on page 95.

9. Unload the plate from the instrument.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the plate temperature can reach 100°C. Allow it to cool to room temperature before handling.

Note: Use Smart Help if the instrument does not eject the plate (see “Request technical support with Smart Help” on page 172).

10. Return the plate to its original packaging.

Note: Do not discard the packaging for the calibration plate. Each calibration plate can be used up to three time if the following conditions are met:

- The plate is stored in its packing sleeve at -25°C to -15°C .
- The plate is used within 6 months after opening.
- The plate is used before the plate expiration date.

View the calibration images

The instrument performs the ROI, uniformity, and background calibrations in sequence. You can view the calibration images after the background calibration is complete.

1. In the home screen, tap  (Settings) ▶ Maintenance and service ▶ Calibration ▶ Status and reminders.

The **Status and Reminders** screen is displayed.



2. Tap a calibration type in the table, then tap **View**.

3. Tap a calibration date in the table, then tap **View**.



The following details are displayed:

• Calibration name	• The operator
• Result	• Reagent name, part number, lot number, and expiration date
• Date	• Comments.
• Whether this calibration is current	

4. Tap **Detail**.

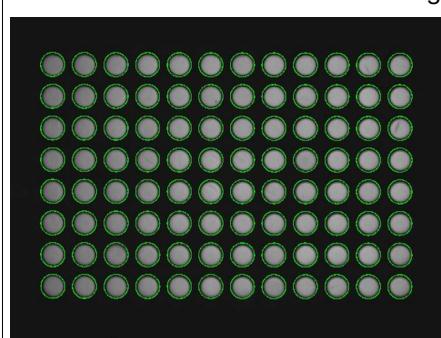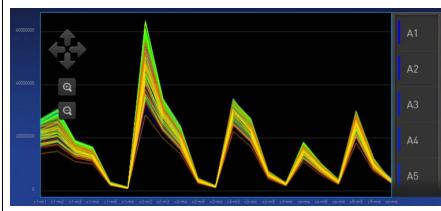Calibration	Example results indicating successful calibration
ROI Note: Select the desired filter combination from the Filter Set dropdown list.	Green circles around all wells and bright well centers. 
Uniformity	Signals from each well following a uniform trend. 

(continued)

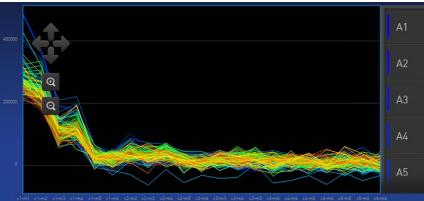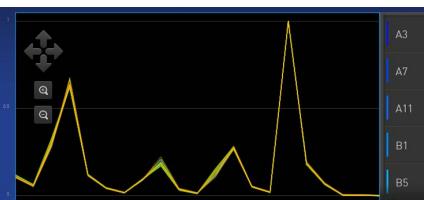Calibration	Example results indicating successful calibration
Background	Few, if any, signals with abnormally high fluorescence. 
Dye	Signals from each well following a uniform trend, and each dye peaks at the correct filter. 

5. (Optional) For Uniformity, Background, and Dye calibration images, tap on an individual well on the right side to see the results for a single well.

Transfer calibration results

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Status and reminders**.
The **Status and Reminders** screen is displayed.
2. Tap a calibration type in the table.
3. Tap **Export**.

Transfer calibration data files

This section describes transferring calibration files if the calibration passed.

If the calibration fails, the calibration data file is not saved. The data file can only be transferred immediately after the calibration (see “Perform calibrations” on page 91).

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Status and reminders**.
The **Status and Reminders** screen is displayed.
2. Tap a calibration type in the table, then tap **View**.
3. Tap a calibration date in the table, then tap **View**.
4. Tap **Transfer EDS**.

Troubleshoot calibration failure

Observation	Possible cause	Recommended action
Calibration failed	The plate was improperly prepared.	<p>Ensure the following:</p> <ul style="list-style-type: none"> • The correct plate was used for the calibration performed. • The plate was properly thawed. • The plate was properly centrifuged. • The plate was properly sealed.
	The plate is damaged or contaminated.	<p>Check for damage, improper plate seal, or contamination.</p> <p>Order a replacement plate. Use Smart Help if the replacement plate fails (see “Request technical support with Smart Help” on page 172).</p>
High fluorescence signal in individual wells	Signals that exceed the limit of normal fluorescence may indicate fluorescent contaminants on the plate or the sample block.	See “Identify contamination” on page 95.
Calibration failed but plate is undamaged	The incorrect plate was used for calibration performed.	Use the plate that matches the calibration performed.
	The plate was improperly prepared.	<p>Repeat the calibration with the plate properly prepared.</p> <p>If the calibration fails again, order a replacement plate. Use Smart Help if the replacement plate fails (see “Request technical support with Smart Help” on page 172).</p>

Identify contamination

Signals that exceed the limit of normal fluorescence may indicate fluorescent contaminants on the calibration plate or the sample block. Common contaminants include ink residue from permanent pens, powder from disposable gloves, and dust.

1. View the calibration data and note the wells that failed the calibration.
2. Remove the plate from the instrument, rotate the plate 180°, then perform the calibration again.

3. Determine the location of the failed wells again as in step 1.

Position of failed wells	Action
Identical	The sample block is contaminated. Decontaminate the sample block (see page 159).
Reversed	The plate is contaminated. Discard the plate, then perform the calibration using a new calibration plate.

4. Use Smart Help if the calibration fails after you decontaminate the sample block and replace the plate (see “Request technical support with Smart Help” on page 172).

Create a background plate (optional)

Whenever possible, use a background plate listed in “Required materials to prepare calibration plates” on page 90. These plates contain a buffer that accurately simulates the reagents used for PCR, and, therefore, produces high-quality calibration data.

If a background plate is not available, you can create one as described below.

Required materials:

- MicroAmp™ optical 96-well reaction plate
- Optical adhesive cover or optical flat caps
- Pipettor, 200-µL (with pipette tips)
- Powder-free gloves
- Safety glasses
- Deionized water

IMPORTANT! Wear powder-free gloves while creating the background plate.

1. Remove a reaction plate from its box and place it on a clean, dry surface.
2. Aliquot 50 µL of deionized water to each well of the reaction plate.
3. Seal the plate using an optical adhesive cover or optical flat caps.
4. Use the plate for background calibration.

Perform ROI/uniformity, background, dye, and normalization calibrations for the TaqMan™ Array Card block

To calibrate the plate blocks, see “Perform ROI/uniformity, background, and dye calibrations for plate blocks” on page 89.

This section covers procedures specific to the TaqMan™ Array Card block. See the following sections for additional information:

- Transfer calibration results (page 94)
- Transfer data files from calibrations (page 94)
- Troubleshoot calibration failure (page 95)

To troubleshoot normalization calibration, see “Troubleshoot normalization calibration failure” on page 102.

- Identify contamination (page 95)

Calibration workflow

Perform an ROI/uniformity calibration

You are automatically prompted to perform background calibration.



Perform a background calibration

Perform any time that ROI/uniformity calibrations are current.



Perform system dye calibrations

Perform any time that ROI/uniformity and background calibrations are current.



Perform normalization calibration

Perform after a dye calibration.



(Optional) Perform custom dye calibrations (see “Calibrate custom dyes” on page 113)

Perform any time that ROI/uniformity and background calibrations are current.

Prepare a calibration TaqMan™ Array Card

Required materials to prepare calibration TaqMan™ Array Card

Unless otherwise indicated, all materials are available through thermofisher.com. "MLS" indicates that the material is available from fisherscientific.com or another major laboratory supplier.

Item	Source
Calibration kit	
ViiA™ 7 Array Card Spectral Calibration Kit, includes the following components: <ul style="list-style-type: none"> Empty TaqMan™ Array Cards FAM™ dye VIC™ dye ROX™ dye ROI dye mix Background buffer FAM™/ROX™ dye VIC™/ROX™ dye 	4432314
Equipment and consumables	
Centrifuge with custom buckets and card holders, one of the following: <ul style="list-style-type: none"> Sorvall™ centrifuge Multifuge™ centrifuge Megafuge™ centrifuge 	Contact your local sales office
TaqMan™ Array Card Sealer	Contact your local sales office
Blank balance TaqMan™ Array Cards	Contact your local sales office
Pipette, 200 µL with tips	MLS
Permanent marker or pen	MLS
Powder-free gloves	MLS
Safety glasses	MLS

Fill the calibration TaqMan™ Array Cards


For detailed instructions to prepare a TaqMan™ Array Card, see Appendix C, "Detailed procedures for preparation of a TaqMan™ Array Card".

IMPORTANT! Wear powder-free gloves while creating the calibration array cards.

This procedure explains how to create *all* of the array cards required to calibrate the block, but not all of them are required for at each calibration. Before preparing array cards for calibration, determine which calibrations are required (see “Calibration and verification schedule” on page 85).

Note: For more information on the array card loading procedure, contact Technical Support.

1. Remove the tubes of calibration solutions from the freezer, allow them to thaw, then vortex the tubes to mix the contents well.
2. Remove the TaqMan™ Array Cards from their box and place them on a clean, dry surface.
3. Mark the side of the empty array cards with the following items:
 - Background
 - FAM™
 - ROI
 - ROX™
 - VIC™
 - FAM™/ROX™
 - VIC™/ROX™

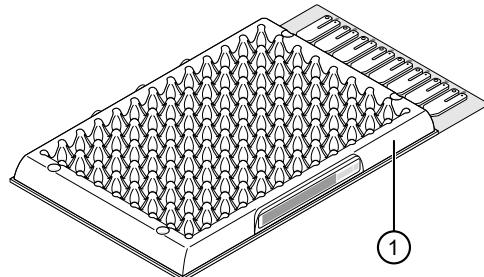

① Location to mark each TaqMan™ Array Card

4. For each array card, pipet 100 µL of the appropriate calibration solution into each of the eight reservoirs in the array card.
See “Load the PCR reaction mix” on page 205.
5. Centrifuge the array cards, then check each array card for the correct fill.
See “Centrifuge the card” on page 205.
6. Seal the array cards.
See “Seal the card” on page 207.

IMPORTANT! If an array card is sealed improperly, it can leak and contaminate the sample block and/or it can cause the associated calibration or RNase P test to fail.

IMPORTANT! After the array cards are sealed, store them in a dark place. Do not expose the array cards to light until you are ready to use them. The dyes in the array cards are photosensitive. Prolonged exposure to light can diminish the fluorescence of the dye.

Perform calibrations

1. In the home screen tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration**.
 The **Calibration** screen is displayed.
2. In the **Calibration** screen, select a calibration type.

Calibration type	Tap
ROI/uniformity	ROI and uniformity
Background	Background
Dye	System dye
Normalization	Instrument normalization

Note: You will be prompted to perform a background calibration after a ROI/Uniformity calibration.

3. *(System dye calibration only)* Select the dye to calibrate.
 - **FAM**
 - **VIC**
 - **ROX**
4. *(Normalization only)* Select the reporter dye and passive reference to calibrate.
 - **FAM/ROX**
 - **VIC/ROX**
5. Follow the instructions on the screen to start the calibration.
6. Load the TaqMan™ Array Card into the instrument.
 See “Load a TaqMan™ Array Card into the instrument” on page 52.



CAUTION! The instrument should be used by trained operators who have been warned of the moving parts hazard.

7. Tap **Start**.
 When the ROI/Uniformity run is complete, you will be prompted to perform a background calibration.
 When the Background calibration is complete, the **Calibration Status** will display. It will indicate whether the calibration passed or failed.

8. (For calibrations that pass) Select an option.

Option	Description
Details	View the calibration images.
Transfer EDS	Transfer the calibration file to Connect, a USB drive, or a network drive.
Reject results	The calibration file is deleted.
Accept results	The calibration file is saved.

If the results are accepted and the calibration file is saved, previous calibration files associated with the combination of the instrument and the block will no longer be used.

9. (For calibrations that fail) Select an option.

Option	Description
Details	View the calibration images.
Transfer EDS	Transfer the calibration file to Connect, a USB drive, or a network drive (recommended for troubleshooting).
Close	The calibration file is deleted.

See “Troubleshoot calibration failure” on page 95.

10. Unload the TaqMan™ Array Card from the instrument.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the TaqMan™ Array Card temperature can reach 100°C. Allow it to cool to room temperature before handling.

Note: Use Smart Help if the instrument does not eject the array card (see “Request technical support with Smart Help” on page 172).

View the calibration images

The instrument performs the ROI, uniformity, and background calibrations in sequence. You can view the calibration images after the background calibration is complete.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Status and reminders**.
The **Status and Reminders** screen is displayed.
2. Tap a calibration type in the table, then tap **View**.
The following details are displayed:
3. Tap a calibration date in the table, then tap **View**.

- Calibration name
- Result
- Date
- Whether this calibration is current
- The operator
- Reagent name, part number, lot number, and expiration date
- Comments.

4. Tap **Detail**.
5. *(Optional)* For Uniformity, Background, and Dye calibration images, tap on an individual well on the right side to see the results for a single well.

Troubleshoot normalization calibration failure

Normalization calibration is only required for the TaqMan™ Array Card. For all other calibration troubleshooting, see “Troubleshoot calibration failure” on page 95.

Observation	Possible cause	Recommended action
Normalization calibration failed	Abnormally low spectra across all wells of the array card.	 CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the TaqMan™ Array Card temperature can reach 100°C. Allow it to cool to room temperature before handling. <ol style="list-style-type: none"> 1. Confirm that you loaded the correct array card into the instrument. If not, perform the calibration again using the correct consumable. 2. If you are using the correct array card, perform the calibration again using a different normalization consumable. 3. If the calibration fails again, contact Support.

Observation	Possible cause	Recommended action
Normalization calibration failed (continued)	One or more wells produced spectra that deviate significantly from the rest of the array card.	<ol style="list-style-type: none">1. Locate the wells with the abnormal signal (see “View the calibration images” on page 101).2. Rotate the array card 180°, then perform the calibration again.3. Determine the location of the contaminated wells again. If the positions of the wells identified in the steps above are:<ul style="list-style-type: none">• Identical: The sample block is contaminated. Decontaminate the sample block.• Reversed: The normalization array card is contaminated. Discard the consumable, then perform the normalization calibration using a new normalization array card.4. If the calibration fails again, contact Support.

Perform verification using RNase P plates

Verification should be performed under the following circumstances:

- After installation and before first use of the instrument and the block
- *(Recommended)* After performing instrument and block calibrations
- As needed to confirm instrument and block performance

The instrument and block require valid ROI/uniformity, background, and dye calibrations to perform verification.

Instrument verification description

Purpose	Description	Pass criteria
Confirms the performance of the instrument.	Quantifies the number of copies of the human RNase P gene in samples with known concentrations of the corresponding DNA template.	<p>The instrument passes performance specifications if the following inequality is true and the instrument successfully distinguishes between unknown populations A and B with a statistical confidence level of 99.7%.</p> $[(C_{qA}) - 3(\sigma_{CqA})] > [(C_{qB}) + 3(\sigma_{CqB})]$ <p>where:</p> <ul style="list-style-type: none"> • C_{qA} = Average C_q of unknown population A • σ_{CqA} = Standard deviation of unknown population A • C_{qB} = Average C_q of unknown population B • σ_{CqB} = Standard deviation of unknown population B

RNase P instrument verification plate

The RNase P plate contains the reagents necessary for the detection and quantitation of genomic copies of the human RNase P gene (a single-copy gene encoding the RNase moiety of the RNase P enzyme). Each well contains the following components:

- PCR Master Mix
- RNase P primers
- A FAM™ dye-labeled probe
- A known concentration of human genomic DNA template

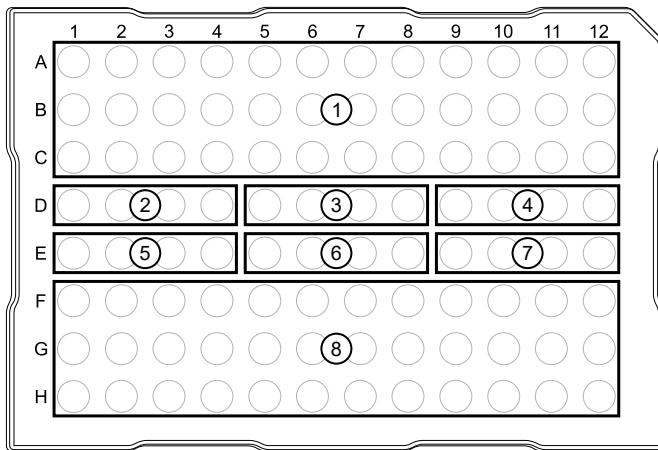


Figure 13 96-well RNase P plate

① Unknown A (5,000)	⑤ STD 5,000 copies
② NTC (no template control)	⑥ STD 10,000 copies
③ STD 1,250 copies	⑦ STD 20,000 copies
④ STD 2,500 copies	⑧ Unknown B (10,000)

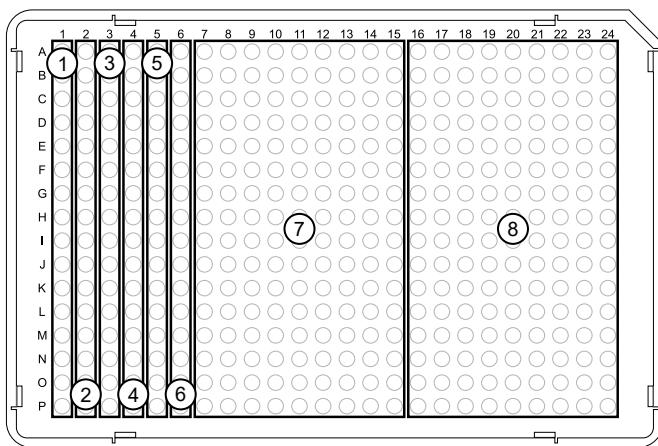


Figure 14 384-well RNase P plate

① NTC (no template control)	⑤ STD 10,000 copies
② STD 1,250 copies	⑥ STD 20,000 copies
③ STD 2,500 copies	⑦ Unknown A (5,000)
④ STD 5,000 copies	⑧ Unknown B (10,000)

Performance specifications pass criteria

After the run, the software calculates average copy number values and standard deviation values. The instrument passes performance specifications if the following inequality is true and the instrument successfully distinguishes between unknown populations A and B with a statistical confidence level of 99.7%.

$$[(C_{qA}) - 3(\sigma_{CqA})] > [(C_{qB}) + 3(\sigma_{CqB})]$$

where:

- C_{qA} = Average C_q of unknown population A
- σ_{CqA} = Standard deviation of unknown population A
- C_{qB} = Average C_q of unknown population B
- σ_{CqB} = Standard deviation of unknown population B

The software automatically adjusts the threshold and omits a defined number of wells from the unknown populations to meet the performance specifications. To view any omitted wells, open the data file for the verification in the desktop software.

Prepare an RNase P plate

Materials required for RNase P plate preparation

- RNase P instrument verification plate
- Centrifuge with plate adapter; buckets cleaned before use
- Powder-free gloves
- Safety glasses


Thaw, vortex, and centrifuge an RNase P plate

IMPORTANT! Expose the RNase P plate to room temperature for no more than 45 minutes, inclusive of thawing and preparation time.

After thawing, the RNase P plate cannot be refrozen.

1. Remove the RNase P plate from the freezer, then thaw the plate in its packaging.
 - Thaw the plate for approximately 5 minutes.
 - Use the plate within 30 minutes of thawing.
2. Confirm that the bench, vortex, and centrifuge are clean. Before use, wipe the vortex and centrifuge using a lint-free tissue.
3. While wearing powder-free gloves, remove the plate from its packaging.
4. Vortex the plate for 5 seconds, then centrifuge at 750–1,000 $\times g$ for 2 minutes.

5. Confirm that the liquid in each well is at the bottom of the well and free of bubbles. If it is not, centrifuge the plate again.

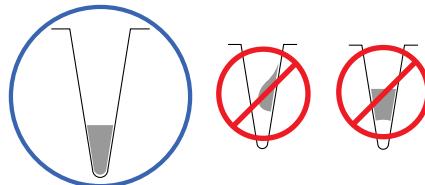

IMPORTANT! Keep the bottom of the plate clean. Fluids and other contaminants on the bottom of the plate can contaminate the sample block and cause an abnormally high background signal.

Perform RNase P verification

1. In the home screen, tap **(Settings)** ▶ **Maintenance and service** ▶ **RNase P verification**.
2. Follow the instructions on the screen to start the verification.
3. Load the plate into the instrument.



CAUTION! The instrument should be used by trained operators who have been warned of the moving parts hazard.

4. Tap **Start**.
5. When the run is complete and the screen displays **Complete**, tap **View results** to confirm the status of the run.

Verification status	Action
Passed	Instrument is ready for use.
Failed	See “Troubleshoot verification failure” on page 108.

6. In the **RNase P Verification Status** screen, tap one of the following options.

Option	Description
Accept results	Save the results to the instrument.
Reject results	Deletes the RNase P verification results.
Export results	Exports the results.

7. Unload the plate from the instrument.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the plate temperature can reach 100°C. Allow it to cool to room temperature before handling.

Note: Use Smart Help if the instrument does not eject the plate (see “Request technical support with Smart Help” on page 172).

The status of the RNase P verification is displayed on the **Status and Reminders** screen (see “View the calibration status on the instrument” on page 87).

IMPORTANT! RNase P plates can only be used one time. Do not reuse the RNase P plate.

Troubleshoot verification failure

Observation	Possible cause	Recommended action
Verification failed	The plate was improperly prepared.	<p>Ensure the following:</p> <ul style="list-style-type: none"> • The correct plate was used for the verification performed. • The plate was properly thawed. • The plate was properly centrifuged. • The plate was properly sealed.
		Open the data file for verification in the desktop software or the Connect platform to view the flags and the troubleshooting details for the failed wells.
	The plate is damaged or contaminated.	<p>Check for damage, improper plate seal, or contamination.</p>
High fluorescence signal	The reaction volume is not correct.	Ensure that reaction volumes in the plate are correct and match the volume that is entered in the Method tab.
	Signals that exceed the limit of normal fluorescence can indicate fluorescent contaminants on the plate or on the sample block.	<p>Examine the bottom of the reaction plate. If there is contamination, prepare a new plate and run the experiment again.</p> <p>Identify the location of contamination on the plate or sample block.</p> <ol style="list-style-type: none"> 1. Obtain or prepare a background plate (see “Create a background plate (optional)” on page 96). 2. Follow the procedures to identify contamination (see “Identify contamination” on page 95).
Verification failed but plate is undamaged	The incorrect plate was used for verification.	Use the correct RNase P plate for verification.

Observation	Possible cause	Recommended action
Verification failed but plate is undamaged <i>(continued)</i>	The plate was improperly prepared.	Repeat the verification with a new properly prepared plate. Note: Each RNase P plate can only be used once.
		Open the data file for verification in the desktop software or the Connect platform to view the flags and troubleshooting details for failed wells.
		If the verification fails again, order a replacement plate. Use Smart Help if the replacement plate fails (see “Request technical support with Smart Help” on page 172).

Perform verification using RNase P with a TaqMan™ Array Card

Verification should be performed under the following circumstances:

- After installation and before first use of the instrument and the block
- *(Recommended)* After performing instrument and block calibrations
- As needed to confirm instrument and block performance

The instrument and block require valid ROI/uniformity, background, normalization, and dye calibrations to perform verification.

For a description of verification, see “Instrument verification description” on page 104.

For a description of the pass criteria, see “Performance specifications pass criteria” on page 106.

RNase P instrument verification TaqMan™ Array Card

The ViiA™ 7 Array Card RNaseP Verification Kit includes one empty array card and eight tubes of solution. Each tube contains reaction mix (PCR master mix, RNase P primers, and FAM™-MGB dye-labeled probe) and a known concentration of human genomic DNA template.

To perform an instrument verification run, each solution is loaded into the empty array card. When complete, the array card contains five replicate groups of standards (200, 400, 800, 1600, and 3200 copies), two of unknown populations (800 and 1600 copies), and one that serves as a no template control (NTC).

See Figure 15 on page 110.

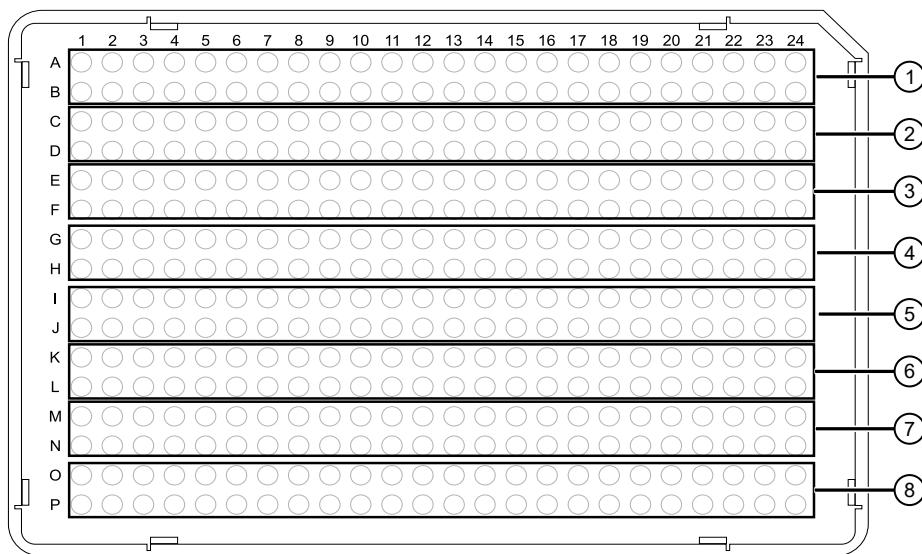


Figure 15 RNase P standards layout

- ① Port 1, No Template Control
- ② Port 2, Unknown A (800 copies)
- ③ Port 3, Unknown B (1,600 copies)
- ④ Port 4, Standard 200 copies
- ⑤ Port 5, Standard 400 copies
- ⑥ Port 6, Standard 800 copies
- ⑦ Port 7, Standard 1,600 copies
- ⑧ Port 8, Standard 3,200 copies

Prepare an RNase P *TaqMan*™ Array Card

Required materials to prepare an RNase P *TaqMan*™ Array Card

Unless otherwise indicated, all materials are available through thermofisher.com. "MLS" indicates that the material is available from fisherscientific.com or another major laboratory supplier.

Item	Source
Calibration kit	
ViiA™ 7 Array Card RNaseP Verification Kit, includes the following components: <ul style="list-style-type: none"> • <i>TaqMan</i>™ Array Card • Eight (8) vials of RNase P detection reagents, one for each port in the <i>TaqMan</i>™ Array Card. Each vial contains the following reagents: <ul style="list-style-type: none"> – PCR master mix – RNase P primers – A FAM™ dye-labeled probe – A known concentration of human genomic DNA template 	4432265

(continued)

Item	Source
Equipment and consumables	
Centrifuge with custom buckets and card holders, one of the following:	
<ul style="list-style-type: none"> • Sorvall™ centrifuge • Multifuge™ centrifuge • Megafuge™ centrifuge 	Contact your local sales office
TaqMan™ Array Card Sealer	Contact your local sales office
Blank balance TaqMan™ Array Cards	Contact your local sales office
Pipette, 200 µL with tips	MLS
Permanent marker or pen	MLS
Powder-free gloves	MLS
Safety glasses	MLS

Prepare an RNase P TaqMan™ Array Card


For detailed instructions to prepare a TaqMan™ Array Card, see Appendix C, “Detailed procedures for preparation of a TaqMan™ Array Card”.

IMPORTANT! Wear powder-free gloves while preparing the array card.

1. Remove the ViiA™ 7 Array Card RNaseP Verification Kit from the freezer, then allow it to warm to room temperature (for approximately 15 minutes).

IMPORTANT! You must fill and run the RNase P array card within 60 minutes of thawing the kit.

2. Remove an array card from its box and place it on a clean, dry surface.
3. Using a permanent marker, mark the side of the empty array card with RNase P.

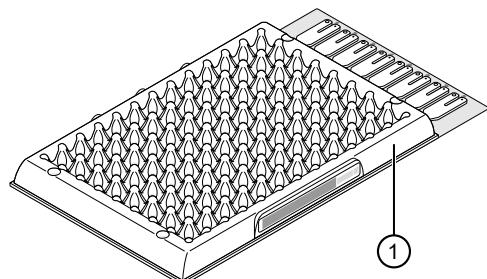

① Location to mark each TaqMan™ Array Card

4. Pipet 100 μ L of each solution into the appropriate port of the array card.

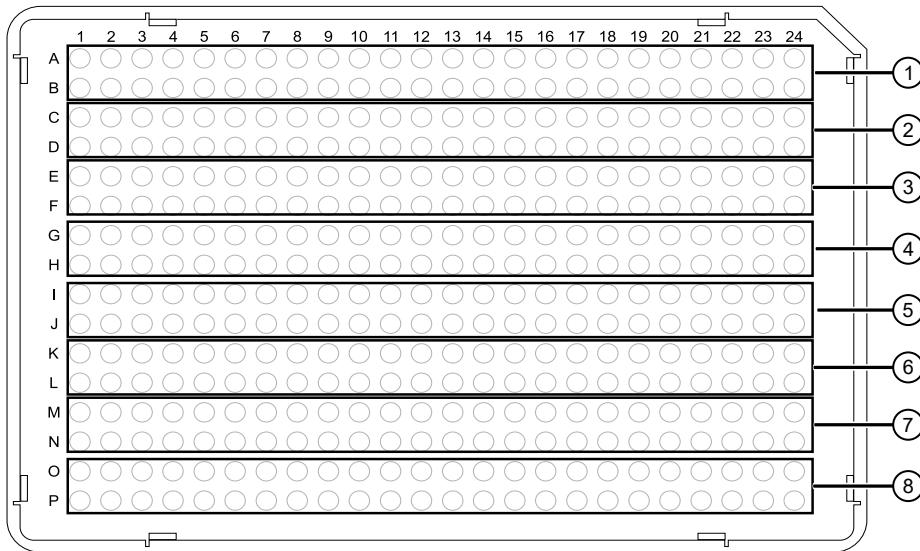


Figure 16 RNase P standards layout

- ① Port 1, No Template Control
- ② Port 2, Unknown A (800 copies)
- ③ Port 3, Unknown B (1,600 copies)
- ④ Port 4, Standard 200 copies
- ⑤ Port 5, Standard 400 copies
- ⑥ Port 6, Standard 800 copies
- ⑦ Port 7, Standard 1,600 copies
- ⑧ Port 8, Standard 3,200 copies

See “Load the PCR reaction mix” on page 205.

5. Centrifuge the array card, then check the array card for the correct fill.

See “Centrifuge the card” on page 205.

6. Seal the array card.

See “Seal the card” on page 207.

IMPORTANT! If an array card is sealed improperly, it can leak and contaminate the sample block and/or it can cause the associated calibration or RNase P test to fail.

7. Run the prepared array card as soon as possible after filling it.

Perform RNase P verification

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **RNase P verification**.
2. Follow the instructions on the screen to start the verification.

3. Load the TaqMan™ Array Card into the instrument.



CAUTION! The instrument should be used by trained operators who have been warned of the moving parts hazard.

4. Tap **Start**.
5. When the run is complete and the screen displays **Complete**, tap **View results** to confirm the status of the run.

Verification status	Action
Passed	Instrument is ready for use.
Failed	See “Troubleshoot verification failure” on page 108.

6. In the **RNase P Verification Status** screen, tap one of the following options.

Option	Description
Accept results	Save the results to the instrument.
Reject results	Deletes the RNase P verification results.
Export results	Exports the results.

7. Unload the TaqMan™ Array Card from the instrument.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the TaqMan™ Array Card temperature can reach 100°C. Allow it to cool to room temperature before handling.

Note: Use Smart Help if the instrument does not eject the plate (see “Request technical support with Smart Help” on page 172).

The status of the RNase P verification is displayed on the **Status and Reminders** screen (see “View the calibration status on the instrument” on page 87).

IMPORTANT! An RNase P TaqMan™ Array Card can only be used one time. Do not reuse the RNase P TaqMan™ Array Card.

Calibrate custom dyes


Overview of custom dyes

The instrument can run assays designed with custom dyes. The following items are defined as custom dyes:

- Dyes that are not manufactured by Thermo Fisher Scientific.
- Dyes or formulations of dyes that are not system dyes for the instrument.

Custom dye calibration workflow

Determine the optimal dye concentration for each custom dye. Use this concentration to prepare all subsequent dye calibration plates.

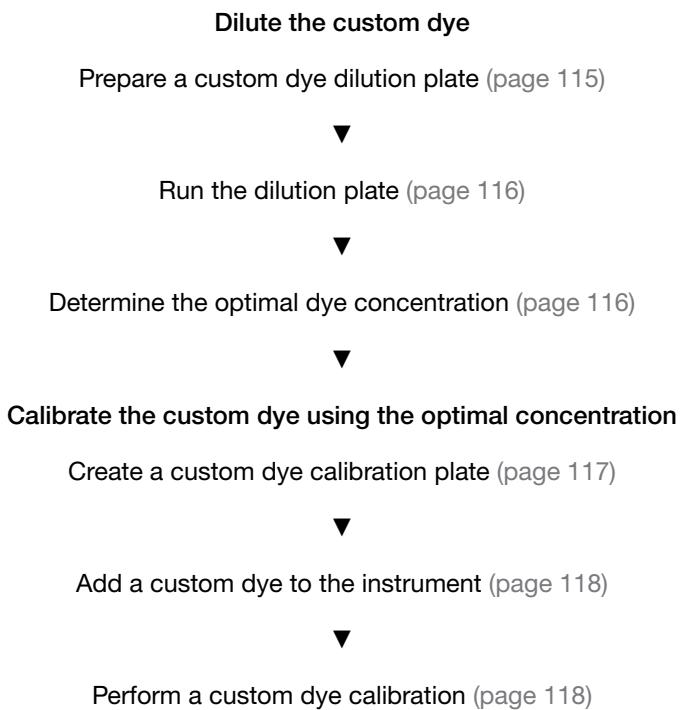

Use a dilution series to determine an optimal custom dye concentration

Custom dye dilution guidelines



Prepare a dilution series for each custom dye.

- Target several dye concentrations within a range of 100–2,000 nM.
- The custom dye for use in calibration must be attached to the 5' end of a short DNA oligonucleotide consisting of the first two bases of the probe sequence without a quencher at the 3' end.
- Choose a 2- or 3-fold difference in dilution points.
- Dispense 10–20 µL per well for 96-well plates.
- Dispense 5 µL per well for 384-well plates.
- Dilute the dye in buffer compatible with your Master Mix.
- (*Intercalating dyes only*) Add the appropriate amount of amplified PCR product to generate fluorescence.


Prepare a custom dye dilution plate

IMPORTANT! Wear powder-free gloves throughout the procedure.

1. Prepare a 2- or 3-fold dilution series of the custom dye.
2. Dispense aliquots of each dilution into the center of a reaction plate, then seal the plate.
A full plate is not needed.

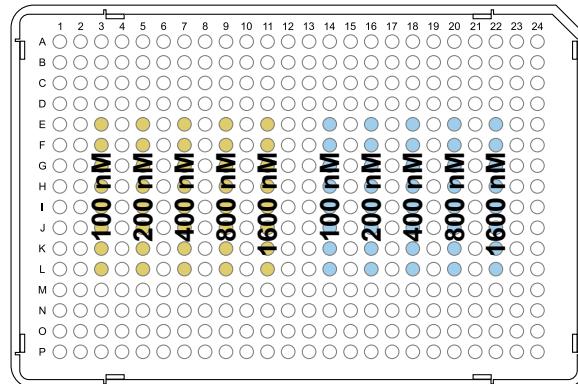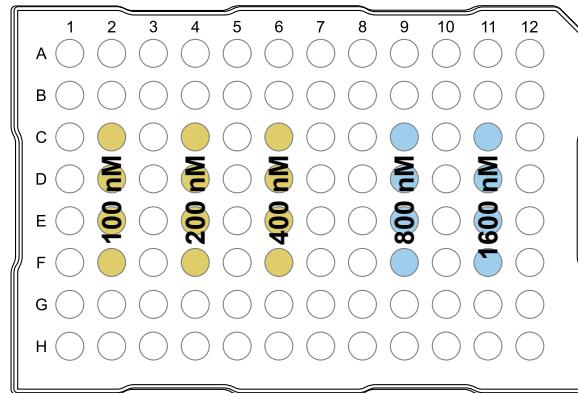

3. Vortex the plate for 5 seconds, then centrifuge at 750–1,000 \times g for 2 minutes.
4. Confirm that the liquid in each well is at the bottom of the well and free of bubbles. If it is not, centrifuge the plate again.

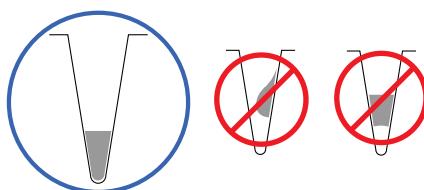

IMPORTANT! Keep the bottom of the plate clean. Fluids and other contaminants on the bottom of the plate can contaminate the sample block and cause an abnormally high background signal.

Run the dilution plate

IMPORTANT! Do not run the dilution plate as a calibration.

1. Load the plate into the instrument.



CAUTION! The instrument should be used by trained operators who have been warned of the moving parts hazard.

2. Set up a genotyping plate file on the instrument (see “Create and run a plate from a system template or a saved plate file” on page 47).
To use a system template for genotyping, see “System template file names” on page 34. Select a genotyping method that is compatible with the Master Mix.
3. In the **Method** tab, set the hold to 60°C for 2 minutes.
4. In the **Plate** tab, enter the dilution series information for the appropriate wells.
5. Tap **Start run**.
6. When the run is complete, transfer the data files for analysis (see “Transfer a data file immediately after a run” on page 81 or “Transfer data files at a later time” on page 81).
7. Unload the plate from the instrument.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the plate temperature can reach 100°C. Allow it to cool to room temperature before handling.

Note: Use Smart Help if the instrument does not eject the plate (see “Request technical support with Smart Help” on page 172).

Determine the optimal dye concentration

Use the Connect platform or the desktop software to review the dye signal data and select the dilution to use for dye calibration.

1. In the **Results** tab, select **Raw data plot**.

This plot displays the raw fluorescence signal of each optical filter, for individual wells.

Note: The raw data plot cannot be viewed on the instrument touchscreen.

2. For each replicate population of dilutions, select the wells in the plate layout to view in the plot.

3. Examine the raw data to identify the wells yielding signals according to the ranges shown in the following table.

Plate type	Acceptable signal range ^[1]
96-well	800,000 to 3,200,000
384-well	400,000 to 2,000,000

^[1] Signal range for the optical filter where the dye is brightest.

4. Examine the raw data to identify the wells yielding signals in the acceptable range for the optical filter where the dye is brightest.
5. (Optional) Export the raw data, then calculate the average fluorescence value for each concentration.
6. Select the lowest optimal dye concentration that falls within the acceptable signal range.


Calibrate the custom dye

Create a custom dye calibration plate

IMPORTANT! Wear powder-free gloves while creating the dye plate.

Create a full plate of the custom dye diluted to the optimal concentration:

1. Dilute the custom dye to the optimal concentration in buffer.
Prepare an adequate volume, using a volume range of 10–20 µL/well for a 96-well plate.
2. Pipet the appropriate volume of the diluted custom dye to the plate wells.
3. Seal the plate.
4. Vortex the plate for 5 seconds, then centrifuge at 750–1,000 \times g for 2 minutes.
5. Confirm that the liquid in each well is at the bottom of the well and free of bubbles. If it is not, centrifuge the plate again.

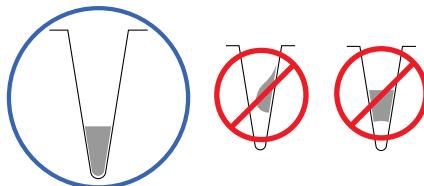

IMPORTANT! Keep the bottom of the plate clean. Fluids and other contaminants on the bottom of the plate can contaminate the sample block and cause an abnormally high background signal.

Add a custom dye to the instrument

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Custom dye**.
The **Custom Dye** screen is displayed.
2. Tap **Add custom dye**.
3. Tap the **Custom Dye Name** field, then enter a name.

IMPORTANT!

- Do not use a system dye name for a custom dye name.
- Dye names are spacing sensitive and cannot contain special characters.

4. Select the type of dye.

Option	Description
Reporter	The dye works in conjunction with a quencher dye to report an increase of PCR product.
Quencher	The dye suppresses the fluorescence of a reporter dye until amplification of PCR product.
Both	The dye can be used as a reporter dye or a quencher dye.

5. Tap **Save**.

IMPORTANT! The custom dye must also be added to the desktop software library before creating, running, or analyzing experiments that use custom dyes.

Perform a custom dye calibration

1. In the instrument home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Custom dye**.

The **Custom Dye** screen is displayed.

2. Tap a custom dye.
3. Review the information, then tap **Update**.
4. Enter the calibration temperature.

Note: The default calibration temperature is set to 60°C.

5. *(Optional)* Tap **Reagents**, then enter the reagent information.
6. Load the plate into the instrument, then tap **Start**.



CAUTION! The instrument should be used by trained operators who have been warned of the moving parts hazard.

Calibration complete is displayed when the run is complete.

7. Tap **View results** ▶ **Details**.

8. Review the plot.

Calibration results that pass show uniform signals with peaks that are aligned with the dye wavelength.

Peak filter	Filter wavelength (nm) ^[1]	
	Excitation	Emission
x1-m1	470 ± 15	520 ± 15
x2-m2	520 ± 10	558 ± 11
x3-m3	550 ± 11	586 ± 10
x4-m4	580 ± 10	623 ± 14
x5-m5	640 ± 10	682 ± 14
x6-m6	662 ± 10	711 ± 12

[1] The central wavelengths are the optimized wavelengths.

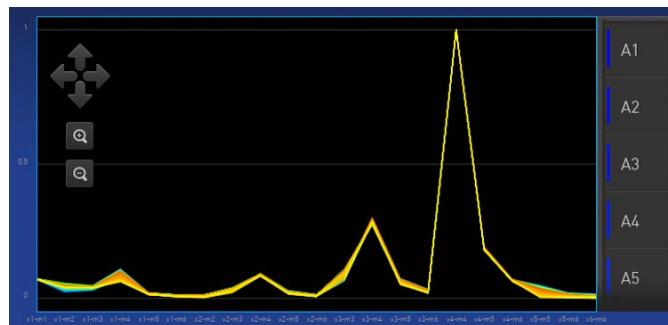


Figure 17 An example of a dye calibration plot

The peaks for your dye may align with a different filter set.

9. Select an action depending on whether the custom dye calibration passed or failed.

Calibration status	Action
Passed	<ol style="list-style-type: none">Tap Accept Results or Reject Results. <p>Note: Accepting the results saves the calibration data to the instrument and overwrites existing data.</p> <ol style="list-style-type: none">(Optional) Tap Transfer to EDS to transfer the calibration data to a USB drive.
Failed	<ol style="list-style-type: none">Perform the calibration again using a new custom dye plate. Create the plate using the next dye concentration greater than the optimal dye concentration (see “Determine the optimal dye concentration” on page 116).For more information about calibration failures, see “Troubleshoot calibration failure” on page 95.

10. Unload the plate from the instrument.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the plate temperature can reach 100°C. Allow it to cool to room temperature before handling.

Note: Use Smart Help if the instrument does not eject the plate (see “Request technical support with Smart Help” on page 172).

Calibrate for a custom melt curve run

Use this calibration to prepare for high resolution melt analysis runs.

Note: A custom melt calibration calibrates a custom dye and a melt calibration at the same time.

- Ensure that all calibrations are current (see “View the calibration status on the instrument” on page 87).
- Add a custom dye (see “Add a custom dye to the instrument” on page 118).

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Calibration** ▶ **Custom melt**.
2. Select the option for the kit that you are using.
 - **PCR + Melt**
 - **Melt only**
3. Select or add a dye, then select a filter set appropriate for your dye's wavelength (see filter-wavelength table below).

Note: See your reagent kit documentation for dye name and wavelength information.

Peak filter	Filter wavelength (nm) ^[1]	
	Excitation	Emission
x1-m1	470 ± 15	520 ± 15
x2-m2	520 ± 10	558 ± 12
x3-m3	550 ± 10	587 ± 10
x4-m4	580 ± 10	623 ± 14
x5-m5	640 ± 10	682 ± 14
x6-m6	662 ± 10	711 ± 12

^[1] The central wavelengths are the optimized wavelengths.

IMPORTANT! If the selected filter set does not match your reagent kit documentation, then the incorrect wavelength may be collected during a run.

4. *(Optional)* Tap **Reagents**, then enter the reagent information.

5. Load the plate into the instrument
6. Tap **Start**.
7. When the run is complete and the screen displays **Calibration complete**, tap **View results** ▶ **Details**.
8. Review the plot.

Passing calibration results show uniform signals with peaks that are aligned with the dye wavelength.

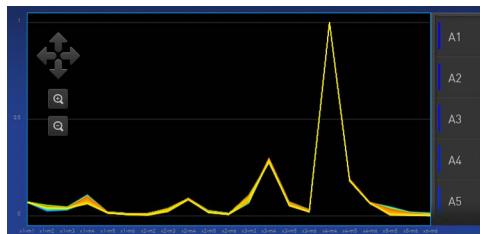


Figure 18 An example dye calibration plot

Note: The peaks for your dye may align with a different filter set.

9. Select an action depending on whether the custom dye calibration passed or failed.

Calibration status	Action
Passed	<ol style="list-style-type: none">1. Tap Accept Results or Reject Results. Note: Accepting the results saves the calibration data to the instrument and overwrites existing data.2. (Optional) Tap Transfer EDS to transfer the calibration data to a USB drive.
Failed	<ol style="list-style-type: none">1. Perform the calibration again using a new custom dye plate. Create the plate using the next dye concentration greater than the optimal dye concentration (see “Determine the optimal dye concentration” on page 116).2. For more information, see “Troubleshoot calibration failure” on page 95.

10. Unload the plate from the instrument.

Note: You must also add the custom dye to the desktop software dye library before you create, run, or analyze experiments that use custom dyes.

Note: To perform a custom melt experiment, you can either create a Standard Curve or a Custom experiment with melt, then specify the data points per degree in the method.

■ Require sign-in	122
■ Recommended order to set up profiles	123
■ Overview of local instrument profiles and Connect profiles	123
■ Connect instrument profile roles and functions	124
■ Test the connection to the Connect platform	124
■ Link the instrument to your Connect account	126
■ Create a new local instrument profile	127
■ Link a local profile to a Connect profile	127
■ Unlink a Connect account	129
■ If you link when you are signed in to the instrument	129
■ If you link when you are not signed in to the instrument	130
■ Manage individual instrument profiles	131
■ Manage all instrument profiles as an administrator	133

Require sign-in

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Sign-in required**.
The **Sign-In Required** screen is displayed.
2. Set the slider.

Option	Description
Off	The instrument will allow guest users to access public folders.
On	The instrument will require a user to sign in.

3. Tap **Done**.

Recommended order to set up profiles

Task	Description
Create a profile for the instrument administrator.	<p>The first profile on the instrument becomes the instrument administrator.</p> <p>The first profile is set up during the initial install with a representative from Thermo Fisher Scientific.</p> <p>See “Create an administrator instrument profile during initial start-up” on page 131.</p>
Select a Connect administrator.	<p>A Connect administrator is distinct from an instrument administrator.</p> <p>The first Connect profile to be linked to the instrument is automatically assigned the role of a Connect administrator. It is recommended that the first Connect profile to be linked is the one that should be a Connect administrator.</p> <p>The roles can be updated at a later time on the Connect platform.</p>
Each person creates a profile.	<p>Create a local instrument profile. Connect-enabled features are not available.</p> <p>(Optional) Link a local instrument profile to a Connect profile.</p>
Assign additional instrument administrators.	<p>Sign in with a Connect profile. Connect-enabled features are available.</p> <p>Any of the profiles can be assigned the role of instrument administrator. The local profile must be created first or the Connect profile must be linked first.</p>

Overview of local instrument profiles and Connect profiles

- Local instrument profile
 - Plate files and data files are stored on the instrument.
 - Plate files and data files can be transferred to the desktop software.
 - Voice commands are not available.
 - Email notifications are not available.
 - The instrument schedule is not available.
 - Smart Help is not available.
 - Separate profiles are required for each instrument if multiple instruments are used.
- Connect profile
 - Plate files and data files are stored on the instrument.
 - Plate files and data files can be transferred to the Connect platform or the desktop software.
 - Voice commands are available.
 - Email notifications are available.

- The instrument schedule is available.
- Smart Help is available.
- The same profile can be used for multiple instruments.


Connect instrument profile roles and functions

The first user who links their local instrument profile to their Connect account is assigned a Connect profile with administrator role.

Instrument profile	Location	Functions allowed
Standard	Connect	<ul style="list-style-type: none"> • Create, save, open, import, and run plate files • Create and modify run settings • View and export data files
Administrator	Connect	<p>All the permissions of a local administrator profile, plus the following functions performed in Connect:</p> <ul style="list-style-type: none"> • See a list of all the Connect profiles that are linked to the instrument • Assign Connect administrator roles to one or more users • Remove a user from an instrument • Disconnect the instrument from Connect • Change the instrument name


Test the connection to the Connect platform

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Connect services**.
 The **Connect Services** screen is displayed.

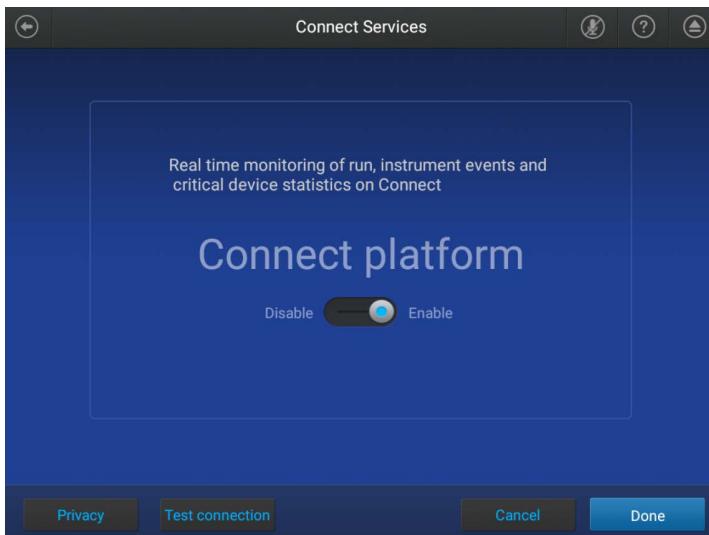


2. Tap **Test connection**.

If the connection can be established, **You are able to connect to the Connect platform** will be displayed.

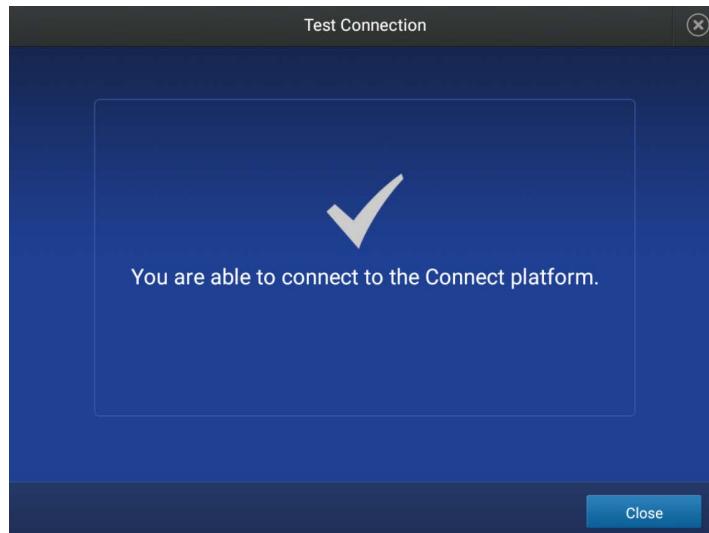

If the connection cannot be established, **Unable to connect** will be displayed.

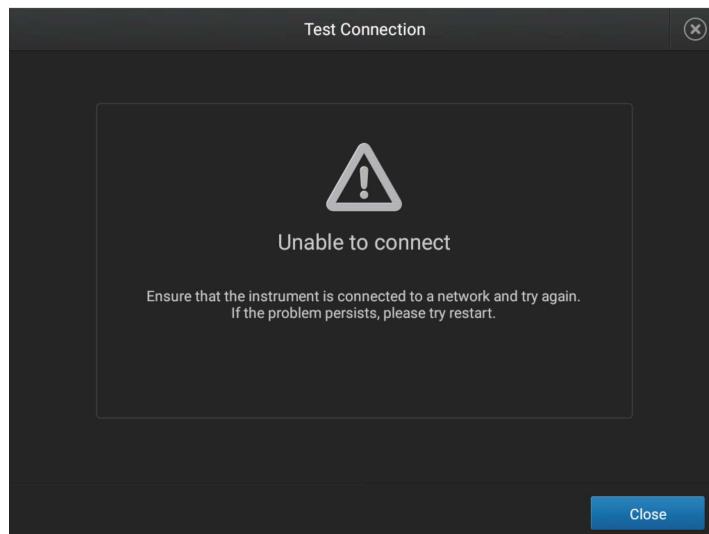

3. Tap **Close**.

Link the instrument to your Connect account

This section describes using your Connect account when you use the instrument for the first time.

The first time you use your Connect account on an instrument, you will be prompted to create a four-digit numerical PIN. This PIN is to use when signing in to the instrument with your Connect account. It will apply to all other instruments when you use your Connect account. This does not change the password when signing in to your Connect account on a browser.

After the PIN is set up on the instrument, it must be changed with your Connect account on a browser (see “Change the PIN for a Thermo Fisher™ Connect profile” on page 133).

The first time the instrument is used with a Connect account, a region must be selected.

For a detailed description of the profile, see “If you link when you are not signed in to the instrument” on page 130.

1. In the **Sign In** screen, tap **Get started ▶ Connect**.
2. *(Optional)* Select the appropriate region.

Option	Description
China	For users in China
U.S.	For users in any country other than China

3. Tap a connection option.

Option	Action
 Mobile devices	<p>Note: Before selecting this option, install and sign in to the Connect application on your mobile device.</p> <p>On the instrument:</p> <ol style="list-style-type: none"> 1. Tap  Mobile devices. 2. Hold the camera on your mobile device over the QR code that is displayed on the touchscreen. 3. Tap Close.
 PC	<p>A link code is displayed on the instrument.</p> <p>On a computer:</p> <ol style="list-style-type: none"> 1. Access the Connect platform. 2. Click Add instrument. 3. Select QuantStudio. 4. Enter the link code.

4. *(Optional)* In the **Enter PIN** screen, tap the **PIN (4 digits required)** field, enter a four-digit numerical PIN, then tap **Enter**.

Tap the **Show PIN** checkbox to show or hide the PIN.

5. *(Optional)* Tap the **Confirm PIN** field, enter the four-digit numerical PIN again, then tap **Enter**.
6. Tap **Done**.

Create a new local instrument profile

A local instrument profile is not compatible with Connect-enabled features. A local instrument profile can be linked to Connect at a later time.

To link a local instrument profile to Connect at a later time, see “Link a local profile to a Connect profile” on page 127.

For a description of linking a local instrument profile to Connect, see “If you link when you are signed in to the instrument” on page 129.

1. In the **Sign In** screen, tap **Get Started**.
2. Tap **Create Profile**.
3. Tap the **User name** field, enter a user name, then tap **Enter**.
4. Tap the **PIN (4 digits required)** field, enter a four-digit numerical PIN, then tap **Enter**.
Tap the **Show PIN** checkbox to show or hide the PIN.
5. Tap the **Confirm PIN** field, enter the four-digit numerical PIN again, then tap **Enter**.
6. Tap **Create Profile**.

Proceed immediately to set up facial authentication (see “Set up facial authentication at the initial creation of an instrument profile” on page 131) or tap **Skip** to set up facial authentication at a later time.

Link a local profile to a Connect profile

This section describes linking a local instrument profile to a Connect profile, if a local instrument profile was created first.

The first time you use your Connect account on an instrument, you will be prompted to create a four-digit numerical PIN. This PIN is to use when signing in to the instrument with your Connect account. It will apply to all other instruments when you use your Connect account. This does not change the password when signing in to your Connect account on a browser.

After the PIN is set up on the instrument, it must be changed with your Connect account on a browser (see “Change the PIN for a Thermo Fisher™ Connect profile” on page 133).

The first time the instrument is used with a Connect account, a region must be selected.

For a detailed description of the profile, see “If you link when you are signed in to the instrument” on page 129.

(Optional) Test the connection (see “Test the connection to the Connect platform” on page 124).

1. In the home screen, tap  **(Profile)**.

The **My Profile** screen is displayed.

2. Tap **Edit ▶ Link**.

The **Connect to Connect Platform** screen is displayed.

3. (Optional) Select the appropriate region.

Option	Description
China	For users in China
U.S.	For users in any country other than China

4. Tap a connection option.

Option	Action
 Mobile devices	<p>Note: Before selecting this option, install and sign in to the Connect application on your mobile device.</p> <p>On the instrument:</p> <ol style="list-style-type: none"> 1. Tap  Mobile devices. 2. Hold the camera on your mobile device over the QR code that is displayed on the touchscreen. 3. Tap Close.
 PC	<p>A link code is displayed on the instrument.</p> <p>On a computer:</p> <ol style="list-style-type: none"> 1. Access the Connect platform. 2. Click Add instrument. 3. Select QuantStudio. 4. Enter the link code.

5. (Optional) In the **Enter PIN** screen, tap the **PIN (4 digits required)** field, enter a four-digit numerical PIN, then tap **Enter**.

Tap the **Show PIN** checkbox to show or hide the PIN.

6. (Optional) Tap the **Confirm PIN** field, enter the four-digit numerical PIN again, then tap **Enter**.

7. Tap **Done**.

Unlink a Connect account

Unlinking a Connect is done from the Connect platform on your computer.

For more information about unlinking a Connect account, see “If you link when you are signed in to the instrument” on page 129 and “If you link when you are not signed in to the instrument” on page 130.

1. Sign in to your account on the desktop Connect platform.
2. In the left pane, click  **(Instrument)**.
3. Select the instrument, then click  **Disconnect**.
4. Tap **Confirm**.

If you link when you are signed in to the instrument

In this scenario, your local instrument profile name is created manually on the instrument before you link. Your local instrument profile name differs from your Connect instrument profile name.

Phase	Steps that occur
Before you link:	<ul style="list-style-type: none">• You enter your local instrument profile name in the Sign In screen.• Your local instrument profile (<i>UserABC</i>) is displayed in the home screen of the instrument.• All plates and results that you create are accessible only when you are signed in with your local instrument profile.
When you link:	<ul style="list-style-type: none">• You link your local instrument profile (see “Link a local profile to a Connect profile” on page 127).• If this is the first time you link, a Connect instrument profile is created using the <i>FirstNameLastInitial</i> of the user name from your thermofisher.com account. Example: <i>User1@thermofisher.com</i> First name is <i>User</i>, Last name is <i>Gray</i>. Connect account username is <i>User G.</i>• Your local instrument profile (<i>UserABC</i>) is linked to your Connect account (<i>User1@thermofisher.com</i>).• Your Connect instrument profile (<i>User G.</i>) replaces your local instrument profile.
After you link:	<ul style="list-style-type: none">• Your Connect instrument profile (<i>User G.</i>) and  is displayed in the home screen of the instrument.• Plate files and data files from your local instrument profile can be transferred to Connect.• New plate files and data files are saved under your Connect instrument profile.• Your Connect instrument profile name (<i>User G. </i>) is available for selection in the Sign In screen.

(continued)

Phase	Steps that occur
If your Connect account is unlinked:	<ul style="list-style-type: none"> Your local instrument profile (<i>UserABC</i>) is displayed in the home screen of the instrument. Plate files and data files that were saved under your Connect instrument profile are accessible under your local instrument profile. Plate files and data files are saved under local instrument profile and can be copied to Connect (see “Copy plate files” on page 77 and “Transfer data files at a later time” on page 81). Your local instrument profile name (<i>UserABC</i>) is available for selection in the Sign In screen.

If you link when you are *not* signed in to the instrument

In this scenario, your local instrument profile name is created automatically at the time that you link. The same user name is used for your local instrument profile and your Connect profile. Plates and results are accessible when you sign in with either profile.

If you link to Connect when you *are not* signed in to the instrument:

Phase	Steps that occur
Before you link:	<ul style="list-style-type: none"> In the Sign In screen, you tap Get Started ▶ Connect.
When you link:	<ul style="list-style-type: none"> You link your profile (see “Link the instrument to your Connect account” on page 126). If this is the first time you link, a Connect instrument and a local instrument profile (with standard role) are created with the same name using the <i>FirstNameLastInitial</i> of the user name from your <i>thermofisher.com</i> account. Example: <i>User1@thermofisher.com</i> First name is <i>User</i>, Last name is <i>Gray</i>. Connect account username is <i>User G</i>. Your local instrument profile (<i>User G.</i>) is linked to your Connect account (<i>User1@thermofisher.com</i>). Your Connect instrument profile (<i>User G.</i>) replaces your local instrument profile.
After you link:	<ul style="list-style-type: none"> Your Connect instrument profile (<i>User G.</i>) and  is displayed in the home screen of the instrument. Plate files and data files from your local instrument profile can be transferred to Connect. New plate files and data files are saved under your Connect instrument profile. Your Connect instrument profile name (<i>User G.</i> ) is available for selection in the Sign In screen.

(continued)

Phase	Steps that occur
If your Connect account is unlinked:	<ul style="list-style-type: none">• Your local instrument profile (<i>User G.</i>) is displayed in the home screen of the instrument.• Plate files and data files that were saved under your Connect instrument profile are accessible under your local instrument profile.• New plates and results are saved under local instrument profile and can be copied to Connect.• Your local instrument profile name (<i>User G.</i>) is available for selection in the Sign In screen.

IMPORTANT! If you sign in with a local profile, without linking to the Connect platform, sign out, then link to Connect (**Get Started ▶ Connect**), you can potentially have two instrument profiles with different names. Plate files and data files from when you are signed in with one instrument profile are not accessible when you are signed in with the other instrument profile.

Manage individual instrument profiles

Create an administrator instrument profile during initial start-up

The first instrument profile that is created during installation is assigned the role of an instrument administrator. Administrators can grant administrative privileges to other users (see “Enable or disable administrator privileges” on page 135).

Note: An instrument administrator is separate from a Connect administrator (see “Recommended order to set up profiles” on page 123).

1. Tap **Name**, enter a user name, then tap **Done**.
2. Tap **PIN**, enter a four-digit numerical PIN, then tap **Enter**.

Note: Tap the **PIN** checkbox to switch the PIN display on or off.

3. Tap **Confirm PIN**, then enter the four-digit numerical PIN again.
4. Tap **Create Profile**.

Set up facial authentication at the initial creation of an instrument profile

When a profile is set up for facial authentication, the photographs should reflect how the user appears in the lab. For example, safety glasses should be worn to set up facial authentication if safety glasses will be worn when signing in.

1. On the **Instrument Profile** page, tap **Set up now**.
2. Prepare for the photographs, then tap **Next**.

3. Center your face within the green frame, then tap **Start**.

Note: You must position yourself close enough to the instrument so that the green frame shows only your face without any of the surrounding background. This position is closer to the instrument than when you sign in with facial authentication.

The instrument will take three separate photos after you tap **Capture**. The instrument will display a countdown before each individual photo is taken.

4. Tap one of the following options.

- **Done**
- **Re-take**

5. Tap **Done** after **Your instrument profile has been created** is displayed.

Add facial authentication to an existing instrument profile

When a profile is set up for facial authentication, the photographs should reflect how the user appears in the lab. For example, safety glasses should be worn to set up facial authentication if safety glasses will be worn when signing in.

1. Sign in.

See “Sign in with a PIN” on page 43.

2. In the home screen, tap  **(Profile)** ▶ **Facial profile** ▶ **Set up now**.

Note: The instrument must be configured for facial authentication (see “Configure the instrument for the hands-free features” on page 185). The **Facial profile** button will not display if the instrument is not configured for facial authentication.

3. Prepare for the photographs, then tap **Next**.

4. Center your face within the green frame, then tap **Start**.

Note: You must position yourself close enough to the instrument so that the green frame shows only your face without any of the surrounding background. This position is closer to the instrument than when you sign in with facial authentication.

The instrument will take three separate photos after you tap **Capture**. The instrument will display a countdown before each individual photo is taken.

5. Tap one of the following options.

- **Done**
- **Re-take**

Change the PIN for an instrument profile

This section describes changing a PIN for a local instrument profile. To change a PIN for a Connect profile, see “Change the PIN for a Thermo Fisher™ Connect profile” on page 133.

1. In the home screen, tap  **(Profile)**.
2. Tap **Edit**.
3. Tap the **Old PIN** field, enter the current four-digit numerical PIN, then tap **Enter**.
4. Tap the **PIN (4 digits required)** field, enter a new four-digit numerical PIN, then tap **Enter**.
Tap the **Show PIN** checkbox to show or hide the PIN.
5. Tap the **Confirm PIN** field, enter the new four-digit numerical PIN again, then tap **Enter**.
6. Tap **Done**.

Change the PIN for a Thermo Fisher™ Connect profile

This section describes changing a PIN for a Thermo Fisher™ Connect profile. To change a PIN for a local instrument profile, see “Change the PIN for an instrument profile” on page 133.

1. Sign in to Thermo Fisher™ Connect on a browser.
2. Click **Update PIN number**.
3. In the **Update PIN number** dialog box, enter a new PIN, then enter the new PIN a second time to confirm it.
4. Click **Send**.

The 4-digit PIN to sign in to instruments with your Thermo Fisher™ Connect profile is updated. This change applies to all of the instruments that you access with your Connect profile.

Manage all instrument profiles as an administrator

View all instrument profiles

1. In the home screen, tap  **(Profile)**.
The **My Profile** screen is displayed.
2. Tap **All accounts**.
A list of all the profiles is displayed. The following information is included for each profile.

• Name	• Role (standard or administrator)
• Date and time created	• Location (local or Connect)

3. *(Optional)* Tap a column in the table to sort by that parameter.
4. Tap **Done**.

Reset the PIN for a profile

1. In the home screen, tap  **(Profile)**.
The **My Profile** screen is displayed.
2. Tap **All accounts**.
A list of all the profiles is displayed.
3. Tap a profile.
The **Manage Account** screen is displayed.
4. Tap **Delete PIN**, then tap **Yes**.
5. Tap **Done** to return to the **My Profile** screen.

The profile must have a new PIN added the next time it is accessed.

Add a profile

1. In the home screen, tap  **(Profile)**.
The **My Profile** screen is displayed.
2. Tap **All accounts**.
3. Tap **Add profile**.
4. Tap the **User name** field, enter a user name, then tap **Enter**.
5. Tap the **PIN (4 digits required)** field, enter a four-digit numerical PIN, then tap **Enter**.
Tap the **Show PIN** checkbox to show or hide the PIN.
6. Tap the **Confirm PIN** field, enter the four-digit numerical PIN again, then tap **Enter**.
7. Tap **Create profile**.
The new profile is displayed on the **My Profile** screen. The new profile is assigned a standard role.

To enable administrator privileges, see “Enable or disable administrator privileges” on page 135.

Enable or disable administrator privileges

Administrator privileges can only be enabled or disabled on the instrument for local instrument profiles. Connect instrument profiles can only have instrument administrator privileges enabled or disabled on the Connect platform.

1. In the home screen, tap  **(Profile)**.
The **My Profile** screen is displayed.
2. Tap **All accounts**.
3. Tap a profile.
The **Manage Account** screen is displayed.
4. Set the **Grant admin rights** slider to **No** or **Yes**.
5. Tap **Done**.
The user role will be updated on the **My Profile** screen.

Delete a profile

1. In the home screen, tap  **(Profile)**.
The **My Profile** screen is displayed.
2. Tap **All accounts**.
3. Tap a profile.
The **Manage Account** screen is displayed.
4. Tap **Delete account**.
5. Tap **Yes** to confirm.
The profile is deleted and it is not listed on the **My Profile** screen.

Use the instrument with the Security, Auditing, and E-signature (SAE) Administrator Console

■ Overview of the SAE Administrator Console components	137
■ Overview of the instrument features when the SAE module is enabled	137
■ SAE Administrator Console functions	138
■ SAE Administrator Console roles	139
■ Dialog boxes in the SAE module	140
■ Profiles when the SAE module is enabled	141
■ SAE-enabled system components	141
■ Enable SAE functions	143
■ Disable SAE on the instrument (administrator only)	145
■ Sign in as a local administrator with SAE enabled	146
■ Sign in with SAE enabled	147
■ Sign out with SAE enabled	148
■ Change your SAE account password on the instrument	149
■ Sign a plate setup	150
■ Enter an audit reason	151
■ View signing records for a plate file	152
■ View audit records	153
■ Use the instrument when the SAE server is offline	154
■ E-signature requirements to start a run	154
■ Signing in after automatic screen locking	155
■ SAE error messages and actions	155

The SAE Administrator Console is only compatible with the QuantStudio™ 7 Pro Real-Time PCR System.

The Security, Auditing, and E-signature (SAE) v2.1 module or later is compatible.

This version converts the URL to **HTTPS** when accessing the SAE Administrator Console. For more information, see the *SAE Administrator Console v2.0 User Guide* (Pub. No. MAN0017468).

The Security, Auditing, and E-signature (SAE) v2.2 includes a language selection.

For more information about the module, including definitions of accounts and roles, see the *SAE Administrator Console v2.0 User Guide* (Pub. No. MAN0017468).

Overview of the SAE Administrator Console components

The SAE Administrator Console includes three components:

- SAE Administrator Console that an administrator uses to configure the module.
- SAE server that stores settings, user accounts, audit records, and e-signature records. By default, the SAE server is installed on the same computer as the SAE Administrator Console.
- SAE screens in an application (sign in and audit that a user interacts with). The QuantStudio™ 7 Pro Real-Time PCR System is an application.

The SAE Administrator Console provides the following SAE functionality on the instrument:

- **System security**—Controls user sign in and access to functions
- **Auditing**—Tracks changes and actions performed by users.
- **E-signature**—Allows users to provide an electronic signature (user name and password) when performing certain functions.

Overview of the instrument features when the SAE module is enabled

The following instrument features are not available when the SAE module is enabled:

- Facial authentication
- Voice commands
- Linking to the Connect cloud-based platform, including using a Connect cloud-based platform account to log in
- System templates

Only SAE-enabled plate files that were created in the QuantStudio™ Design and Analysis Software v2 can be opened on the instrument.

SAE Administrator Console functions

Functions that are controlled

The following functions are controlled, depending on the user role:

- Plate setup, including the following functions:
 - Create a new template
 - Edit the run method
 - Edit the analysis settings
 - Define, edit, or delete the targets or the assays
 - Assign the target or the assay
 - Define, edit, or delete the sample
 - Assign the samples
 - Define, edit, or delete the reagents
 - Assign the reagents
- Add and delete a custom dye
- Start and stop an instrument run
- Calibration runs and RNase P verification runs
- Instrument configuration, including the following functions:
 - Software update
 - Network configuration
 - Instrument name change
 - Date and time change
 - SAE module setting modification
 - Sleep setting modification
 - Instrument access modification
 - Disk management setting modification
 - Instrument back-up and restore a back-up
- Log into timed out sessions
- Perform e-signatures

Functions that can be audited

Certain instrument functions can be audited. This depends on how the SAE administrator has configured the audit settings.

The following functions can be audited:

- Changes to a plate file, including the following items:
 - Plate layout
 - Analysis settings
 - Run method
 - Samples
 - Assays
- Sign in success and sign in failure
- Sign out
- Enable and disable the SAE Administrator Console
- Perform a calibration run
- Perform an RNase P verification run
- Start and stop an instrument run
- Pause and resume an instrument run

Functions that can be signed

Certain instrument functions can be signed. This depends on how the SAE administrator has configured the e-signature settings.

The following functions can be signed with an e-signature:

- Start an instrument run
- Plate setup
- Plate results



SAE Administrator Console roles

Role name	Description
Administrator	Has all the available privileges.
Scientist	Has all the privileges, except for privileges related to instrument configuration and maintenance.
Technician	Has access to the instrument to perform a run.

Dialog boxes in the SAE module


Some of the features and functions described might not be accessible to you. The features and functions that are available depend on the way that the SAE administrator has configured them.

When you use the software, the dialog boxes and prompts shown in the following table can appear.

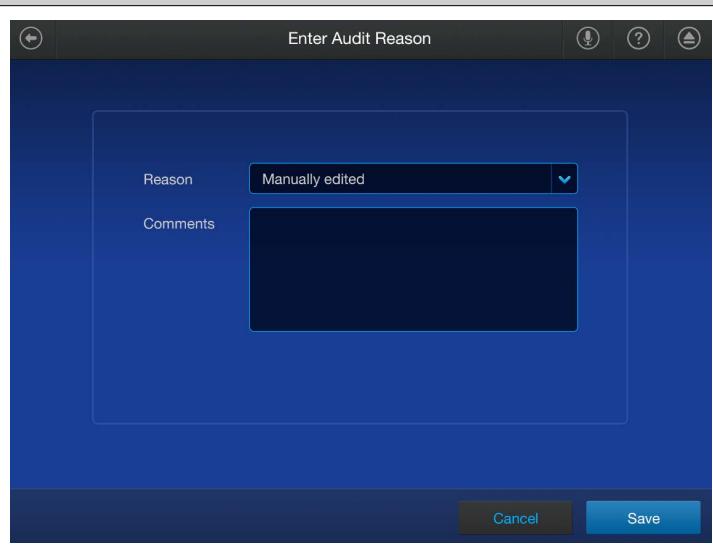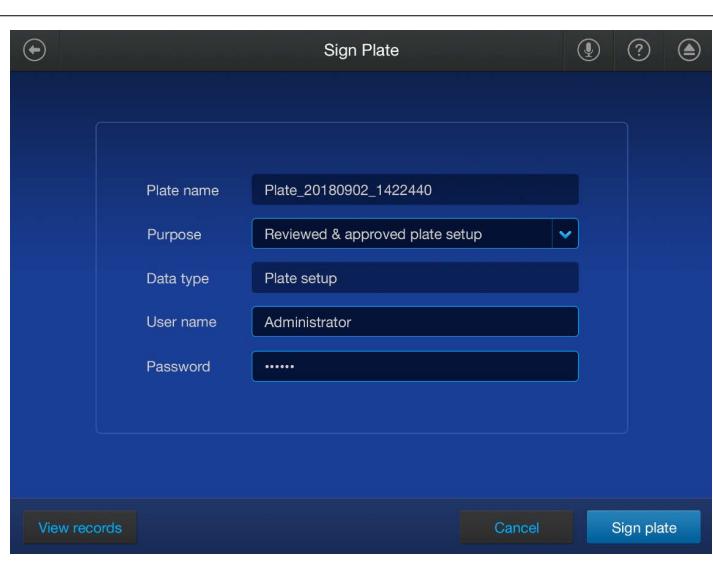Dialog box	Description
	<p>An action is set up for auditing and requires you to specify a reason for the action.</p>
	<p>An action is set up for electronic signature and allows you to enter your password to allow the action.</p>

If you do not have access to a function, you will not be able to select it.

For example, the buttons to start any of the calibrations are inactive if your role is not able to perform a calibration.

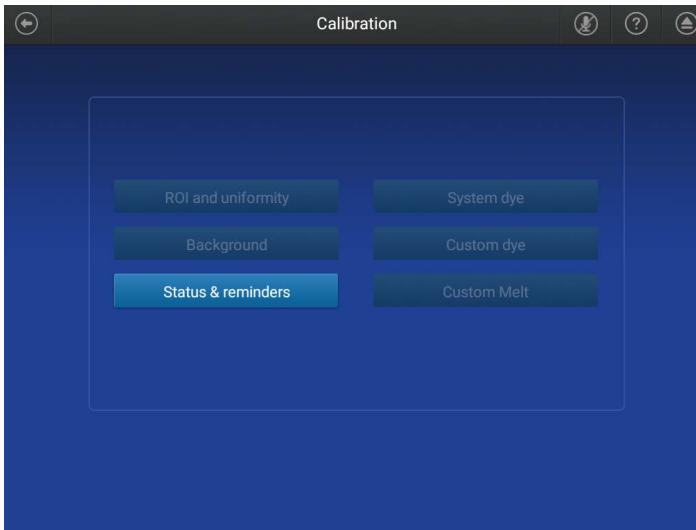

Profiles when the SAE module is enabled

After the SAE module is enabled on the instrument, the local instrument profiles and the Connect cloud-based platform profiles will not be available.

An account from the SAE Administrator Console must be used to sign in to the instrument when the SAE module is enabled. A local instrument administrator can sign in to the instrument to perform limited administrator functions.

The local instrument profiles and the Connect cloud-based platform profiles are available if the SAE module is disabled.

SAE-enabled system components

The following system components can be used with SAE functions enabled:

- QuantStudio™ Design and Analysis Software v2
- QuantStudio™ 7 Pro Real-Time PCR System
- QuantStudio™ 7 Pro Real-Time PCR System plate file—SAE-enabled plate files are created in the QuantStudio™ Design and Analysis Software v2 with SAE enabled.
 - If SAE is enabled in the QuantStudio™ Design and Analysis Software v2, you can only create plate files for the QuantStudio™ 7 Pro Real-Time PCR System.
 - If SAE is enabled on the QuantStudio™ 7 Pro Real-Time PCR System, you cannot create or edit a plate file from the instrument touchscreen.
- QuantStudio™ 7 Pro Real-Time PCR System data file—SAE-enabled data files are created on the QuantStudio™ 7 Pro Real-Time PCR System with SAE enabled.

We recommend enabling SAE for all system components (see “Enable SAE functions” on page 143). If one or more of the components have conflicting SAE status, some functions may not be available (see “Compatibility between SAE-enabled and SAE-disabled components” on page 142).

Compatibility between SAE-enabled and SAE-disabled components

We recommend enabling SAE for all system components (for more information, see “Enable SAE functions” on page 143). If one or more of the components have conflicting SAE status, some functions may not be available. See the following table for more information.

Component	Functionality with an SAE-enabled plate or data file	Functionality with an SAE-disabled plate or data file
QuantStudio™ Design and Analysis Software v2 with SAE enabled	<ul style="list-style-type: none"> The file can be edited depending on SAE configuration. The audit trail is continued. 	<ul style="list-style-type: none"> The file is opened in read-only mode. The file cannot be edited or saved.
QuantStudio™ Design and Analysis Software v2 with SAE disabled	<ul style="list-style-type: none"> SAE-disabled files allowed—The file is opened and can be edited. The file can be saved as an invalid SAE file only.^[1] QuantStudio™ 7 Pro Real-Time PCR System forbidden—The file cannot be opened. 	The file can be opened, edited, and saved.
QuantStudio™ 7 Pro Real-Time PCR System with SAE enabled	<ul style="list-style-type: none"> The file can be opened from the run queue, a USB drive, or other sources.^[2] The file cannot be edited. The audit record is continued. 	The file cannot be opened.
QuantStudio™ 7 Pro Real-Time PCR System with SAE disabled	<ul style="list-style-type: none"> The plate file can be opened and edited. The file can be saved as an invalid SAE file only.^[1] The file can be used to start a run, but the data file will be an invalid SAE file.^[1] 	The file can be opened, edited, and saved.

^[1] Invalid SAE files contain incomplete audit records.

^[2] You cannot import plate files from the Connect cloud-based platform when the instrument has SAE enabled.

Enable SAE functions

Workflow: Enable SAE functions

Configure SAE Administrator Console application profiles (page 143)



Enable SAE on the instrument and specify the SAE server (administrator only) (page 143)



Connect to the SAE server (page 144)



Enable SAE functions in QuantStudio™ Design and Analysis Software v2 (page 145)

Configure SAE Administrator Console application profiles

Note: Configuring application profiles in the SAE Administrator Console requires an SAE administrator account.

In the SAE Administrator Console, an application profile contains default settings for an application. Before using the SAE Administrator Console, an administrator must install, then configure profiles for the following applications:

1. QuantStudio™ 7 Pro Real-Time PCR System
2. QuantStudio™ Design and Analysis Software v2

Note: The QuantStudio™ 7 Pro Real-Time PCR System profile must be installed before the QuantStudio™ Design and Analysis Software v2 profile.


For information on configuring application profiles, see *SAE Administrator Console v2.0 User Guide* (Pub. No. MAN0017468).

Enable SAE on the instrument and specify the SAE server (administrator only)

This procedure requires a local administrator profile on the instrument and an SAE administrator account in the SAE Administrator Console.

1. In the home screen, tap  **(Settings)** ▶ **SAE**.
The **SAE Mode** screen is displayed.
2. In the **SAE Mode** screen, set the **SAE Mode** slider to **Enable**.

3. Tap the **Server IP** field, then enter the IP address of the SAE server.

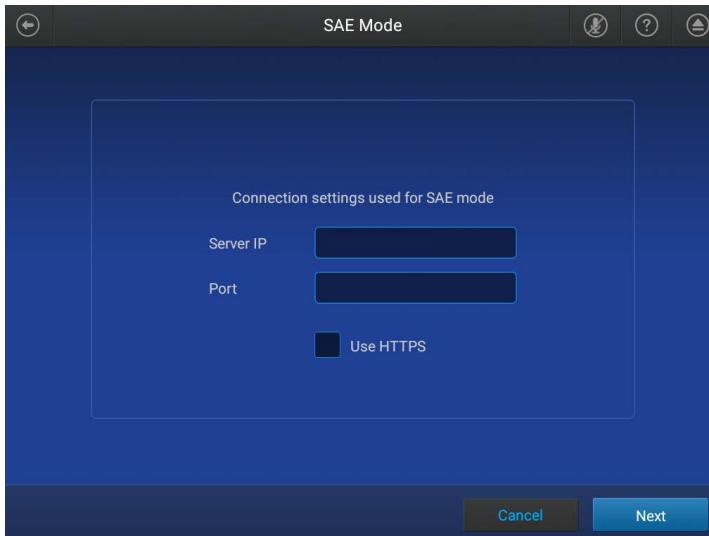

4. Tap the **Port** field, enter the port, then tap **Next**.
5. Enter the SAE administrator user name and password when prompted, then tap **Enable**.

The home screen is displayed. The SAE administrator is signed in.

Connect to the SAE server

- Install the SAE Administrator Console and the SAE server on a computer with a static IP address (*recommended*) or a dynamic IP address.
- Before you connect, close all plate files and data files.

Note: Connect the software and any instruments to the same SAE server to ensure that audit records are maintained across system components.

1. In the menu bar, click **System** ▶ **SAE Connection Settings**.

2. Enter the IP address and port number of the SAE server.

Note: If using a dynamic IP address, enter the hostname instead of the IP address to prevent the loss of a connection.

3. *(Optional)* Click **Test Connection** to confirm that the connection information is correct.

4. Click **Save**.

Enable SAE functions in QuantStudio™ Design and Analysis Software v2

This procedure requires an SAE administrator account.

Before you enable SAE functions in the QuantStudio™ Design and Analysis Software v2, you must complete the following tasks:

- Connect to the SAE server (see “Connect to the SAE server” on page 144).
- Close all plate files and data files.

1. In the QuantStudio™ Design and Analysis Software v2, select **System** ▶ **Enable Security**.
2. Enter your SAE administrator account username and password, then click **Sign In**.


The SAE administrator account is automatically signed into the software after SAE is enabled. The SAE username is displayed in the upper-right corner of the software menu bar.

Disable SAE on the instrument (administrator only)

This procedure requires a local administrator profile and an SAE administrator account.

Sign in with a local administrator account (see “Sign in as a local administrator with SAE enabled” on page 146).

1. In the home screen, tap **⚙ (Settings)** ▶ **SAE**.
The **SAE Mode** screen is displayed.
2. In the **SAE Mode** screen, set the **SAE Mode** slider to the **Disable** position, then tap **Done**.
3. Enter the password for the SAE administrator account, then tap **Disable**.

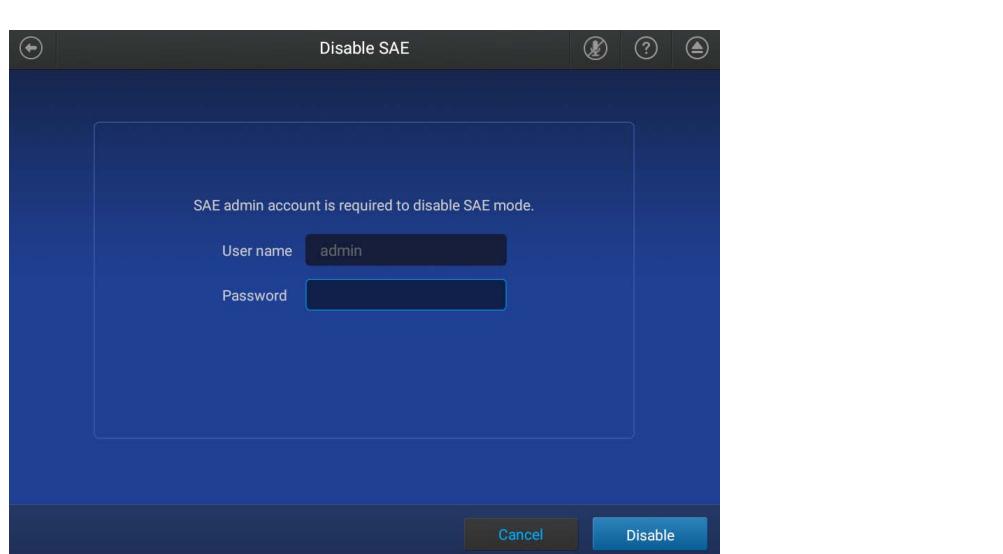


The **Sign In** screen is displayed.

Sign in as a local administrator with SAE enabled

Sign in as a local administrator to access the instrument settings. Plate files are not accessible if you are signed in as a local administrator.

1. In the **Sign In** screen, tap **Sign in** under **Local sign in**.

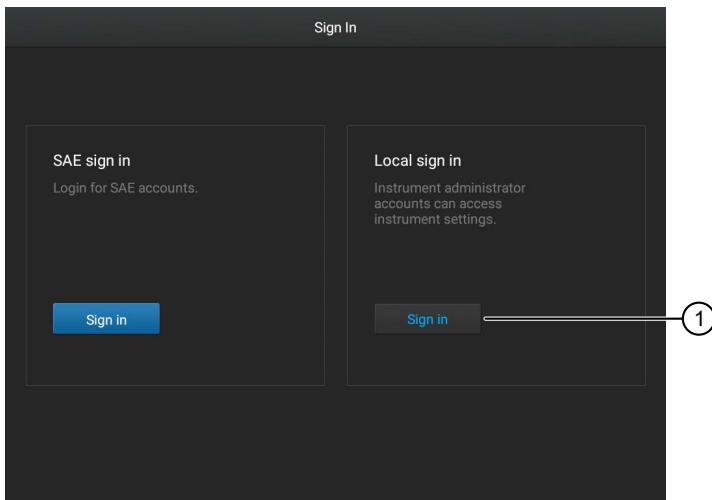

① Sign in for administrator

The **Local Administrator Sign In** screen is displayed.


2. In the **Local Administrator Sign In** screen, select your local administrator profile.
3. Enter your PIN, then tap **Enter**.

The **Settings** screen is displayed.

Sign in with SAE enabled

You must create a profile on the SAE Administrator Console before you sign in when SAE is enabled on the instrument.

1. In the **Sign In** screen, tap **Sign in** under **SAE sign in**.

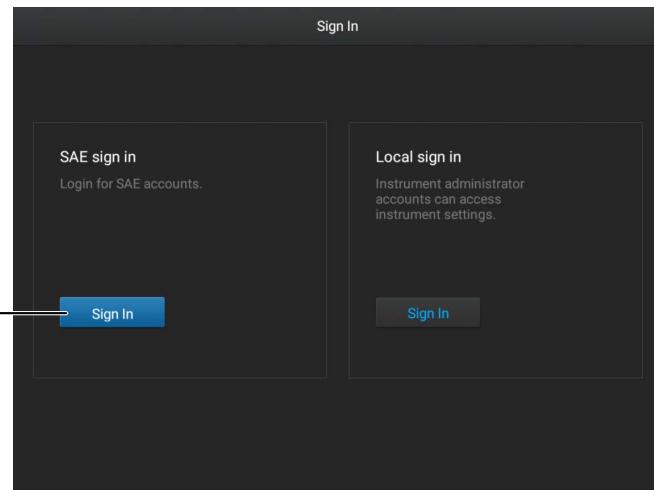

① Sign in to an SAE account

The **SAE Sign In** screen is displayed.

2. In the **SAE Sign In** screen, enter your SAE user name and your SAE password, then tap **Sign in**.

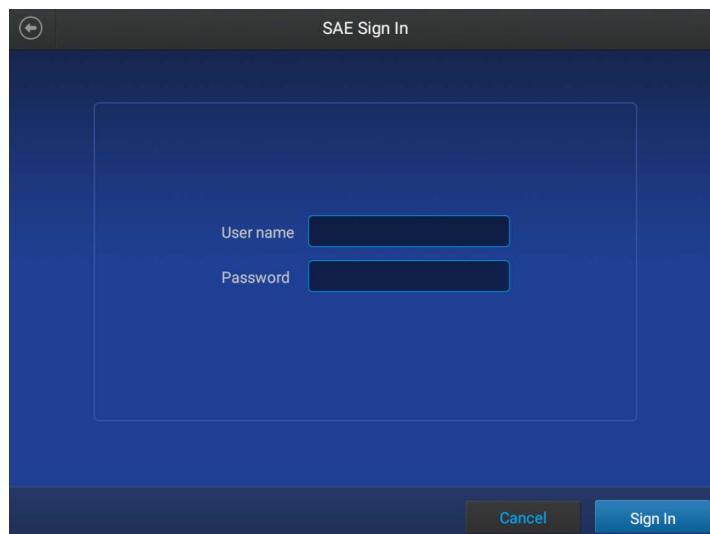


Figure 19 SAE Sign In screen

The home screen is displayed.

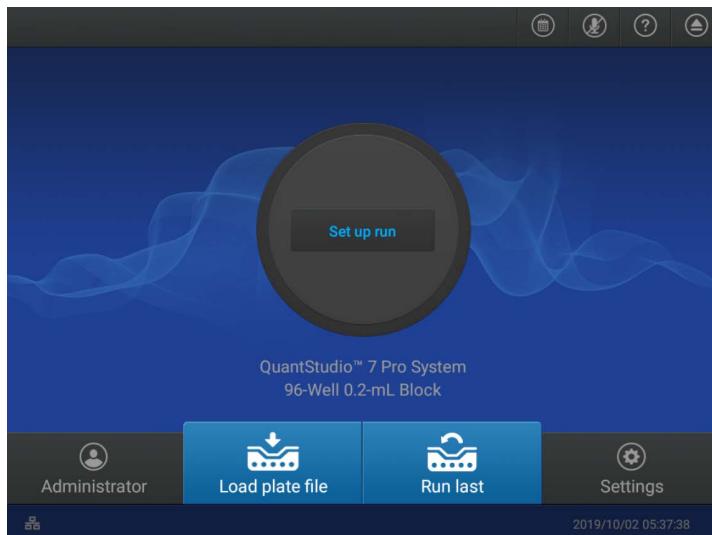


Figure 20 Home screen

Sign out with SAE enabled

1. In the home screen, tap  **(Profile)**.

The **My Profile** screen is displayed. The **SAE** button is selected and the profile that is signed in is listed.

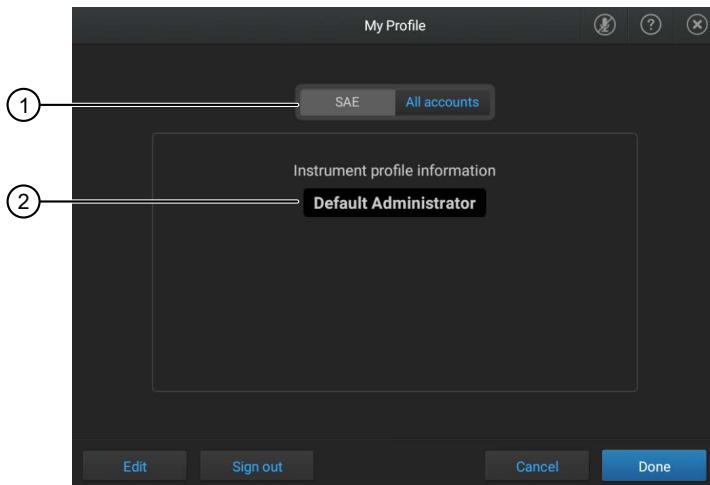


Figure 21 My Profile screen

- ① **SAE** button is active
- ② Profile that is signed in

2. Tap **Sign out**, then tap **Yes** to confirm.

You are signed out and the **Sign in** screen is displayed.

Change your SAE account password on the instrument

The QuantStudio™ Design and Analysis Software v2 and the instrument must be connected to the same SAE server in order for the SAE account password to be updated for both the software and the instrument.

You must be signed in with the instrument in SAE mode.

See “Sign in with SAE enabled” on page 147.

1. In the home screen, tap  **(Profile)**.

The **My Profile** screen is displayed. The **SAE** button is selected, and the profile that is signed in is listed.

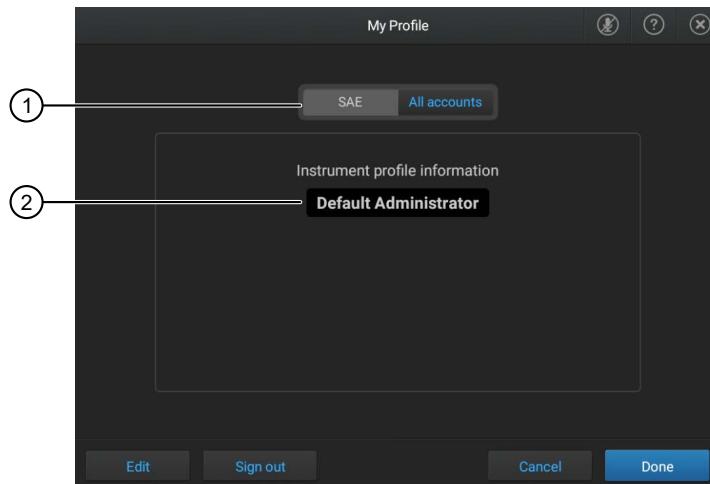


Figure 22 My Profile screen

- ① **SAE** button is active
- ② Profile that is signed in

2. Tap **Edit**.

The **Edit My Profile** screen is displayed.

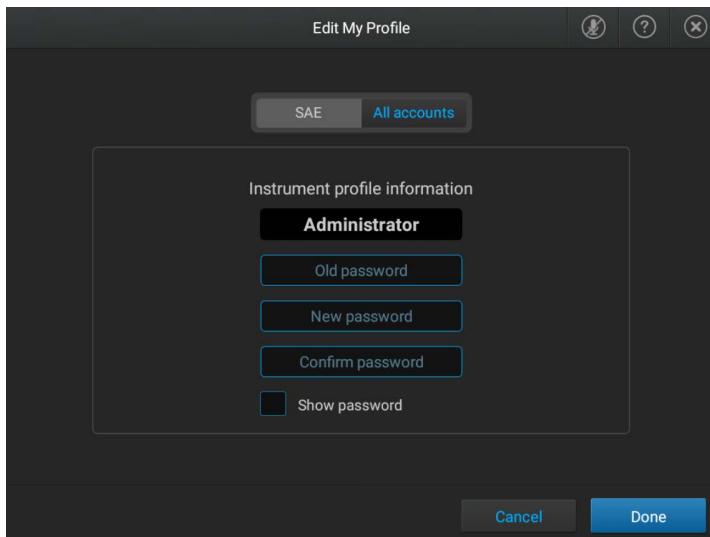


Figure 23 Edit My Profile screen

3. Tap the **Old password** field, enter the current SAE account password, then tap **Enter**.
4. Tap the **New password** field, enter a new SAE account password, then tap **Enter**.
(Optional) Tap the **Show password** checkbox to show or hide the password.
5. Tap the **Confirm password** field, enter the new SAE account password again, then tap **Enter**.
6. Tap **Done**.

Your SAE password will also be changed on the QuantStudio™ Design and Analysis Software v2 if the instrument and the software are connected to the same SAE server.

Sign a plate setup

1. In the home screen, tap one of the following options:
 -  **(Load plate file)**
 - **Set up run**
2. Tap the location of the system template or the plate file in the left column.
 -  **Run Queue**
 -  **Public**
 -  **My Instrument**
 -  **Network Drive**
 -  **USB Drive**
3. Tap the template file name in the right column.


The **Plate Properties** screen is displayed.

4. In the **Plate Properties** screen, tap **Actions**.

The **Actions** screen is displayed.

5. Tap **Sign plate**.

The **Sign Plate** screen is displayed.

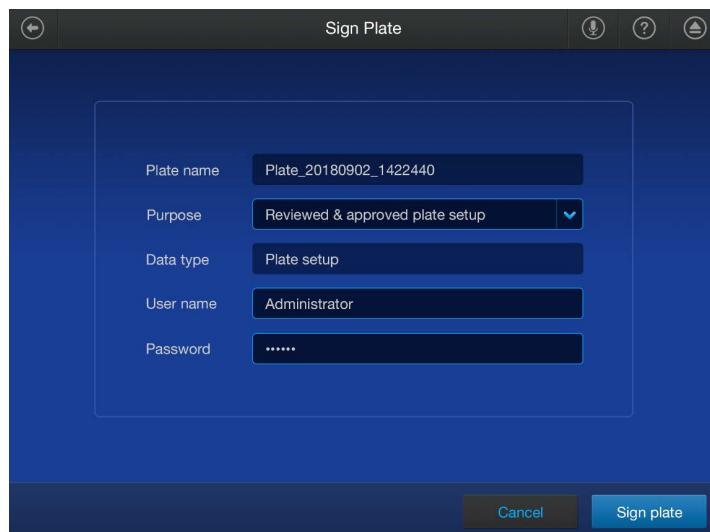

6. Select the purpose from the **Purpose** dropdown list.


7. Enter your SAE user name and password.

8. Tap **Sign plate**.

Enter an audit reason

Depending on the way that your SAE administrator configures audit settings, the **Enter Audit Reason** screen may be displayed when you make changes to one of the following items:

- Plate file
- Calibration
- RNase P run

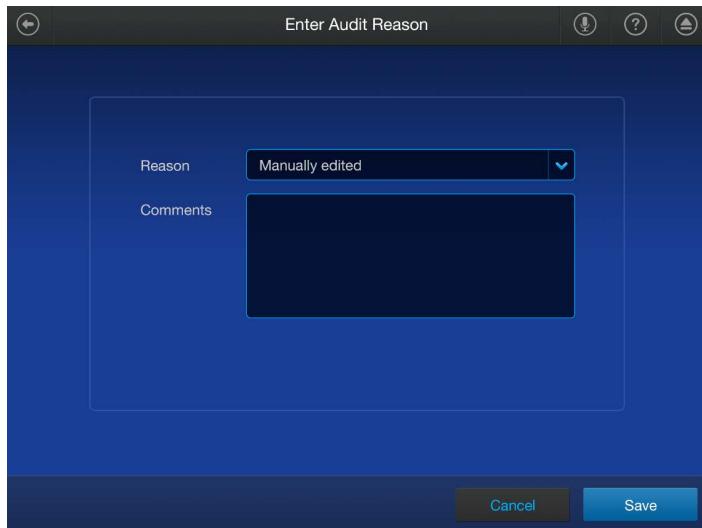


View signing records for a plate file

1. In the home screen, tap one of the following options:
 - **(Load plate file)**
 - **Set up run**
2. Tap the location of the system template or the plate file in the left column.
 - **Run Queue**
 - **Public**
 - **My Instrument**
 - **Network Drive**
 - **USB Drive**
3. Tap the template file name in the right column.
The **Plate Properties** screen is displayed.
4. In the **Plate Properties** screen, tap **Actions**.
The **Actions** screen is displayed.
5. Tap **View signing records**.
The **Signing records** screen is displayed.



6. Tap a row to view the details.

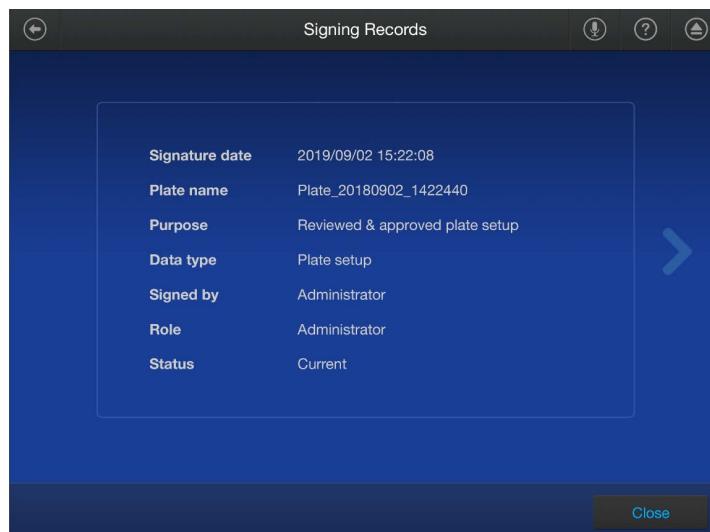

7. Tap **Close** to return to the **Signing records** screen.

View audit records

The audit records cannot be viewed on the instrument touchscreen. The audit records are viewed on the QuantStudio™ Design and Analysis Software v2 or on the SAE Administrator Console.

Use the instrument when the SAE server is offline

If your SAE administrator has configured your instrument to allow use when the SAE server is offline (**Client offline login** System setting in the SAE module), you can use the instrument for the period of time specified by the SAE administrator for **Client offline login**.

Note: If you have not previously signed in to the instrument with your SAE account, you cannot sign in when the SAE server is offline.


All SAE records are retained if the instrument is disconnected from an SAE server. When the instrument is reconnected to the SAE server, SAE records are uploaded to the server.

The following functions are not available when the SAE server is offline:

- Account lockout, password reminder, mandatory password change
- Disable SAE
- Change Password

E-signature requirements to start a run

Depending on the way your SAE administrator has configured your instrument, E-signatures may be required when you start a run. If starting a run is configured to require signatures, a message is displayed when you start a run.

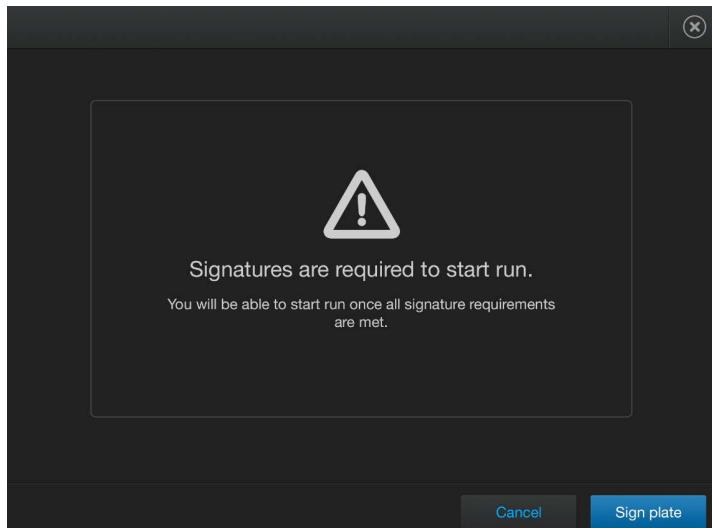

Tap **Sign plate** to sign the plate and start the run.

Signing in after automatic screen locking

Depending on the way your SAE administrator has configured your instrument, the instrument touchscreen may automatically lock after a specified duration.

An Administrator role or a Scientist role can sign in after automatic screen locking.

SAE error messages and actions

Message	Possible cause	Action
Unable to connect to SAE server. Check current connections.	The SAE server connection settings are incorrect.	<ol style="list-style-type: none">1. Check the SAE server IP address.2. In the instrument Sign In screen, sign in with a local administrator account.3. Set the correct IP address (see “Enable SAE on the instrument and specify the SAE server (administrator only)” on page 143).
	There is a problem with the computer on which the SAE Administrator Console is installed or a problem with the network.	Troubleshoot computer or network problems.
	The computer on which the SAE Administrator Console has a dynamic IP address that is disconnecting the server when the computer is restarted.	Set a static IP address on the computer.

■ Install a block	156
■ Change the block	157
■ Remove the block	158
■ Decontaminate the sample block	159
■ Replace the instrument fuses	161
■ Power on or off, store, and move	163

IMPORTANT! This chapter contains user maintenance procedures for the instrument. Procedures other than those described in this document must be performed by a qualified Thermo Fisher Scientific service representative.

Install a block

This section describes installing a block when there is no block in the instrument. To change a block, see “Change the block” on page 157.

A block can be installed by any user. This function is not restricted to administrators.

Each combination of a block and an instrument has its own calibration status. To determine the calibration status of a block on an instrument, see “View the calibration status and set reminders” on page 87. To calibrate a block on a specific instrument, see Chapter 7, “Calibrate and verify instrument performance”.

1. Tap  (Settings) ▶ Change block.
2. Insert the block and heated cover into the instrument drawer.
The instructions are displayed on the screen.
3. Tap **Install**.

The instrument displays one of the following options, depending on the calibration status:

- Confirmation that the block was installed successfully and the expiration date of the current calibration.
- A warning that the block is not calibrated on the instrument.
- A warning that the calibration of the block on the instrument is expired.

Change the block



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the sample block temperature can reach 100°C. Allow it to cool to room temperature before handling.

This section describes changing a block and heated cover when there is already a block installed. To install a block when one is not installed, see “Install a block” on page 156.

A block can be changed by any user. This function is not restricted to administrators.

Each combination of a block and an instrument has its own calibration status. To determine the calibration status of a block on an instrument, see “View the calibration status and set reminders” on page 87. To calibrate a block on a specific instrument, see Chapter 7, “Calibrate and verify instrument performance”.

1. Tap  **(Settings)** ▶ **Change block**.

2. In the **Change Block** screen, tap **Change**, then tap **Ok**.


The instrument unlocks the block and the heated cover. This step takes approximately one minute. The drawer opens with the block and the heated cover.

3. Use the handles to remove the block and the heated cover.

The instructions are displayed on the screen.

The block and the heated cover are removed together.

IMPORTANT! Only touch the block and the heated cover at the handles.

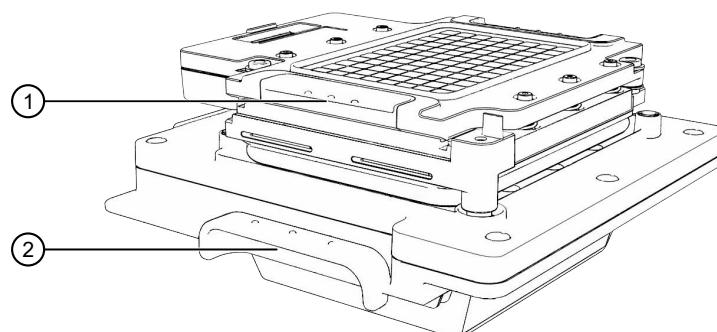

① Heated cover handle

② Block handle

4. Insert the new block and the heated cover into the instrument drawer, then tap **Install**.

The **Install** button is inactive until the instrument detects a block and heated cover that are positioned correctly in the drawer.

The instrument displays one of the following options, depending on the calibration status:

- Confirmation that the block was installed successfully and the expiration date of the current calibration.
- A warning that the block is not calibrated on the instrument.
- A warning that the calibration of the block on the instrument is expired.

It is recommended to store the block and the heated cover in the provided protective box when they are not installed on an instrument.

Remove the block



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the sample block temperature can reach 100°C. Allow it to cool to room temperature before handling.

This section describes removing the block and heated cover without installing another block and heated cover into the instrument.

1. Tap  (Settings) ▶ Change block.
2. In the **Change Block** screen, tap **Change**.


3. Tap **OK**.

The instrument unlocks the block and the heated cover. This step takes approximately one minute. The drawer opens with the block and the heated cover.

4. Use the handles to remove the block and the heated cover.

The block and the heated cover are removed together.

IMPORTANT! Only touch the block and the heated cover at the handle.

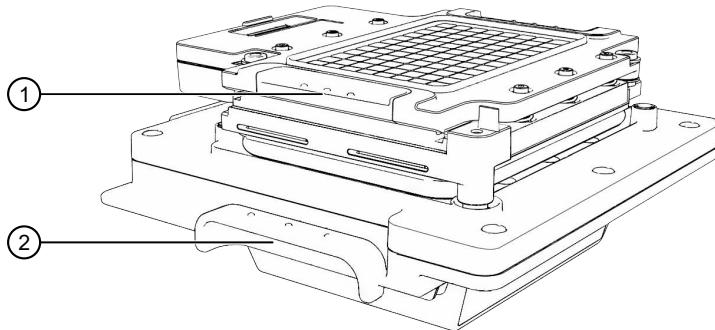

- (1) Heated cover handle
- (2) Block handle

5. Tap  (Eject) to close the instrument drawer.

It is recommended to store the block and the heated cover in the provided protective box when they are not installed on an instrument.

Decontaminate the sample block

Perform this procedure to eliminate fluorescent contaminants from the instrument sample block. Contamination is generally evident in failed background calibrations where one or more wells consistently exhibit abnormally high signals.



CAUTION! PHYSICAL INJURY HAZARD. Do not remove the instrument cover. There are no components inside the instrument that you can safely service yourself. If you suspect a problem, contact Support.



CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the sample block temperature can reach 100°C. Allow it to cool to room temperature before handling.



CAUTION! Before using a cleaning or decontamination method other than those recommended by Thermo Fisher Scientific, confirm with Thermo Fisher Scientific that the proposed method will not damage the equipment.

Materials required

- Safety glasses
- Powder-free gloves
- Tissue, lint-free
- Cotton or nylon swabs and lint-free cloths
- Pipette (100- μ L) with pipette tips
- Deionized water
- Ethanol, 95% solution
- Bleach, 10% solution

Clean the sample block




CAUTION! PHYSICAL INJURY HAZARD. During instrument operation, the sample block temperature can reach 100°C. Allow it to cool to room temperature before handling.

IMPORTANT! Wear powder-free gloves when you perform this procedure.

IMPORTANT! Always use deionized water to rinse wells after cleaning with bleach or ethanol solution.

1. Identify the contaminated wells of the sample block (see “Identify contamination” on page 95).
2. Remove the sample block and heated cover (see “Remove the block” on page 158).
3. Place the sample block and heated cover on a clean surface.

4. Rinse the contaminated wells with deionized water (see “Detailed procedures for cleaning the sample block” on page 160).

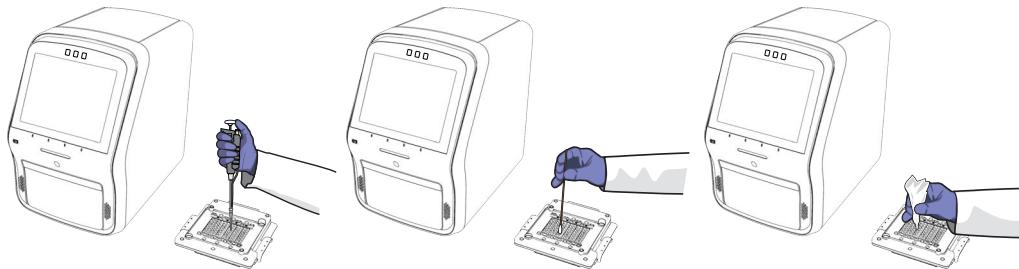

5. Install the block and heated cover (see “Install a block” on page 156).
6. Perform a background calibration to confirm that you have eliminated the contamination (see “Perform calibrations” on page 91).
7. If the contamination remains, clean the contaminated wells with a 95% ethanol solution.

Repeat step 2 to step 6 with a 95% ethanol solution (see “Detailed procedures for cleaning the sample block” on page 160).

IMPORTANT! Always use deionized water to rinse wells after cleaning with bleach or ethanol solution.

8. If the contamination remains, clean the contaminated wells with a 10% bleach solution.

Repeat step 2 to step 6 with a 10% bleach solution (see “Detailed procedures for cleaning the sample block” on page 160).

IMPORTANT! Always use deionized water to rinse wells after cleaning with bleach or ethanol solution.

9. Use Smart Help if the contamination remains (see “Request technical support with Smart Help” on page 172).

Detailed procedures for cleaning the sample block

IMPORTANT! Use these cleaning procedures *only* in conjunction with the complete decontamination procedure (see “Clean the sample block” on page 159).

- Rinse the sample block with deionized water.
 - Pipet a small volume of deionized water into each contaminated well.
 - In each well, pipet the water up and down several times to rinse the well.
 - Pipet the water to a waste beaker.
 - Use a cotton swab to scrub inside of each contaminated well.
 - Use a lint-free cloth to absorb the excess deionized water.

- Clean the sample block with 95% ethanol.
 - a. Pipet a small volume of 95% ethanol solution into each contaminated well.
 - b. In each well, pipet the solution up and down several times to rinse the well.
 - c. Pipet the ethanol solution to a waste beaker.

IMPORTANT! Always use deionized water to rinse wells after cleaning with bleach or ethanol solution.

- Clean the sample block with 10% bleach.
 - a. Pipet a small volume of 10% bleach solution into each contaminated well.
 - b. In each well, pipet the solution up and down several times to rinse the well.
 - c. Pipet the bleach solution to a waste beaker.

IMPORTANT! Always use deionized water to rinse wells after cleaning with bleach or ethanol solution.

Replace the instrument fuses



CAUTION! FIRE HAZARD. For continued protection against the risk of fire, replace fuses only with listed and certified fuses of the same type and rating as those currently in the instrument.


Materials required

- Fuses (2) – 10A, Time-Lag T, 250VAC, 5 × 20 mm
- Safety glasses
- Powder-free gloves
- Screwdriver, flathead

Replace the fuses

1. Power off, then unplug the instrument.
2. Allow the instrument cool for 15 minutes.

3. Using a flat-head screwdriver, unscrew and remove the fuse holder.

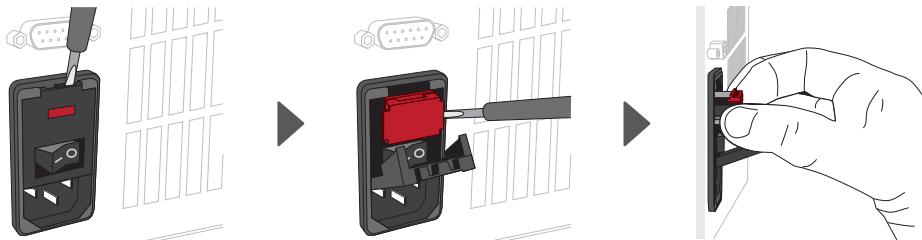

4. Remove each fuse from its fuse holder, then inspect it for damage.


Carbon typically coats the inside of failed fuses.

Good	Failed

5. Replace each failed fuse.

Note: The voltage and amperage ratings are on the fuse holder.

6. Align each new fuse in the fuse holder as shown in the figure below.

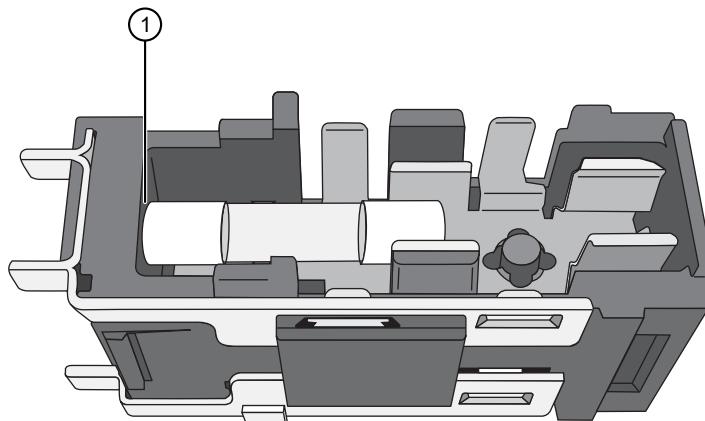

① Fuse alignment in the fuse holder

7. Install the fuse holder back into the instrument.

8. Plug in, then power on the instrument.

The installation is successful if the instrument powers on.

Note: Fuse failure can result from fluctuations in the supplied power to the system. To prevent further failures, consider installing an electrical protective device, such as a UPS or a surge protector. If issues with the fuse persist, use Smart Help (see “Request technical support with Smart Help” on page 172).

Power on or off, store, and move

Enable and edit sleep mode

Sleep mode allows the instrument to use less energy when not in use. The length of time that the instrument is inactive before going into sleep mode can be customized.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Sleep mode**.
2. Set the slider.

Option	Description
Off	The instrument will not go into sleep mode.
On	The instrument will go into sleep mode after a period of inactivity.

3. Tap the **Idle time before sleep mode** field, then enter a time.
4. Tap **Done**.

Power on the instrument

If left unattended, the instrument automatically enters sleep mode to conserve power. The default time that an instrument is inactive before going into sleep mode is 15 minutes.

1. Tap anywhere on the touchscreen to determine if the instrument is in sleep mode.
If the screen becomes active, the instrument is already powered on.
2. If the home screen does not display, power on the instrument by pressing the ON/OFF switch on the rear panel.

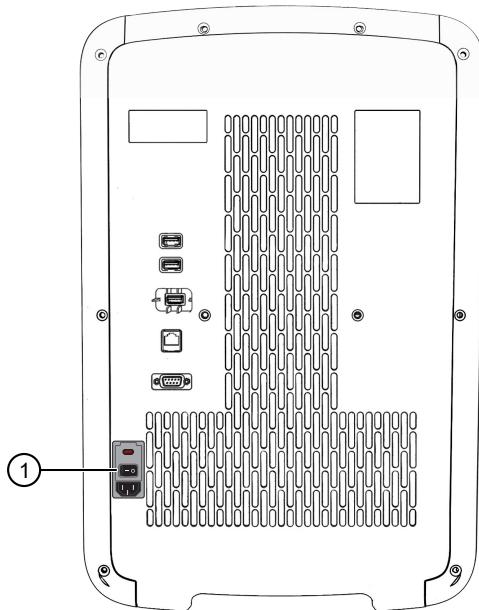


Figure 24 Location of the ON/OFF switch

① ON/OFF switch

Note: The length of time that the instrument is inactive before going into sleep mode can be customized (see “Enable and edit sleep mode” on page 163).

Power off the instrument

1. Power off the instrument using the ON/OFF switch on the back of the instrument.
For the location of the ON/OFF switch, see Figure 6 on page 43.
2. Power off the computer if there is a computer directly connected to the instrument.

Prepare the instrument to store, move, or ship

Remove the block and heated cover to store, move, or ship the instrument.

It is recommended to store the block and heated cover in the provided protective box when they are not installed on an instrument.

Ship the instrument and the block and heated cover separately.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Ship prep mode** ▶ **Next**.
The **Ship Prep Mode** screen is displayed.
2. In the **Ship Prep Mode** screen, tap **Next**.
3. Follow the instructions on the touchscreen to remove the block and heated cover.

The instrument is now ready to store, move, or ship.

Move the instrument



CAUTION! PHYSICAL INJURY HAZARD. Do not attempt to lift the instrument or any other heavy objects unless you have received related training. Incorrect lifting can cause painful and sometimes permanent back injury. Use proper lifting techniques when lifting or moving the instrument. At least two people are required to lift it.

IMPORTANT! Moving your instrument can create subtle changes in the alignment of the instrument optics. Recalibrate the instrument if necessary.

- Ensure that the surface on which you place the instrument can support at least 35.85 kg (79 lbs.).
- Ensure that the path to transport the instrument is clear of obstructions.
- At least two people are needed to lift and carry the instrument.
- Keep your spine in a good neutral position.
- Bend at the knees and lift with your legs.
- Do not lift an object and twist your torso at the same time.
- Coordinate your intentions with your assistant before lifting and carrying.

IMPORTANT! After moving the instrument, perform an RNase P instrument verification run. If the run fails, perform ROI/ uniformity, background, and dye calibrations.

■ View the instrument details	166
■ View the license agreement	167
■ View the privacy policy for the Connect cloud-based platform	167
■ Configure the instrument settings	168
■ Configure maintenance and service settings	172
■ Request remote support	183
■ Configure the instrument for the hands-free features	185
■ Manage the run history	185
■ Manage the file name convention of the data file	185
■ Enable or disable the reminder to insert a plate	186
■ Select the language	186
■ View the notifications	187
■ Enable and disable home screen notifications	187
■ Set up email notifications	187
■ Clear the notifications	188

View the instrument details

1. In the home screen, tap  **(Settings)** ▶ **About instrument**.

The **About Instrument** screen displays the following information:

• Model name	• Block serial number
• Wired IP address	• Heated cover serial number
• Wireless IP address	• Block type
• Instrument serial number	• Software version

2. *(Optional)* Tap **Details**.

The following information is displayed:

- Instrument model name
- Wired IP address
- Wired MAC address
- Wireless IP address
- Wireless MAC address
- Instrument serial number
- Block serial number
- Heated cover serial number
- Block type
- Software version
- Firmware version
- User interface version
- Instrument server version
- Heated cover firmware version
- Block firmware version


3. Tap  **(Back)** to return to the **Settings** screen.

View the license agreement

1. In the home screen, tap  **(Settings)** ▶ **About instrument**.
The **About Instrument** screen is displayed.
2. Tap **EULA**.
The end user license agreement (EULA) is displayed.
3. *(Optional)* Tap **Export**.
4. Tap **Close**.

View the privacy policy for the Connect cloud-based platform

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Connect services**.
The **Connect Services** screen is displayed.

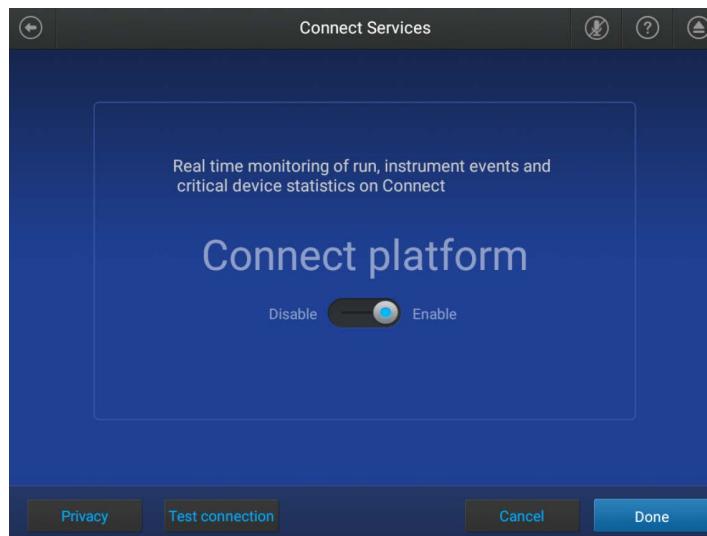

2. Tap **Privacy**.

The privacy policy is displayed.

3. Tap **Close**.

Configure the instrument settings

Enable or disable demonstration mode

Note: The instrument restarts when demonstration mode is turned on or off.

When the instrument is in demonstration mode, the following features apply:

- The instrument simulates a run but does not start a run.
- The instrument drawer will not open.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Demo mode**.

Turn demo mode off is displayed if the instrument is in demo mode.

2. In the **Confirmation** screen, tap **Ok**.

You will be prompted to restart the instrument.

Demo mode is displayed on the instrument when it is in demonstration mode.

Edit the instrument name (administrator only)

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Instrument name**.2. Tap the **Instrument name** field, then enter or edit the instrument name.

The instrument name must be unique if it is connected to a network.

3. Tap **Done**.

Set the date and the time

The date and time can be set manually or automatically. The instrument must be connected to the internet to set the date and time automatically.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Date and time**.

The **Date and Time** screen is displayed.

2. Set the slider to **Off** or **On** to determine if the date and time are set automatically.3. *(Optional)* Select the time zone from the dropdown list.

4. Select the date format and the time format from the dropdown lists.

5. Tap **Done**.

Configure the heated cover settings

The heated cover can be heated when the instrument is idle or it can be left without being heated when the instrument is idle. To save power, configure the settings so the heated cover is left without heating. It will take longer to bring the heated cover up to the correct temperature when the instrument is brought back into use.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Cover temperature**.
The **Heated Cover** screen is displayed.
2. Set the slider.

Option	Description
On	The heated cover remains heated when the instrument is idle.
Off	The heated cover is not heated when the instrument is idle.

3. Tap the **Cover temperature** field to edit the heated cover temperature.
4. Tap **Done**.

Adjust the speaker volume

The selected speaker volume is saved for each user.

When a new Connect profile is linked to the instrument, the default speaker volume is set to the maximum.

1. In the home screen tap,  **(Settings)** ▶ **Instrument settings** ▶ **Speaker volume**.
The **Volume Control** screen is displayed.
2. Slide the control to the adjust the volume.
The speakers are muted when the slider is at the far left side of the bar.
3. Tap **Done**.

Configure the network

For detailed instructions, see Appendix B, “Connect the instrument to a network”.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Network configuration**.
The **Network Configuration** screen is displayed.
2. Tap **Edit**.
The fields for **Wired** and **Wireless** are now editable.
3. Enter the information that is needed for a wireless or wired connection.
 - Tap the **Network** field to enter network information for a wireless connection.
 - Tap the **IP address** field to enter the IP address for a wired connection.

4. (Optional) Tap **Proxy server** to connect the instrument to a proxy server (see “Connect a proxy server” on page 170).
5. Tap **Done**.

Connect a proxy server

A proxy server sits between the instrument and the main server. The proxy server intercepts requests to the main server and evaluates if it can complete the request. If it cannot fulfill the request, it will forward the request to the main server.

Consult your Network or IT specialist before linking the instrument to a proxy server.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Proxy server**.
The **Network Configuration** screen is displayed.
2. Tap each field to enter the required information.
 - **Proxy server**
 - **Proxy port**
 - **User name**
 - **Password**
3. Tap **Done**.

The instrument is connected to the specified proxy server.

Select the region for the Connect platform

Note: The instrument restarts when the region for the Connect platform is changed.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Connect region**.
The **Select Region** window opens.
2. Select the region.

Option	Description
China	For users in China
U.S.	For users in any country other than China

3. Tap **Done**.

Select a disk storage management mode

This setting determines how the instrument manages data files when the disk space is full.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Disk storage management**.
2. Select a mode for disk storage management.

Option	Description
Automatic	<p>The instrument automatically deletes data files until there is disk space available to complete a run.</p> <p>Note: The instrument deletes data files in chronological order, starting with the oldest files. Only data files that have been transferred to another location are deleted.</p> <p>IMPORTANT! The system does not notify a user when data files are deleted.</p>
Manual	<p>The instrument will not delete data files when there is no disk space available to complete a run.</p> <p>Note: The system notifies a user when data files must be deleted (see “Delete data files from the instrument” on page 82).</p>

3. Tap **Done**.

Enable and edit automatic sign-out

The automatic sign-out feature signs out a user after a period of inactivity.

A run in progress is not considered inactivity. The period of inactivity begins after the run is complete.

The period of inactivity before a user is signed out automatically can be customized.

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Auto sign-out**.
2. Set the slider.

Option	Description
Off	The instrument will not automatically sign out a user after a period of inactivity.
On	The instrument will automatically sign out a user after a period of inactivity.

3. Tap the **Idle time before auto sign-out** field, then enter a time.
4. Tap **Done**.

Set an access key for external software to connect to the instrument

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Instrument access**.
2. In the **Instrument Access** screen, set the **Enable access** slider to the **On** position.
The instrument displays an access code.

3. *(Optional)* Set the **Visible to all users** slider.

The **Visible to all users** slider is only displayed if the **Enable access** slider is set to the **On** position.

4. Tap **Close**.

5. In the external software that is being used to connect to the instrument, enter the remote access key to add the instrument.

Reset the access key for external software to connect to the instrument

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Instrument access**.

2. Tap **Generate code**.

The instrument displays a new access code.

3. *(Optional)* Set the **Visible to all users** slider.

4. Tap **Close**.

In the external software that is being used to connect to the instrument, enter the new remote access key to add the instrument.

Configure maintenance and service settings

Request technical support with Smart Help

Request technical support with Smart Help for any technical problem with the instrument. A Connect profile is required.

Note: A recording to describe the technical problem is not stored on the instrument.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Smart Help**.

The **Smart Help** screen is displayed.

2. Tap **Edit**, then enter a description in the **Description** field.

3. Fill out the **Email** field and the **Phone** field.

4. Select a country from the dropdown list.

5. Select a preferred contact method from the dropdown list.

6. *(Optional)* Select the **Include instrument log file** checkbox.

The log file is not required but it is recommended for troubleshooting.

7. *(Optional)* Tap **Attach run files**, select run files, then tap **Attach files**.

8. *(Optional)* Tap  to record a description of the technical issue.
9. Tap **Done**.

A confirmation email is sent. It is sent even if email was not selected as the preferred contact method.

Request instrument service with Smart Help

Request instrument service with Smart Help for any instrument service request. A Connect profile is required.

Note: A recording to describe the technical problem is not stored on the instrument.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Smart Help**.
The **Smart Help** screen is displayed.
2. Tap **Edit**, then enter a description in the **Description** field.
3. Fill out the **Email** field and the **Phone** field.
4. Select a country from the dropdown list.
5. Select a preferred contact method from the dropdown list.
6. *(Optional)* Select the **Include instrument log file** checkbox.
The log file is not required but it is recommended for troubleshooting.
7. *(Optional)* Tap **Attach run files**, select run files, then tap **Attach files**.
8. *(Optional)* Tap  to record a description of the technical issue.
9. Tap **Done**.

A confirmation email is sent. It is sent even if email was not selected as the preferred contact method.

Perform and view self-verification tests

Perform a hardware test

There will be elevated noise levels while the fans are tested.

IMPORTANT! Ensure that there is no obstruction at the front of the instrument. The instrument drawer opens and closes as part of the test.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Self verification test**.
The **Self Verification Test** screen is displayed.
2. Tap **Hardware test**.
3. Tap **Start test**.
The test takes approximately 10 minutes.

The **Hardware Test** screen is displayed when the self-verification test is complete. The **Hardware Test** screen displays the functions that pass and fail.

4. *(Optional)* Tap **Export** to export the results.
5. Tap **Close**.

Use Smart Help if any hardware functions fail (see “Request technical support with Smart Help” on page 172).

View the last hardware test result

The self verification test checks the instrument hardware functions. You can view the results of the last self-verification test that was run.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Self verification test**.
The **Self-Verification Test** screen is displayed.
2. Tap **Hardware test**.
3. Tap **Last test**.
The **Self-Verification Test** screen displays the functions that pass and fail.

Use Smart Help if any hardware functions fail (see “Request technical support with Smart Help” on page 172).

Perform a peripheral test


The peripheral test takes approximately 5 minutes. Tap  **(Close)** on any screen to exit the peripheral test.

The peripheral test checks the following functions:

- Touchscreen accuracy
- Speaker
- Microphone
- Camera
- Proximity sensor


1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Self verification test**.
The **Self Verification Test** screen is displayed.
2. Tap **Peripheral test**.
3. Tap **Start**.
The peripheral test begins with the touch accuracy test.

4. In the **Touch Accuracy Test** screen, tap the screen in several locations.

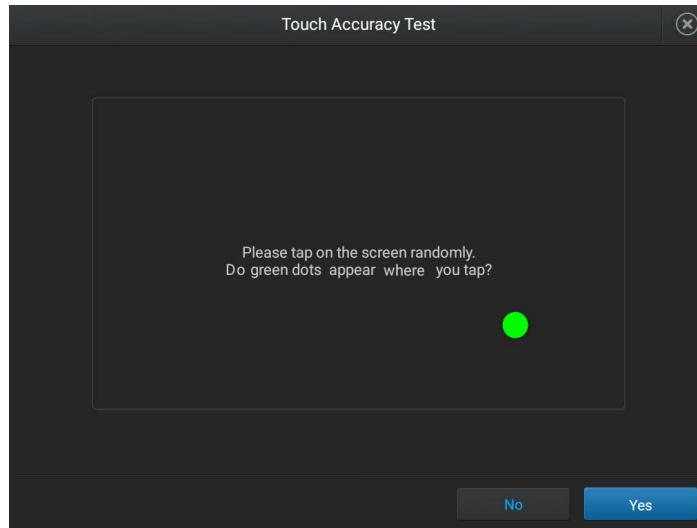

Green dots appear where you tap.


5. Tap **Yes** to proceed.
6. In the **Speaker Test** screen, tap **(Play)**.

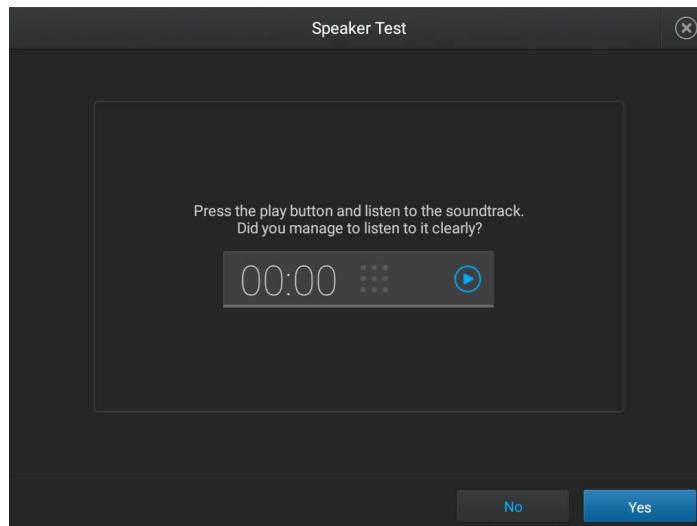


A soundtrack is played.

7. Tap **Yes** to proceed.


8. In the **Microphone Test** screen, tap  to record.

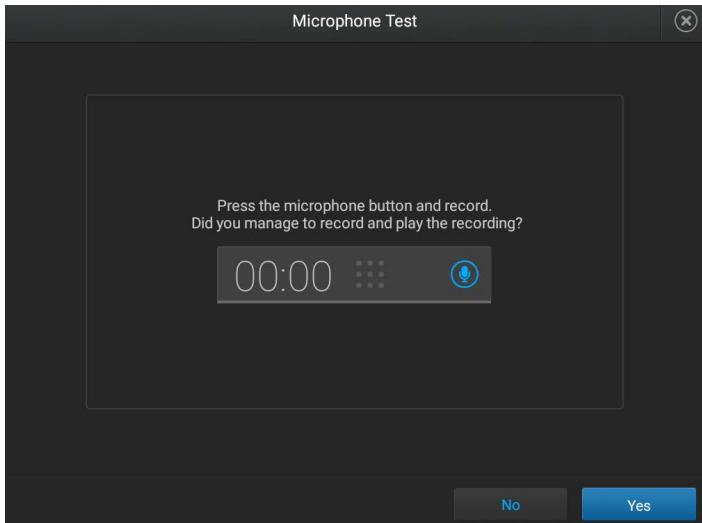


9. Tap **Yes** to proceed.
10. In the **Camera Test** screen, tap **Yes** to confirm that the camera is on.
A preview is displayed if the camera is on.
11. In the **Camera Test** screen, tap **Yes** to confirm that the camera detects a face.

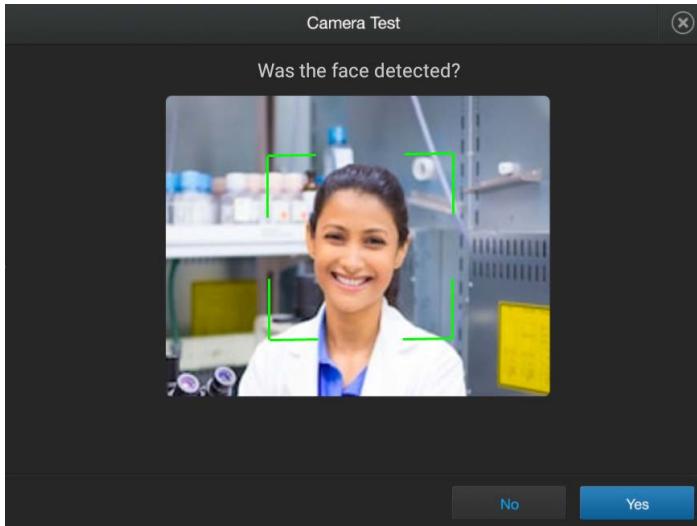


12. In the **Camera Test** screen, tap **Capture** to capture an image, then tap **Yes** to confirm that an image was captured.
The image should appear clear and in focus.

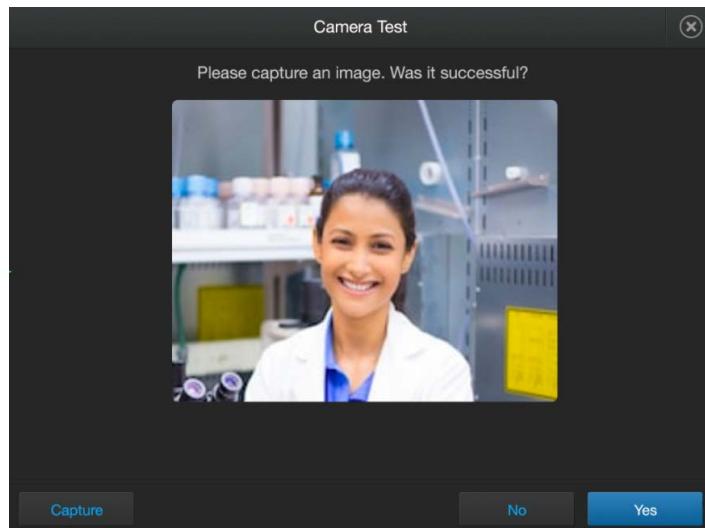

13. In the **Camera Test** screen, tap **Yes** to confirm that the depth map is displayed on the screen.

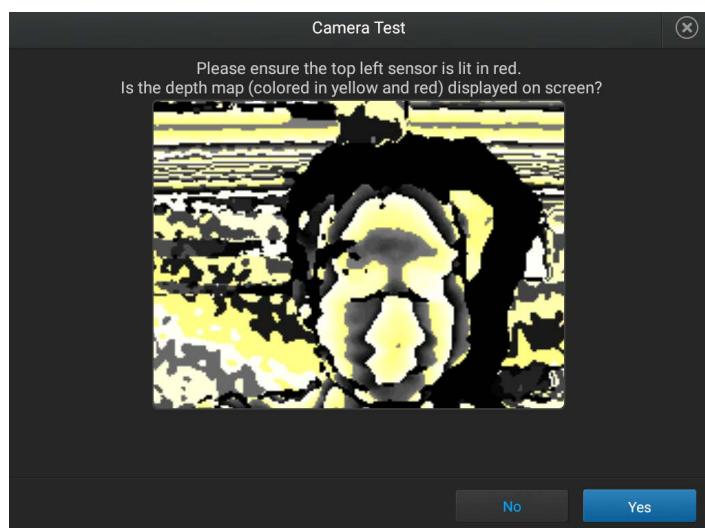


14. In the **Proximity Test** screen, move forwards towards the instrument and backwards away from the instrument.

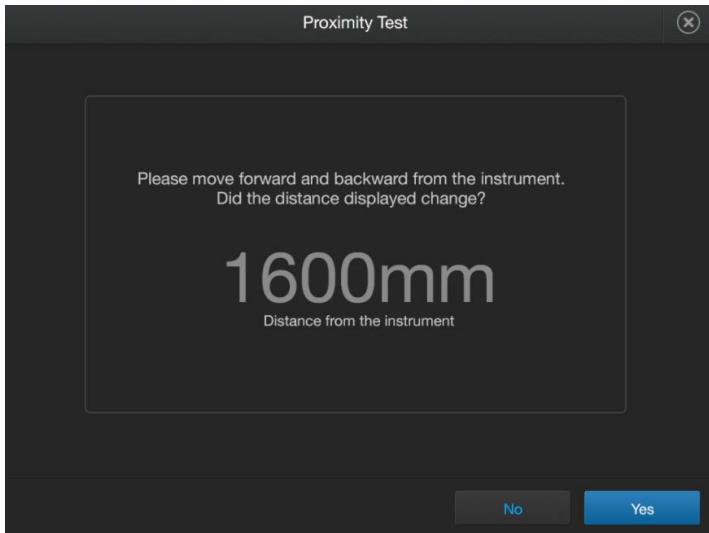

The distance on the instrument screen changes to show your distance from the instrument.

15. Tap **Yes** to complete the test.

The **Peripheral Test** screen is displayed. It shows a **Pass** or **Fail** result for each test.

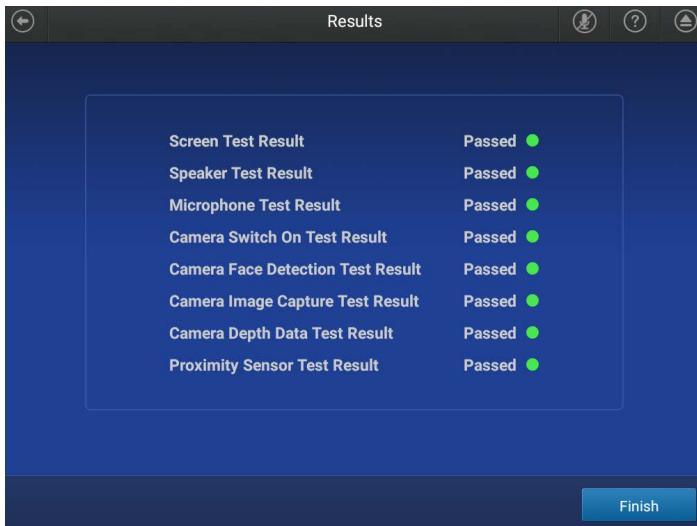

16. (Optional) Tap **Export** to export the results.

View the instrument log

1. In the home screen, tap **(Settings)** ▶ **Maintenance and service** ▶ **Instrument log**.
The **Instrument Log** screen lists the instrument event types with a time stamp.

- Run
- Change Block
- Error
- Power
- Calibration
- Self-Test
- Update
- User

2. *(Optional)* Tap any of the table headings to sort by that parameter.
3. *(Optional)* Tap an instrument event type in the left pane to filter the list that is displayed by that event type.
4. Tap **Cancel**.

Export the instrument log

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Instrument log**.
2. *(Optional)* Tap a type of instrument event to export the log for only that type of event.
If an instrument event is not selected, the logs for all the instrument events will be exported.
3. Tap **Export**.
4. Select a destination to save the instrument log files.
 -  **Connect**
 -  **Network Drive**
 -  **USB Drive**

Note: You must be signed in with a Connect profile in order to export the instrument log to Connect.

5. Tap **Done**.

View the instrument statistics

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Instrument statistics**.
The **Instrument Statistics** screen is displayed.
 - **Disk Space Remaining (MB)**
 - **Block Cycle Count**
 - **Block Degree Climbed**
 - **LED Usage (hours)**
 - **RNase P Status**
 - **Block Serial Number**
 - **Heated Cover Serial Number**
 - **Block Type**
2. Tap **Cancel**.

Enable or disable remote monitoring

The following items can be monitored remotely:

- The progress of a run
- Instrument events
- Instrument statistics

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Connect services**.
2. Set the slider to the **On** or **Off** position.
3. Tap **Done**.

Update the software

Software updates are displayed on the **Notifications** button in the **Settings** screen when they are available.



① Notifications, including software updates

To enable notifications to be displayed on the home screen, see “Enable and disable home screen notifications” on page 187.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Software update**.
The **Software Update** screen is displayed. It indicates whether a software update is available.
2. Tap **Update now**.
The software will update.

The **Software Update** screen will display **Update is complete** when the software update is complete. The instrument will restart.

The software can be repaired if the update failed (see “Repair the software” on page 181).


Repair the software

If a software update failed, the software can be repaired.

The software version installed during a repair is the same version that was previously installed on the instrument, even if a newer version is available.

Instrument settings, calibrations, profiles, and data files are preserved during a repair.

If the software update failed, the **Software Update** screen displays **Failed to upgrade**.

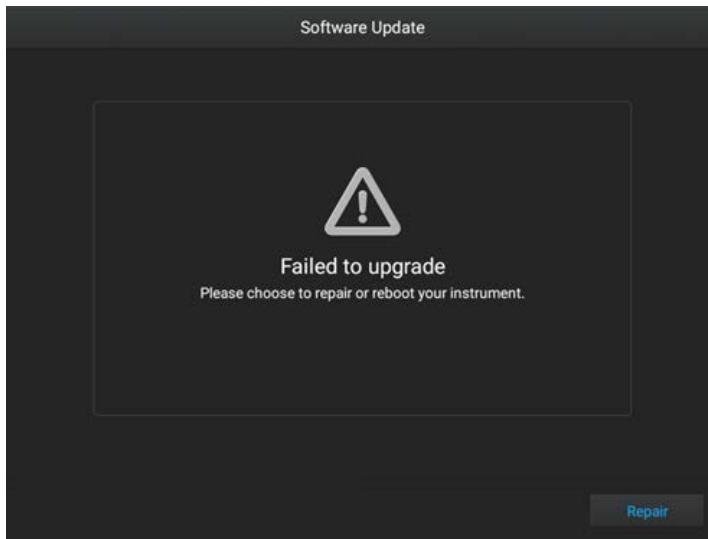

In the **Software Update** screen, tap **Repair**.

Back up the instrument

The instrument can be backed up to a USB drive or to Connect.

A back-up of the instrument can only be performed by an administrator.

IMPORTANT! A back-up of the instrument includes instrument profiles and settings. It does not include plate files or data files. Plate files and data files must be transferred separately. See Chapter 6, “View and manage files”.

1. In the home screen, tap  **(Settings)** ▶ **Maintenance and service** ▶ **Back up/Restore** ▶ **Back up instrument**.
The **Back Up/Restore** screen is displayed.
2. Select a location to store the backup of the instrument.
 -  **USB Drive**
 -  **Connect**
3. *(Optional)* If you are backing up the instrument to a USB drive, insert a USB drive into the USB port on the front of the instrument.

4. Enter a backup file name.
5. Select the elements to back up.
 - Instrument settings
 - User profiles
 - Calibration records
6. Tap **Back up**.
Instrument back up is complete is displayed on the **Back Up Instrument** screen when the back up is complete.
7. Tap **Done**.

Restore a backup of the instrument

IMPORTANT! A back-up of the instrument includes instrument profiles and settings. It does not include plate files or data files. Plate files and data files must be transferred separately. See Chapter 6, “View and manage files”.

Restoring a backup of the instrument can only be performed by an administrator.

System templates are included when a backup of the instrument is restored.

1. In the home screen, tap **⚙ (Settings)** ▶ **Maintenance and service** ▶ **Back up/Restore** ▶ **Restore a backup**.
The **Restore a Backup** screen is displayed.
2. Select the backup to restore.
3. Tap **Restore**.
Your instrument has been restored is displayed on the **Restore a Backup** screen.
4. Tap **Done**.

Upload a certificate for a TaqMan™ Array Plate with RFID

1. In the home screen, tap **⚙ (Settings)** ▶ **Maintenance and service** ▶ **RFID certificate**.
The TaqMan™ Array Plate with RFID **Certificate** screen is displayed.
2. Tap **Upload**.
3. Tap the location of the certificate.
 -  **Connect**

Note: You must be signed in with a Connect profile.

 -  **Network**
 -  **USB**



Overview of an instrument reset (restore factory default)

IMPORTANT! Back up the instrument before resetting the instrument (see “Back up the instrument” on page 181). Transfer plate files and data files (see Chapter 6, “View and manage files”). Plate files and data files are not included in a backup of the instrument.

Resetting the instrument deletes the following information:

- Instrument profiles
- User-created plate files and data files
- Custom dyes
- Custom melt calibrations

The following information is retained if the instrument is reset:

- System templates
- The last known calibrations, including expired calibrations

Reset the instrument (restore factory default)

IMPORTANT! Back up the instrument before resetting the instrument (see “Back up the instrument” on page 181). Plate files and data files are not backed up. Transfer plate files and data files (see Chapter 6, “View and manage files”).

For a description of the information that is deleted and the information that is retained during an instrument reset, see “Overview of an instrument reset (restore factory default)” on page 183.

1. In the home screen, tap **(Settings) ▶ Maintenance and service ▶ Restore factory default.**
2. Restart the instrument when prompted.

Request remote support

Remote support grants a support representative the ability to access the instrument remotely.


A Support representative can perform the following functions during a remote support session:

- Use a pointer that displays on the touchscreen
- Draw on the touchscreen
- Capture an image of the touchscreen
- Capture a recording of the touchscreen
- Control the instrument via the touchscreen
- Send keyboard inputs
- Annotate on the touchscreen
- View instrument disk space and memory information

The instrument must be connected to the internet to use remote support.

Contact Support to receive a PIN prior to beginning a remote support session.

1. In the home screen, tap **⚙ (Settings) ▶ Remote support**.
The **Remote Support** screen is displayed.
2. Tap **Start remote session**, then enter the PIN provided by Support.

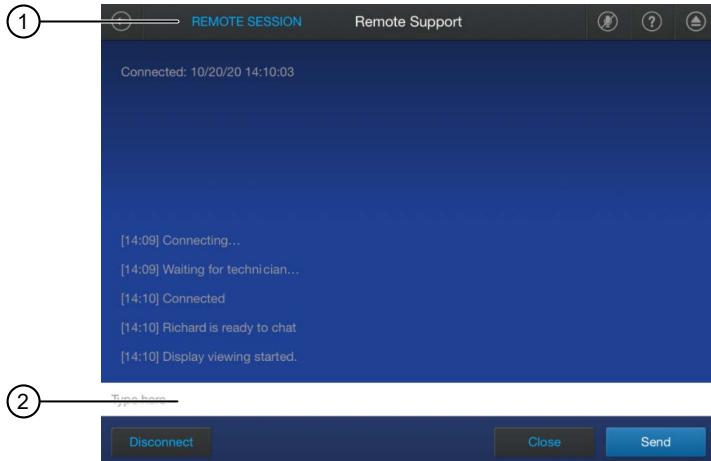

① **Remote Session**

② **Text field**

Remote Session is displayed at the top of the touchscreen when a successful connection is established. A chat is displayed on the touchscreen.

3. Tap the text field to type a message, then tap **Send**.
4. *(Optional)* Navigate to another screen on the instrument.
Remote Session is displayed at the top of the touchscreen to indicate that the session is active.
5. Tap **Remote Session** to return to the remote support session.
The **Remote Support** screen is displayed.
6. In the **Remote Support** screen, tap the active session listed in the table, then tap one of the following options.
 - **Disconnect** to end the remote support session.
 - **View** to return to the chat.

The remote session can also be disconnected from the chat screen. A Support representative can end the remote support session. **Remote Session** is not displayed on the touchscreen after a session has been disconnected.

Configure the instrument for the hands-free features

The administrator must be signed in with a Connect profile in order to enable voice commands.

1. In the home screen, tap  **(Settings)** ▶ **Hands-free operation**.
2. Set the sliders to the **Off** position or the **On** position.
 - Enable facial authentication
 - Enable voice command
3. Tap **Done**.


Manage the run history

- To view the run history, see “View the list of data files on the instrument” on page 82.
- To export the run history, see “Transfer data files at a later time” on page 81.
- To delete the run history, see “Delete data files from the instrument” on page 82.

Manage the file name convention of the data file

1. In the home screen, tap  **(Settings)** ▶ **Run settings** ▶ **File name convention**.
The **Data File Naming Convention** screen is displayed.
2. Tap **Attributes**.
The list of available attributes for the data file name is displayed.
3. Select any of the following attributes to add to the data file name format, then tap **Done**.
 - **Plate file name** (default)
 - **Time stamp** (default)
 - **User name**
 - **Barcode**
 - **Custom text**

4. Tap the lines beside an attribute, then drag the attribute above or below the other attributes in the table.

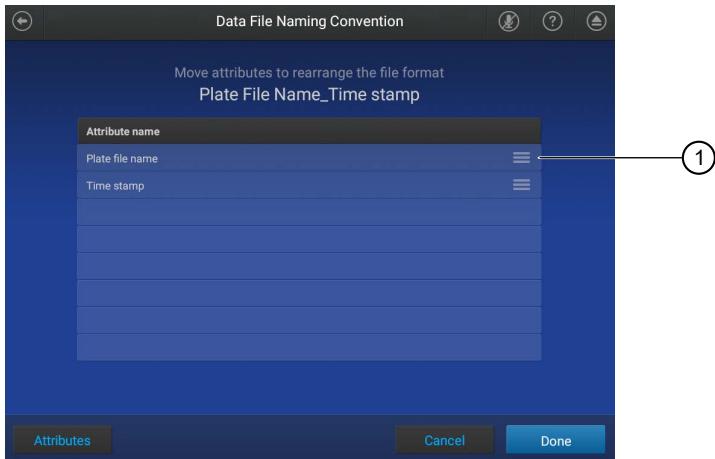

① Location to tap and drag an attribute

The order of the attributes in the table is the order that they will appear in the data file name.

5. Tap **Done**.

Enable or disable the reminder to insert a plate

1. In the home screen, tap  **(Settings)** ▶ **Run settings** ▶ **Insert plate reminder**.
The **Insert Plate Reminder** screen is displayed.
2. Set the slider to the **On** or the **Off** position.
3. Tap **OK**.

Select the language

1. In the home screen tap  **(Settings)** ▶ **Language**.
The **Language** screen is displayed.
2. Tap the language of interest.
3. Tap **Done**.



View the notifications

1. In the home screen, tap **(Settings)** ▶ **Notifications**.

The **Notifications** screen is displayed. It lists the notification and the date and the time that it was received.

2. Tap **Close**.

Enable and disable home screen notifications

When home screen notifications are enabled, the number of notifications appears on the **Settings** button.

1. In the home screen, tap **(Settings)** ▶ **Notifications**.

The **Notifications** screen is displayed.

2. Set the slider to the **On** or the **Off** position.
3. Tap **Close**.

Set up email notifications

Email notifications are only available when a user is signed in with a Connect account. The emails notifications will be sent to the email address associated with the Connect account.

1. In the home screen, tap **(Settings)** ▶ **Email notifications**.

Note: If a user is not signed in with a Connect account, the **Email notifications** button will not be displayed.

2. Select the event type.

- Run started
- PCR cycling started
- Run paused
- Run stopped
- Run is about to end
- Run is complete
- Instrument error

Note: Deselect all of the event types in order to not receive email notifications.

3. Tap **Done**.

Clear the notifications

1. In the home screen, tap  **(Settings)** ▶ **Notifications**.

The **Notifications** screen is displayed.

2. Clear the notifications.

Option	Description
Clear a single notification	Tap a single notification in the table, then tap Clear .
Clear all the notifications	Tap Clear all .

3. Tap **Close** to return to the **Settings** screen.




Instrument specifications and layout

■ Instrument dimensions	189
■ Electrical requirements	192
■ Environmental requirements	193


Instrument dimensions

The instrument, including a block, has a total weight of 35.85 kg.

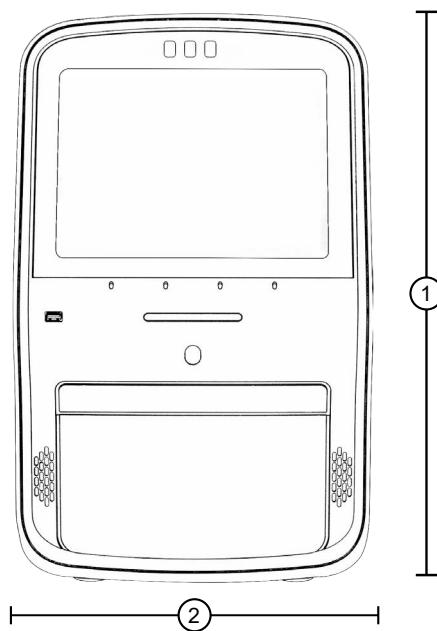

① Height: 54.7 cm

② Width: 33.7 cm

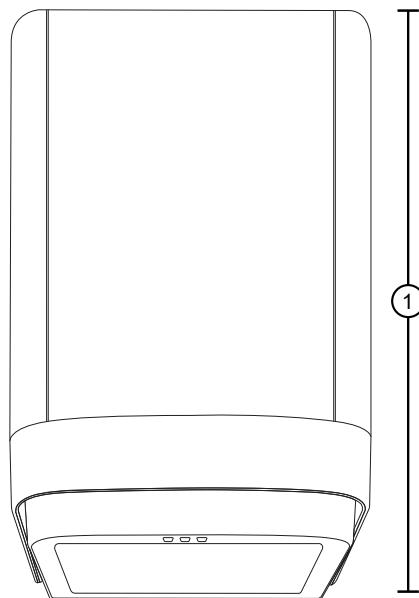

① Depth: 52.5 cm

Instrument and computer clearances

During instrument installation and maintenance, it is necessary to access the back of the instrument. If the back of the instrument faces a wall, ensure that there is sufficient clearance on the bench to rotate the instrument for access.

IMPORTANT! For safety, the power outlet for the instrument must be accessible.


Component	Top	Front	Sides	Back
Instrument	30.5 cm (12 in.)	30.5 cm (12 in.)	—	15.25 cm (6 in.)
Computer ^[1]	—	15.25 cm (6 in.)	—	15.25 cm (6 in.)

^[1] Co-locating the computer with the instrument is optional.

Configured system dimensions


Allow space for the configured instrument.

It is not a requirement to locate the computer next to the instrument. The clearances for the instrument shown in the figures also apply to an instrument without a computer.

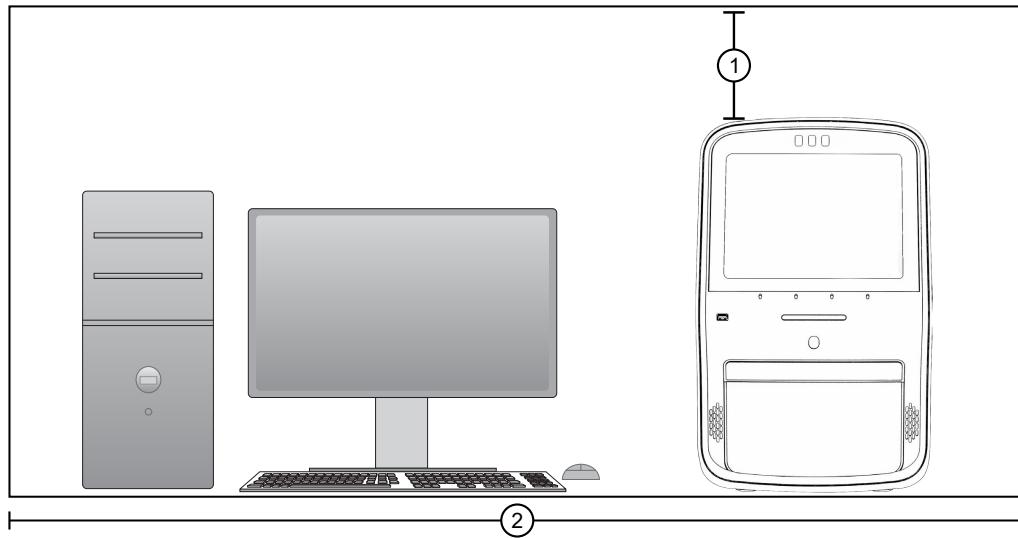

① Top clearance for the instrument: 30.5 cm

② Width of system: 185 cm

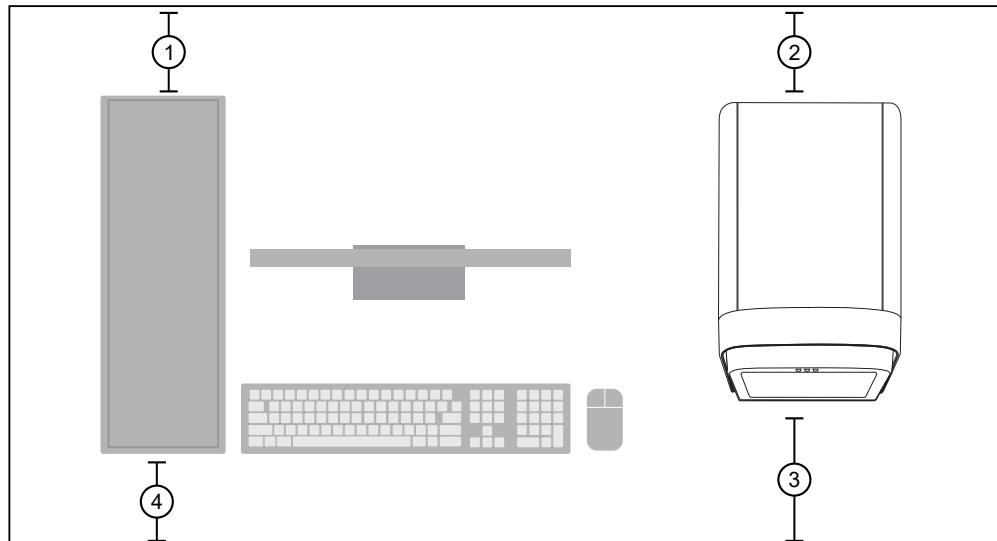

① Back clearance for the instrument: 15.25 cm

② Back clearance for the computer: 15.25 cm

③ Front clearance for the instrument: 30.5 cm

④ Front clearance for the computer: 15.25 cm

Light guidelines

Use the following guidelines to ensure that there is proper lighting for facial authentication.

- Avoid direct light above or behind the user.
- Place the instrument in a position that minimizes backlight.
- For optimal results, place the instrument between two ceiling lights.

Electrical requirements



CAUTION! Do not unpack or plug in any components until they are configured for the proper operating voltage by the service representative.



WARNING! For safety, the power outlet for the instrument must be accessible at all times. See “Instrument dimensions” on page 189 for information about the space needed between the wall and the instrument. In case of emergency, you must be able to immediately disconnect the main power supply to all the equipment. Allow adequate space between the wall and the equipment so that the power cords can be disconnected in case of emergency.

- Electric receptacle required: Grounding capability required
- Maximum power dissipation: ~417 VA, 371 W (not including computer and monitor)
- Mains AC line voltage tolerances must be up to ± 10 percent of nominal voltage

Device	Rated voltage	Circuit required	Rated frequency	Rated power
Instrument	100–240 $\pm 10\%$ VAC ^[1]	10 A	50/60 Hz	960 W
Computer (desktop)	100–240 $\pm 10\%$ VAC	10 A	50/60 Hz	125 VA
Monitor				65 VA
Computer (laptop)	100–240 $\pm 10\%$ VAC	10 A	50/60 Hz	90 VA

^[1] If the supplied power fluctuates beyond the rated voltage, a power line regulator may be required. High or low voltages can adversely affect the electronic components of the instrument.

Environmental requirements

Condition	Acceptable range
Installation site	Indoor use only
Electromagnetic interference	<p>Do not use this device in close proximity to sources of strong electromagnetic radiation (for example, unshielded intentional RF sources). Strong electromagnetic radiation may interfere with the proper operation of the device.</p> <p>This equipment has been designed and tested to CISPR 11 Class A. In a domestic environment it may cause radio interference. You may need to take measures to mitigate the interference.</p>
Altitude	Located between sea level and 2000 m (6500 ft.) above sea level
Humidity (instrument and computer)	Operation: 15%–80% (noncondensing)
Temperature (instrument and computer)	15°C to 30°C (60°F to 85°F) Note: The room temperature must not fluctuate more than 2°C over a 2-hour period.
Transient category	Installation categories II
Overvoltage category	Installation categories II
Vibration	The instrument is not adjacent to strong vibration sources, such as a centrifuge, pump, or compressor. Excessive vibration will affect instrument performance.
Pollution degree	II Install the instrument in an environment that has nonconductive pollutants such as dust particles or wood chips. Typical environments with a Pollution Degree II rating are laboratories and sales and commercial areas.
Liquid waste collection	Dispose of the polymer, buffer, reagents and any liquid waste as hazardous waste in compliance with local and national regulations.
Other conditions	<p>Ensure the room is away from any vents that could expel particulate material on the components.</p> <p>Avoid placing the instrument and computer adjacent to heaters, cooling ducts, or in direct sunlight.</p>

■ Connect the computer to the instrument directly or to a LAN	194
■ Connect the instrument to a wired network	195
■ Connect the instrument to a wireless network	196
■ Instrument and computer connections	197
■ Download and install the desktop software	198
■ Networking	199

Connect the computer to the instrument directly or to a LAN

This section describes direct wired connection of the computer provided by Thermo Fisher Scientific to the instrument or to a LAN.

Do not connect a customer-provided computer to the instrument.

1. Connect an Ethernet cable from the instrument or a LAN to the computer.
2. Power on the computer, then log in using a Windows™ Administrator account.
3. In the Windows™ desktop, right-click **My Network Places** ▶ **Properties**.
4. Right-click **Local Area Connection**, then select **Properties**.
5. Select **Internet Protocol (TCP/IP)** ▶ **Properties**.



6. Set the Internet Protocol (TCP/IP) Properties for either DHCP or Static IP communication:

Network configuration	Action
DHCP	<ol style="list-style-type: none"> 1. Select Obtain an IP address automatically. 2. Set the DNS address. If the computer obtains DNS addresses: <ul style="list-style-type: none"> • Automatically – Select Obtain DNS server address automatically. • Statically – Select Use the following DNS address, then enter the address of the preferred and alternate DNS servers (if available).
Static IP	<ol style="list-style-type: none"> 1. Select Use the following IP address. 2. In the IP Address field, enter the static IP address. 3. If necessary, enter a subnet mask. 4. If necessary, enter a static gateway address in the Default Gateway field.

7. If your network requires advanced TCP/IP setup (such as WINS), define the settings:

- a. Click **Advanced** in the Internet Protocol (TCP/IP) Properties dialog box.
- b. Define the IP Settings, DNS, and WINS tabs as instructed by your systems administrator, then click **OK**.

8. Close all dialog boxes by clicking **OK**, then re-start the computer.

The computer is now visible to other computers on the network.

Connect the instrument to a wired network

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Network configuration**.
The **Network Connections** screen is displayed.
2. Tap **Edit**.
3. Tap one of the active fields found under the **Wired** option.
4. Select a wired network connection:

Wired network connection	Action
DHCP	No further action is required.
Static IP	<ol style="list-style-type: none"> 1. In the IP Address field, enter the static IP address. 2. If necessary, enter a subnet mask. 3. If necessary, enter a static gateway address in the Default Gateway field.

5. Tap **Done**.

The **Network Connections** screen is displayed. The **Status** and **IP Address** fields under the **Wired** option have the information from the selected wired network.

6. Tap **Done**.

This instrument is now connected to the selected wired network.

Connect the instrument to a wireless network

1. In the home screen, tap  **(Settings)** ▶ **Instrument settings** ▶ **Network configuration**.

The **Network Connections** screen is displayed.

2. Tap **Edit**.

3. Tap one of the active fields found under the **Wireless** option.

4. Select a wireless network to connect to the instrument.

5. *(Optional)* Tap **Refresh** to prompt the instrument to search for available wireless networks.

6. Tap **Done**.

The **Network Connections** screen is displayed. The **Status**, **Network**, and **IP Address** fields under the **Wireless** option are filled with the information from the selected wireless network.

7. Tap **Done**.

This instrument is now connected to the selected wireless network.

Instrument and computer connections

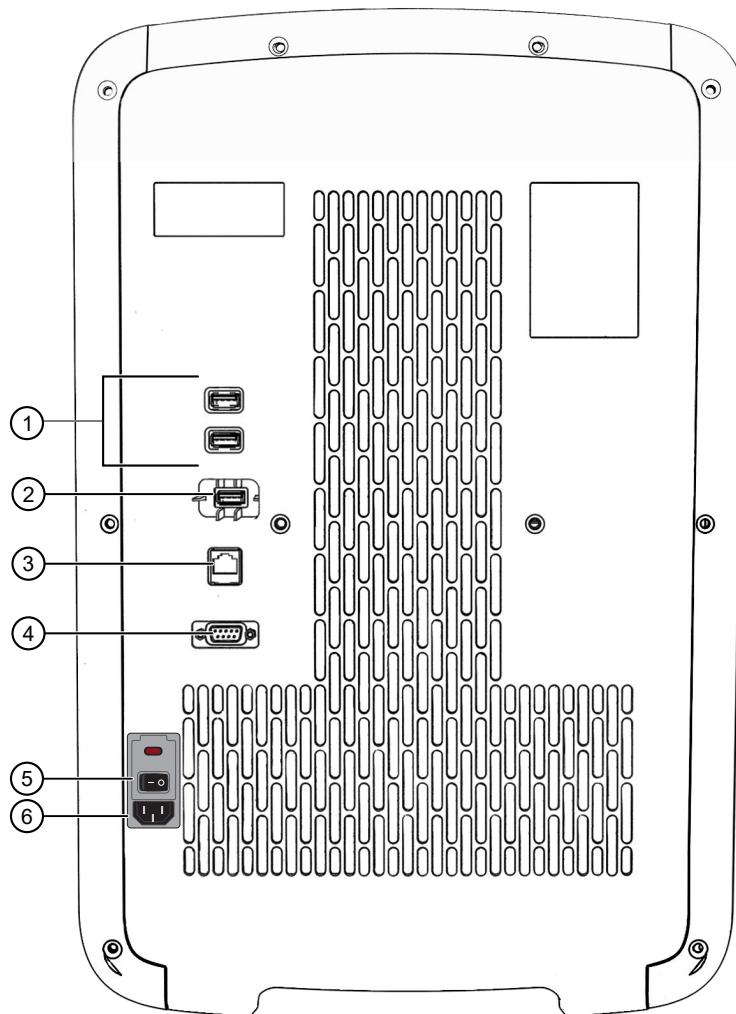


Figure 25 Instrument back panel

- ① USB ports
- ② Wi-Fi dongle port
- ③ Ethernet port
- ④ RS-232 port (service use only)
- ⑤ ON/OFF switch
- ⑥ Power inlet receptacle

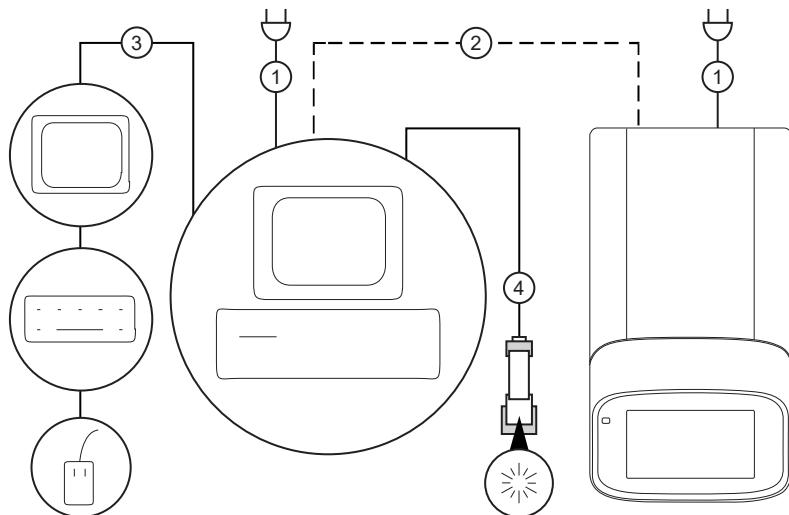


Figure 26 Instrument-to-computer connections (minitower configuration)

- ① Detachable power supply cord compatible with local power supply receptacle.
- ② Connection between the computer and the instrument.
- ③ Connection between the computer and the monitor, keyboard, and mouse.
- ④ Connection between the computer and the (*optional*) handheld barcode scanner.

Download and install the desktop software

Computer requirements for the desktop software

The desktop software can be installed on a customer-provided computer. The following list contains the minimum software requirements for a customer-provided computer.

- Operating system—Windows™ 10 (64-bit) or Macintosh™ OS 10.01
- Processor—Pentium® 4 processor or comparable
- Memory—4 GB RAM
- Hard drive—10 GB
- Monitor—1280 × 1024 resolution

Download the desktop software

1. Go to thermofisher.com/qpcrsoftware.
2. Select the link for the QuantStudio™ 6 Pro and 7 Pro Real-Time PCR Systems.
3. Download the software.

The example files are embedded in the software.

Install the desktop software

If you ordered a computer supplied by Thermo Fisher Scientific, the Field Service Engineer will configure the computer and install the desktop software during system installation.

1. Use an administrator account to log in to the computer on which you are installing the desktop software.
2. Unzip the downloaded software.
The example files are embedded in the software.
3. Double-click **setup.exe**.
4. Follow the **InstallShield Wizard** prompts to install the software.
5. Accept the License Agreement.
6. Select **Typical** as the setup preference, then click **Next**.
7. Click **Finish**.


Networking

IMPORTANT! This section provides general networking information. It *does not* provide adequate detail to integrate the instrument into all possible network architectures. Because a network may contain advanced features (such as a firewall or network domains), we recommend that you consult a network administrator before connecting the instrument to your laboratory network.




Supported options for instrument and computer connections

We support the following direct, networked (LAN—local area network), or Connect cloud-based platform configurations. Configurations other than those listed are not recommended. Select a configuration that meets the needs of your laboratory's instrument, software, and workflow requirements.

Note: For detailed information about networking your instrument, see the *QuantStudio™ 6 Pro and 7 Pro Real-Time PCR Systems IT Checklist* (Pub. No. MAN0018160).

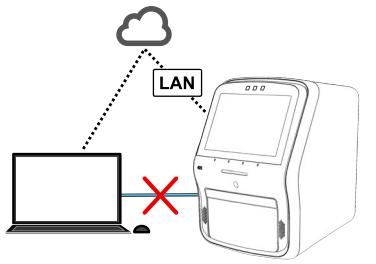Connect cloud-based platform connection	
<ul style="list-style-type: none">• Internet access and a computer with the Chrome™ web browser to access the Connect platform.• Computer-to-Connect platform connection (<i>select an option</i>):<ul style="list-style-type: none">– Wired connection to the internet using an Ethernet cable –<i>or</i>–– Wireless connection to the internet	

(continued)

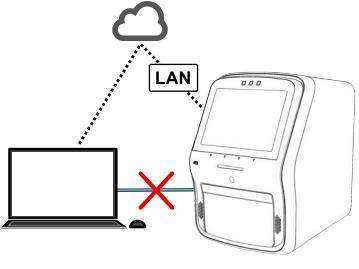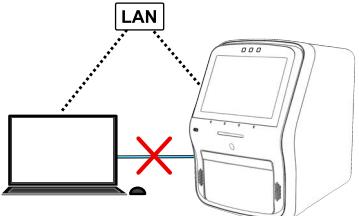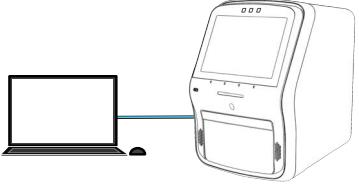Connect cloud-based platform connection	
<ul style="list-style-type: none">Instrument-to-Connect platform connection (<i>select an option</i>):<ul style="list-style-type: none">Wired connection to the network using an Ethernet cable –or–Wireless connection to the network using the instrument Wi-Fi module	
Local area network (LAN) connection	
<ul style="list-style-type: none">A computer provided by Thermo Fisher Scientific with the QuantStudio™ Design and Analysis Desktop Software.The computer and instrument must be on the same subnet mask.	
<ul style="list-style-type: none">Computer-to-LAN connection (<i>select an option</i>):<ul style="list-style-type: none">Wired connection to the network using an Ethernet cable –or–Wireless connection to the network	
<ul style="list-style-type: none">Instrument-to-LAN connection (<i>select an option</i>):<ul style="list-style-type: none">Wired connection to the network using an Ethernet cable –or–Wireless connection to the network using the instrument Wi-Fi module	
Direct connection	
<ul style="list-style-type: none">A computer provided by Thermo Fisher Scientific with the QuantStudio™ Design and Analysis Desktop Software.	
<ul style="list-style-type: none">Computer-to-instrument connection: Direct, wired connection between the computer and the instrument using an Ethernet cable.	
<p>IMPORTANT! A direct instrument-to-computer connection <i>cannot</i> be combined with the LAN or the Connect cloud-based platform configurations.</p>	



Control and monitor networked instruments

The following items apply when the instrument is connected to a network:

- Computers on the network that are running the desktop software can control the instrument. Networked instruments can be controlled by only one computer at a time.
- Instruments linked to the Connect cloud-based platform cannot be controlled remotely. The following functions can be performed:
 - Remotely access the Connect cloud-based platform to create plate files and analyze data files.
 - Send plate files to the instrument run queue, then start a run from the instrument.
 - Monitor a run in real-time from the Connect cloud-based platform.

Networking guidelines and best practices

- Consult a network administrator before connecting the instrument to a network.
- To enable the full functionality of the software, the computer requires a network connection.
- Open the firewall port for the instruments to be discovered. See “Firewall ports that must be open” on page 201.
- Observe the restrictions to mDNS and Autodiscovery.

The instrument supports mDNS but only when the instrument and computer share a direct network connection and are within the same subnet. Network computers that are separated from the instrument by a router, hub, or another network device may not be able to access the instrument by its host name.

- Confirm the uniqueness of the instrument name.
 - The instrument name must be unique within the subnet. The desktop software can automatically discover instruments on the link-local network.
 - The instrument does not test the uniqueness of the instrument name within the subnet when it is set.

Firewall ports that must be open

Ports	Condition
80/443	Standard ports for instrument-to-Connect platform and computer-to-Connect platform connections
mDNS, 7443	Instrument-to-computer connection
mDNS, 5353	Instrument discovery

Ethernet port overview

The Ethernet port of the instrument supports:

- Static IP network service with subnet mask, primary and secondary data network service (DNS), and default gateway settings, or dynamic host configuration protocol (DHCP) network service.
- mDNS/DNS for local domains.

Note: Because mDNS is limited to direct network connections, an instrument configured for mDNS may not be visible to other nodes that are separated by a router, hub, or another network device.

- IPv4 linklocal (IPV4LL) in the RFC (also known as Automatic Private IP Addressing [APIPA] or Internet Protocol Automatic Configuration [IPAC]).

Note: When an instrument is set for DHCP, APIPA is automatically enabled, and the instrument provides an IP address when no address is supplied by the DHCP server.

Third-party software

Before installing third-party software on the computer running the desktop software, confirm that the third-party software will not do the following:

- Restrict Ethernet communication.
- Interfere with instrument or computer operation.



Detailed procedures for preparation of a TaqMan™ Array Card


Guidelines for preparation of a card

- Keep the card protected from light and stored as indicated until ready for use. Excessive exposure to light may affect the fluorescent probes of the dried-down assays in the card.
- Before removing the card from its packaging:
 - Prepare each sample-specific PCR reaction mix.
 - Allow the card to reach room temperature.
- Load each fill reservoir with 100 µL of sample-specific PCR reaction mix.
 - Each fill reservoir contains a single sample as determined by the card layout.
 - The 100-µL volume ensures adequate filling of each reaction well. Volumes smaller than 100 µL result in insufficiently filled cards.
- Do not allow the micropipette tip to contact the coated foil beneath the fill port.


- Load the card with PCR reaction mix *before* centrifuging the card. During centrifugation, the PCR reaction mix resuspends the dried-down assays in each well of the card. Adding sample after centrifuging disrupts the assay layout of the card.
- Run the card within the time allowed by the PCR master mix. See the user guide for the assay or the master mix.
- If the card is not run immediately, protect it from light and store at 2-8°C.

TaqMan™ Array Card diagram

A TaqMan™ Array Card includes 8 fill reservoirs and 384 reaction wells.

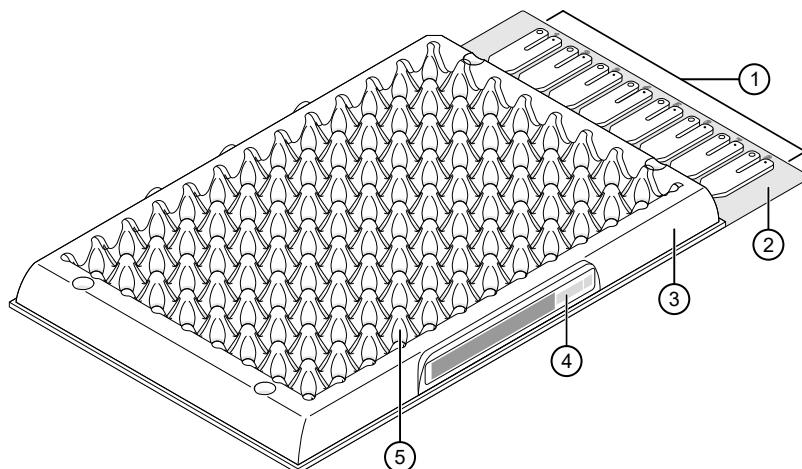


(1) Fill reservoirs (8 total)
 (2) Fill reservoir strip
 (3) Array card carrier
 (4) Array card barcode
 (5) Reaction well (384 total)

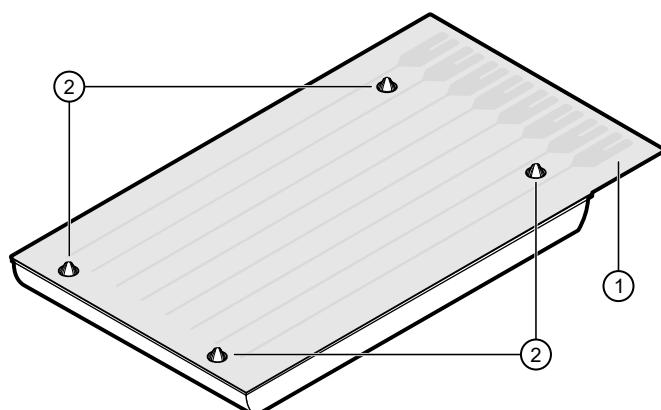

(1) Foil
 (2) Array card feet

The fill reservoir includes a fill port and a vent port. Use the fill port to load PCR reaction mix into the card.

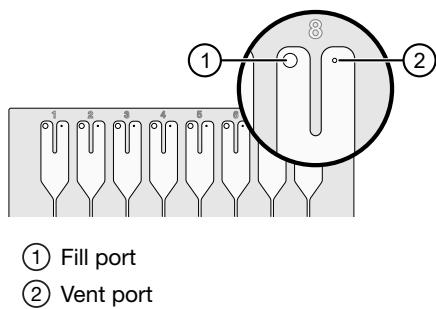


Load the PCR reaction mix

Before removing the card from its packaging:

- Prepare each sample-specific PCR reaction mix.
- Allow the card to reach room temperature.


1. Carefully remove the card from its packaging.
2. Place the card on the benchtop with its foil-side down.
3. Load 100 μ L of the sample-specific PCR reaction mix into a micropipette.
4. Hold the micropipette in an angled position, then place the tip into a fill port of the card.
5. Slowly dispense the entire volume of reaction mix so that it sweeps in and around the fill reservoir toward the vent port.

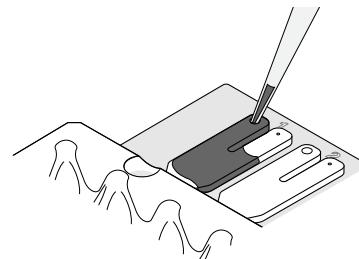

Centrifuge the card

1. Load the cards into the centrifuge buckets.
 - a. Place the bucket on the benchtop with its label facing the front of the bench.
 - b. Insert the cards into the card holder, ensuring that:
 - The fill reservoirs extend upwards out of the card holder.
 - The reaction wells face the label-side of the card holder.
 - c. Insert blank balance cards into any empty positions of the card holder. All three positions in the card holder must be filled.

d. Place the loaded card holder into the bucket so that the card holder label faces the front of the bucket.

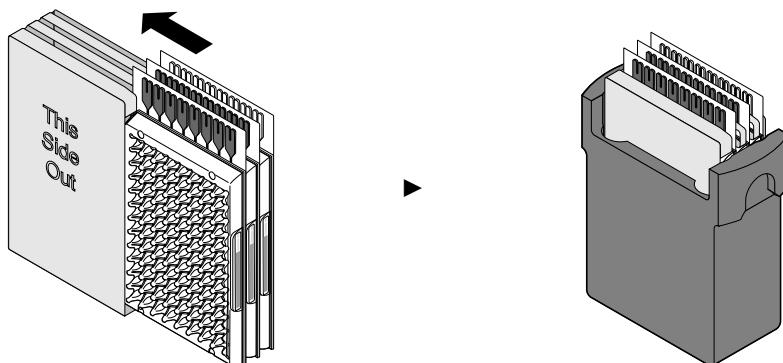

2. Configure the centrifuge using its front-panel controls.

- Set the bucket type to **15679**.
- Set the following parameters according to the control panel type.


Parameter	EASYSet (touchpad-operated)	QUIKSet (knob-operated)
Increasing ramp rate	9	3
Decreasing ramp rate	9	—
Rotational speed	1,200 rpm (331 \times g)	1,200 rpm
Centrifugation time ^[1]	1 minute	1 minute

^[1] You will centrifuge the cards twice, each time for 1 minute (see step 4).

IMPORTANT! A speed that is set too high can deform the card.

3. Load the buckets into the centrifuge.

- Press  on the centrifuge control panel to open the centrifuge cover.
- Place each loaded bucket onto an open rotor arm of the centrifuge.
Ensure that each bucket can swing easily within its slotted position on the rotor arm.
- If there are empty rotor arms, prepare buckets with blank balance cards as described in step 1. Place the balance buckets onto the rotor arms.
The rotor must be evenly loaded and opposing buckets must hold the same weight.
- Close the centrifuge cover.

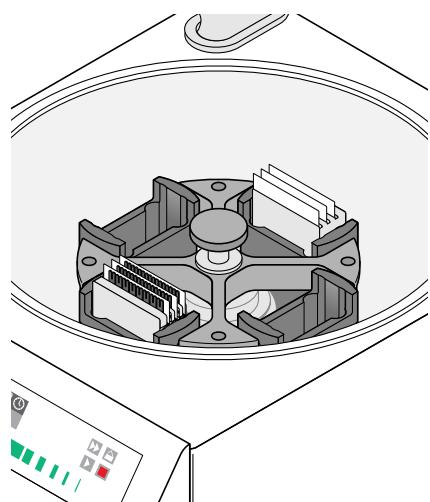

Centrifuge is properly loaded and balanced.



4. Run the centrifuge.

a. Press ►.

The centrifuge will start, then automatically stop after 1 minute.

b. Repeat substep 4a so that the cards are centrifuged for a total of two, consecutive, 1-minute centrifugations.

IMPORTANT! Do not centrifuge the cards continuously for 2 minutes. The ramping up in speed during the two, consecutive 1-minute centrifugations is important for proper filling.

5. Remove the cards from the centrifuge.


a. Press ▲.

b. Remove the buckets from the centrifuge, then remove the card holders from the buckets.

c. Remove each card from the card holder by lifting it gently by the card carrier sides.

6. Examine the cards for proper filling.

When properly filled, the remaining volumes of PCR reaction mix are consistent reservoir to reservoir.

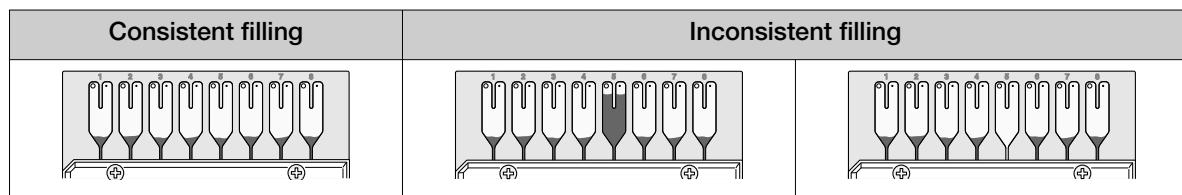

Seal the card

The TaqMan™ Array Card Sealer isolates the wells of an array card after it is loaded with PCR reaction mix and centrifuged. The sealer uses a precision stylus assembly (under the carriage) to seal the main fluid distribution channels of the array card.


Note: In some documents, the TaqMan™ Array Card Sealer is referred to as a Stylus Staker.

1. Position the TaqMan™ Array Card Sealer and its carriage before inserting a card.

a. Place the sealer on a benchtop that is approximately waist-high so that the carriage can be easily maneuvered.

b. Position the sealer with the carriage starting position toward the front of the bench.

Ensure that the engraved arrows on the sealer point away from you.

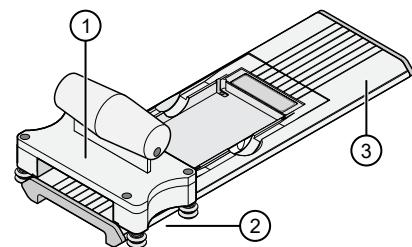

① Carriage

② Carriage starting position


③ Carriage ending position

- c. Ensure that the carriage for the sealer is at the starting position.

IMPORTANT! Do not insert a filled card into the sealer if the carriage is *not* in its starting position. The card will be irreparably damaged if the carriage is moved backwards across the card towards the starting position.

2. Insert a filled, centrifuged card into the sealer.


- a. Hold the card with its foil-side up.
- b. Orient the card over the sealer with the fill reservoirs of the card toward the ending position.
- c. Align the rear pin grooves of the card to the alignment pins of the sealer.

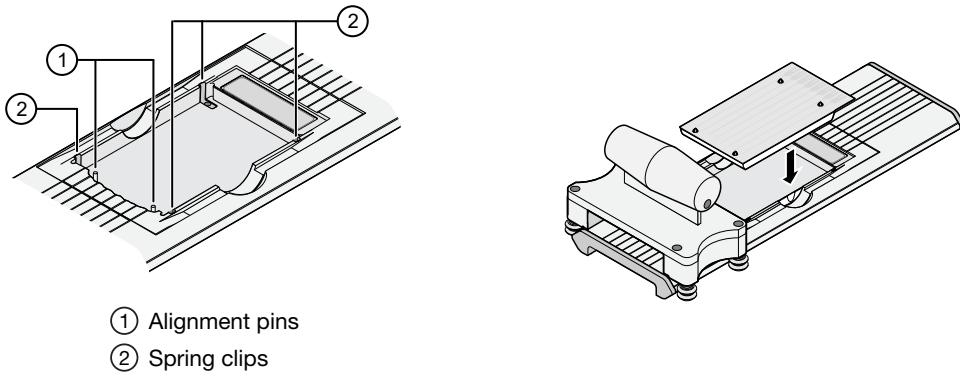

- d. Gently place the card on top of the sealer.


- e. Gently push the card until it is seated securely in the sealer.

When properly seated, the foil surface of the card is level with the base of the sealer and the spring clips hold the card securely in place.

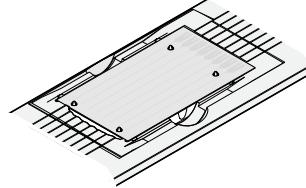

3. Slowly and steadily push the carriage across the sealer in the direction of the engraved arrows.

Push the carriage across the entire length of the card until the carriage meets the mechanical stops at the ending position.

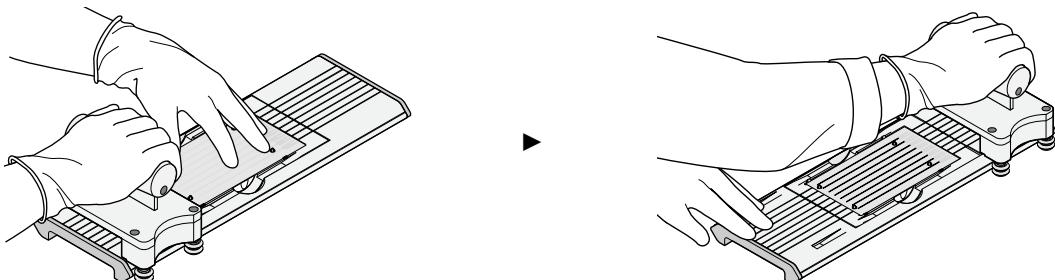



IMPORTANT!

- Do not use excessive force or speed when pushing the carriage across the card.
- Do not move the carriage back across the card. Leave the carriage at the ending position while removing the card from the sealer.

4. Remove the sealed card from the sealer by grasping the sides of the card and lifting it off.


Use the thumb slot in the middle of the sealer to access the card.

5. Examine the card for proper sealing.

Note: When properly sealed, the indentations from the sealer carriage align with the main channels of the card.

If the indentations do not align or if the foil is damaged, do not use the card.

6. Use scissors to trim the fill reservoir strip from the card. Use the edge of the card carrier as a guide.

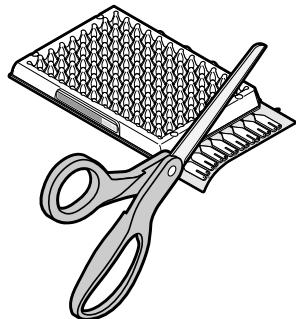

IMPORTANT! Completely remove the fill reservoir strip. Any remaining plastic that extends beyond the card edge can prevent the card from seating properly on the sample block and can affect amplification.

Correct trim	Incorrect trim

The card is now ready to run on the instrument.

(Optional) Store the sealed card, protected from light, at 2–8°C. The length of time the card can be stored depends on the master mix. See the user guide for the assay or the master mix.



Parts and materials

Kits, consumables, accessories, and reagents

Unless otherwise indicated, all materials are available through thermofisher.com.

Store all calibration and RNase P plates at -20°C. All other items can be stored at 15–30°C. Use all materials by the expiration date on the packaging.

Catalog numbers that appear as links open the web pages for those products.

96-well 0.2 mL consumables

Consumable	Amount	Cat. No.
MicroAmp™ Optical 8-Cap Strips	300 strips	4323032
MicroAmp™ Optical 8-Tube Strip, 0.2 mL	125 strips	4316567
MicroAmp™ Optical Tube without Cap, 0.2 mL	2,000 tubes	N8010933
MicroAmp™ Optical 96-Well Reaction Plate with Barcode	20 plates	4306737
	500 plates	4326659
MicroAmp™ EnduraPlate™ Optical 96-Well Clear GPLE Reaction Plates with Barcode	20 plates	4483348
	500 plates	4483351

Instrument verification or calibration plate	Cat. No.
TaqMan™ RNase P Instrument Verification Plate, 96-Well 0.2-mL	4432382
Region of Interest (ROI) and Background Plates, 96-Well 0.2-mL (2 plates)	4432364
QuantStudio™ 3/5 10-Dye Spectral Calibration Kit, 96-Well 0.2-mL (contains all 3 spectral calibration plates listed below)	A26343
QuantStudio™ 3/5 Spectral Calibration Plate 1 (FAM™, VIC™, ROX™, and SYBR™ dyes), 96-Well 0.2-mL	A26331
QuantStudio™ 3/5 Spectral Calibration Plate 2, 96-Well 0.2-mL (ABY™, JUN™, and MUSTANG PURPLE™ dyes)	A26332
QuantStudio™ 3/5 Spectral Calibration Plate 3, 96-Well 0.2-mL (TAMRA™, NED™, and Cy®5 dyes)	A26333

96-well 0.1 mL consumables

Consumable	Contents	Cat. no.
MicroAmp™ Optical 8-Cap Strips	300 strips	4323032
MicroAmp™ Fast 8-Tube Strip, 0.1 mL	125 strips	4358293
MicroAmp™ Fast Reaction Tube with Cap, 0.1 mL	1000 tubes	4358297
MicroAmp™ 96-Well Tray for VeriFlex™ Block	10 trays	4379983
MicroAmp™ Fast Optical 96-Well Reaction Plate, 0.1 mL	10 plates	4346907
MicroAmp™ EnduraPlate™ Optical 96-Well Fast Clear Reaction Plates with Barcode	20 plates	4483485
	500 plates	4483494

Calibration or instrument verification plate	Cat. no.
Region of Interest (ROI) and Background Plates, Fast 96-Well 0.1-mL (2 plates)	4432426
QuantStudio™ 3/5 10-Dye Spectral Calibration Kit, 96-well, 0.1-mL	A26342
QuantStudio™ 3/5 Spectral Calibration Plate 1, (FAM™, VIC™, ROX™, and SYBR™ dyes), 96-well, 0.1 mL	A26336
QuantStudio™ 3/5 Spectral Calibration Plate 2 (ABY™, JUN™, MUSTANG PURPLE™ dyes), 96-well Fast (0.1-mL) Plate	A26337
QuantStudio™ 3/5 Spectral Calibration Plate 3 (TAMRA™, NED™ and Cy®5 dyes) 96-well Fast (0.1-mL) Plate	A26340
TaqMan™ RNase P Instrument Verification Plate, Fast 96-Well	4351979

384-well consumables

Consumable	Contents	Cat. no.
MicroAmp™ Optical 384-Well Reaction Plate with Barcode	50 plates	4309849
	500 plates	4326270
	1000 plates	4343814
MicroAmp™ EnduraPlate™ Optical 384-Well Clear Reaction Plate with Barcode	20 plates	4483285
	500 plates	4483273

Calibration or instrument verification plate	Cat. no.
TaqMan™ RNase P Instrument Verification Plate, 384-well	4455280
Region of Interest (ROI) and Background Plates, 384-well	4432320
QuantStudio™ 5 10-Dye Spectral Calibration Kit, 384-well (contains the 2 spectral calibration plates listed below)	A26341
QuantStudio™ 5 Spectral Calibration Plate 1, (FAM™, VIC™, ROX™, TAMRA™, and SYBR™ dyes), 384-well	A26334
QuantStudio™ 5 Spectral Calibration Plate 2 (ABY™, JUN™, MUSTANG PURPLE™, NED™, and Cy®5 dyes), 384-well	A26335

TaqMan™ Array Card consumables

Go to thermofisher.com/taqmanarrays to select assays and order TaqMan™ Array Cards.

Item	Source
Calibration kit	
ViiA™ 7 Array Card Spectral Calibration Kit, includes the following components: <ul style="list-style-type: none"> Empty TaqMan™ Array Cards FAM™ dye VIC™ dye ROX™ dye ROI dye mix Background buffer FAM™/ROX™ dye VIC™/ROX™ dye 	4432314
Verification kit	
ViiA™ 7 Array Card RNaseP Verification Kit, includes the following components: <ul style="list-style-type: none"> TaqMan™ Array Card Eight (8) vials of RNase P detection reagents, one for each port in the TaqMan™ Array Card. Each vial contains the following reagents: <ul style="list-style-type: none"> PCR master mix RNase P primers A FAM™ dye-labeled probe A known concentration of human genomic DNA template 	4432265

Accessories

Item	Amount	Cat. No.
MicroAmp™ 96-Well Tray/Retainer Set	10 trays	4381850
MicroAmp™ Multi Removal Tool	1 tool	4313950
MicroAmp™ Cap Installing Tool (handle style)	1 tool	4330015
MicroAmp™ Optical Adhesive Film	25 films	4360954
	100 films	4311971
MicroAmp™ Adhesive Film Applicator	5 applicators	4333183
RT-PCR Grade Water	10 x 1.5 mL tubes	AM9935
Handheld Barcode Scanner	1 scanner	4488442

The following accessories are required for TaqMan™ Array Cards. Contact your local sales office for details.

- Centrifuge with custom buckets and card holders, one of the following:
 - Sorvall™ centrifuge
 - Multifuge™ centrifuge
 - Megafuge™ centrifuge
- TaqMan™ Array Card Sealer (referred to as the Stylus Staker in some documents)
- Blank balance TaqMan™ Array Cards (included with the instrument block upgrade / installation kit)



Safety

■ Symbols on this instrument	214
■ Safety information for instruments not manufactured by Thermo Fisher Scientific	221
■ Instrument safety	221
■ Safety and electromagnetic compatibility (EMC) standards	225
■ Chemical safety	230
■ Biological hazard safety	231



WARNING! GENERAL SAFETY. Using this product in a manner not specified in the user documentation may result in personal injury or damage to the instrument or device. Ensure that anyone using this product has received instructions in general safety practices for laboratories and the safety information provided in this document.

- Before using an instrument or device, read and understand the safety information provided in the user documentation provided by the manufacturer of the instrument or device.
- Before handling chemicals, read and understand all applicable Safety Data Sheets (SDSs) and use appropriate personal protective equipment (gloves, gowns, eye protection, and so on). To obtain SDSs, see the “Documentation and Support” section in this document.

Symbols on this instrument

Symbols may be found on the instrument to warn against potential hazards or convey important safety information. In this document, the hazard symbol is used along with one of the following user attention words.

- **CAUTION!**—Indicates a potentially hazardous situation that, if not avoided, may result in minor or moderate injury. It may also be used to alert against unsafe practices.
- **WARNING!**—Indicates a potentially hazardous situation that, if not avoided, could result in death or serious injury.
- **DANGER!**—Indicates an imminently hazardous situation that, if not avoided, will result in death or serious injury.

Standard safety symbols

Symbol and description
 CAUTION! Risk of danger. Consult the manual for further safety information.
 CAUTION! Risk of electrical shock.
 CAUTION! Hot surface.
 CAUTION! Potential biohazard.
 CAUTION! Ultraviolet light.

Symbole et description
 MISE EN GARDE ! Risque de danger. Consulter le manuel pour d'autres renseignements de sécurité.
 MISE EN GARDE ! Risque de choc électrique.
 MISE EN GARDE ! Surface chaude.
 MISE EN GARDE ! Danger biologique potentiel.
 MISE EN GARDE ! Rayonnement ultraviolet.

Additional safety symbols

Symbol and description
 CAUTION! Moving parts.
 MISE EN GARDE ! Parties mobiles.

Location of safety labels

The instrument contains warnings at the locations shown in the figures below.

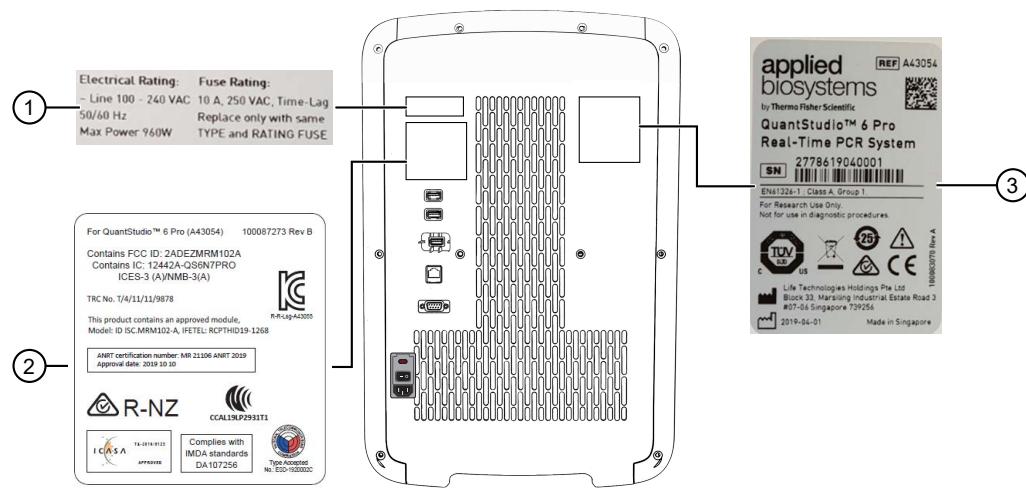


Figure 27 Labels on the back of the instrument (QuantStudio™ 6 Pro Real-Time PCR System)

- ① Electrical rating label
- ② RFID label
- ③ Safety label

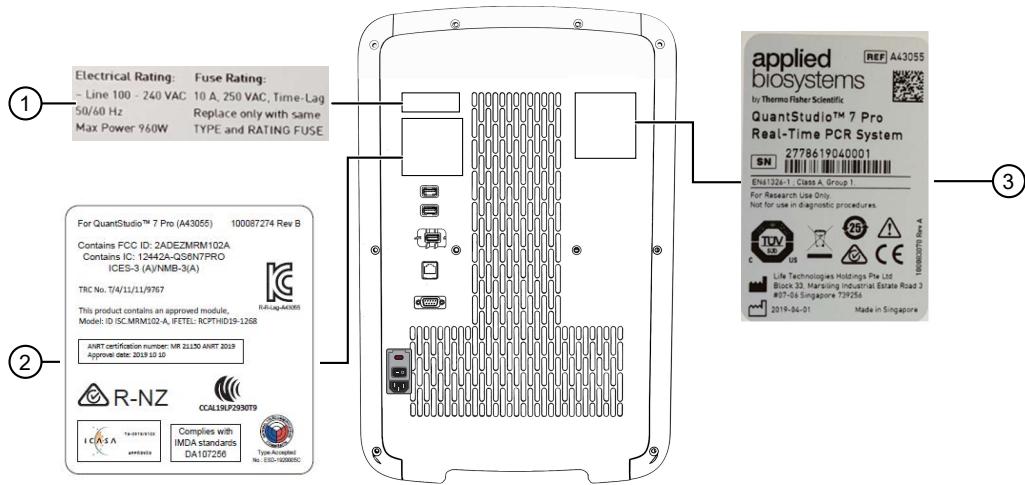


Figure 28 Labels on the back of the instrument (QuantStudio™ 7 Pro Real-Time PCR System)

- ① Electrical rating label
- ② RFID label
- ③ Safety label

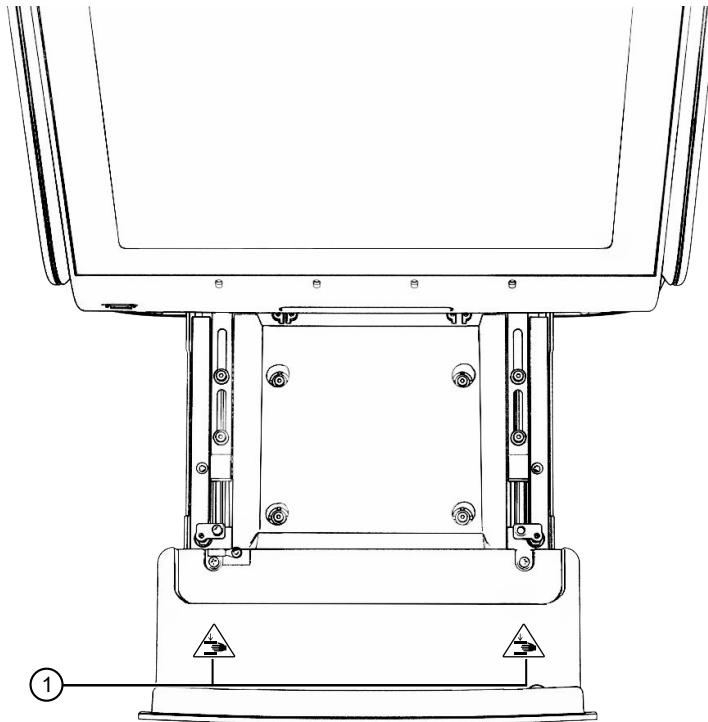


Figure 29 Labels on the drawer

- ① Moving parts label

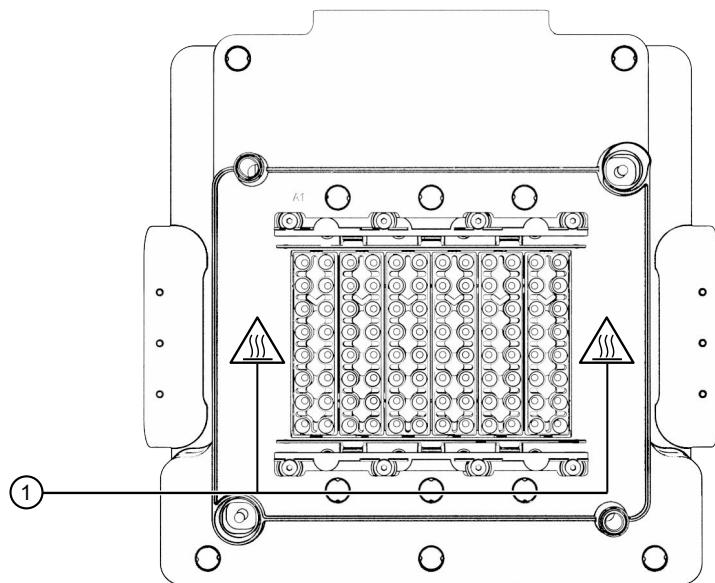


Figure 30 Labels on the top of the block

(1) Hot surface label

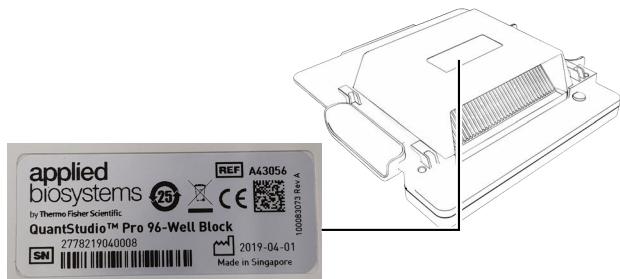


Figure 31 Labels on the bottom of the block

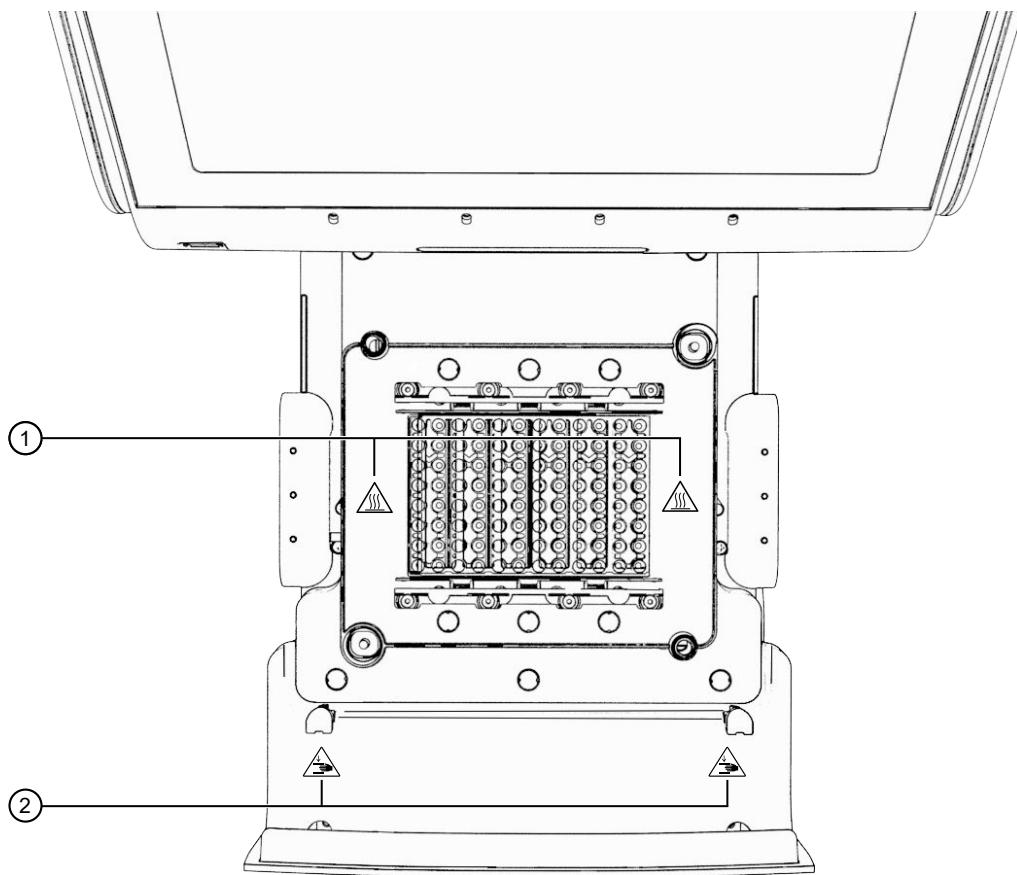


Figure 32 Labels on drawer and the installed block

- ① Hot surface label
- ② Moving parts label

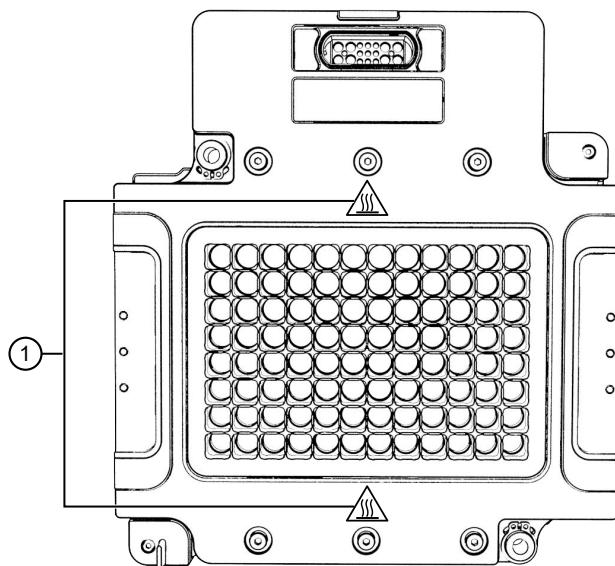


Figure 33 Labels on the top of the heated cover

- ① Hot surface label

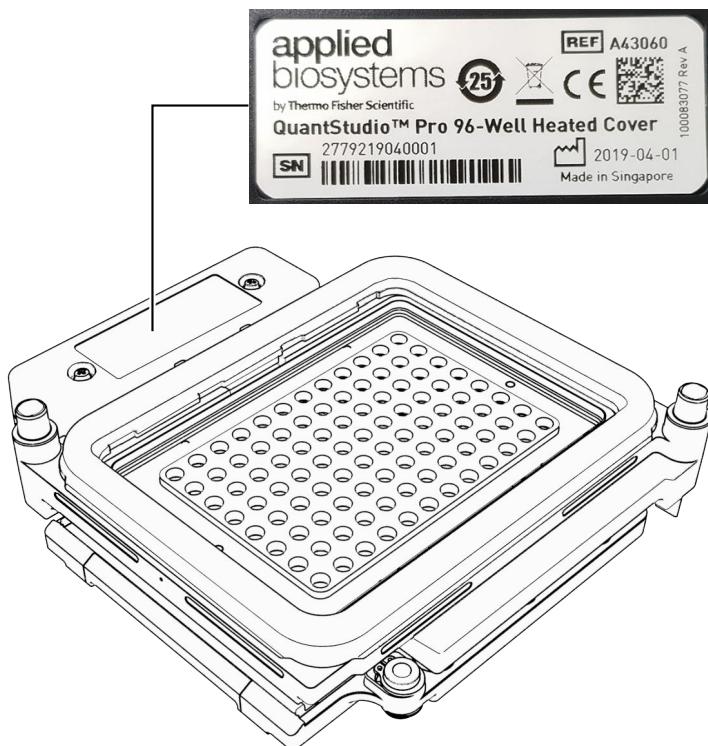


Figure 34 Labels on the bottom of the heated cover

Control and connection symbols

Symbols and descriptions	
	On (Power)
○	Off (Power)

Conformity symbols

Conformity mark	Description
	Indicates conformity with safety requirements for Canada and U.S.A.
	Indicates conformity with China RoHS requirements.
	Indicates conformity with European Union requirements.
	Indicates conformity with Australia/New Zealand standards for electromagnetic compatibility and radio frequency requirement.

(continued)

Conformity mark	Description
	Indicates conformity with the WEEE Directive 2012/19/EU.  CAUTION! To minimize negative environmental impact from disposal of electronic waste, do not dispose of electronic waste in unsorted municipal waste. Follow local municipal waste ordinances for proper disposal provision and contact customer service for information about responsible disposal options.
	Indicates conformity with US radio frequency standard.

Safety information for instruments not manufactured by Thermo Fisher Scientific

Some of the accessories provided as part of the instrument system are not designed or built by Thermo Fisher Scientific. Consult the manufacturer's documentation for the information needed for the safe use of these products.

Instrument safety

General



CAUTION! Do not remove instrument protective covers. If you remove the protective instrument panels or disable interlock devices, you may be exposed to serious hazards including, but not limited to, severe electrical shock, laser exposure, crushing, or chemical exposure.



CAUTION! Solvents and Pressurized fluids. Wear eye protection when working with any pressurized fluids. Use caution when working with any polymeric tubing that is under pressure:

- Extinguish any nearby flames if you use flammable solvents.
- Do not use polymeric tubing that has been severely stressed or kinked.
- Do not use polymeric tubing with tetrahydrofuran or nitric and sulfuric acids.
- Be aware that methylene chloride and dimethyl sulfoxide cause polymeric tubing to swell and greatly reduce the rupture pressure of the tubing.
- Be aware that high solvent flow rates (~40mL/min) may cause a static charge to build up on the surface of the tubing and electrical sparks may result.

Physical injury



CAUTION! Moving and Lifting Injury. The instrument is to be moved and positioned only by the personnel or vendor specified in the applicable site preparation guide. Improper lifting can cause painful and permanent back injury.

Things to consider before lifting or moving the instrument or accessories:

- Depending on the weight, moving or lifting may require two or more persons.
- If you decide to lift or move the instrument after it has been installed, do not attempt to do so without the assistance of others, the use of appropriate moving equipment, and proper lifting techniques.
- Ensure you have a secure, comfortable grip on the instrument or accessory.
- Make sure that the path from where the object is to where it is being moved is clear of obstructions.
- Do not lift an object and twist your torso at the same time. Keep your spine in a good neutral position while lifting with your legs.
- Participants should coordinate lift and move intentions with each other before lifting and carrying.
- For smaller packages, rather than lifting the object from the packing box, carefully tilt the box on its side and hold it stationary while someone else slides the contents out of the box.



CAUTION! Moving Parts. Moving parts can crush, pinch and cut. Keep hands clear of moving parts while operating the instrument. Disconnect power before servicing.



Electrical safety



WARNING! Fuse Installation. Before installing the instrument, verify that the fuses are properly installed and the fuse voltage matches the supply voltage. Replace fuses only with the type and rating specified for the unit. Improper fuses can damage the instrument wiring system and cause a fire.



AVERTISSEMENT ! Installation des fusibles. Avant d'installer l'instrument, vérifier que les fusibles sont correctement insérés et que leur tension correspond à celle fournie par le circuit d'alimentation. Ne remplacer les fusibles que par des modèles du type et de la puissance spécifiés pour l'appareil. L'utilisation de fusibles inadaptés peut endommager le circuit électrique de l'instrument et provoquer un incendie.



WARNING! Ensure appropriate electrical supply. For safe operation of the instrument:

- Plug the system into a properly grounded receptacle with adequate current capacity.
- Ensure the electrical supply is of suitable voltage.
- Never operate the instrument with the ground disconnected. Grounding continuity is required for safe operation of the instrument.



AVERTISSEMENT ! Veiller à utiliser une alimentation électrique appropriée. Pour garantir le fonctionnement de l'instrument en toute sécurité :

- Brancher le système sur une prise électrique correctement mise à la terre et de puissance adéquate.
- S'assurer que la tension électrique est convenable.
- Ne jamais utiliser l'instrument alors que le dispositif de mise à la terre est déconnecté. La continuité de la mise à la terre est impérative pour le fonctionnement de l'instrument en toute sécurité.



WARNING! Power Supply Line Cords. Use properly configured and approved line cords for the power supply in your facility.



AVERTISSEMENT ! Cordons d'alimentation électrique. Utiliser des cordons d'alimentation adaptés et approuvés pour raccorder l'instrument au circuit électrique du site.



WARNING! Disconnecting Power. To fully disconnect power either detach or unplug the power cord, positioning the instrument such that the power cord is accessible.



AVERTISSEMENT ! Déconnecter l'alimentation. Pour déconnecter entièrement l'alimentation, détacher ou débrancher le cordon d'alimentation. Placer l'instrument de manière à ce que le cordon d'alimentation soit accessible.

Laser safety



WARNING! LASER HAZARD. The QuantStudio™ 6 Pro System and the QuantStudio™ 7 Pro System are compatible with an optional Handheld Barcode Scanner. Lasers can burn the retina, causing permanent blind spots. To ensure safe laser operation:

- Never look directly into the laser beam.
- Do not remove safety labels, instrument protective panels, or defeat safety interlocks.
- The system must be installed and maintained by a Thermo Fisher Scientific Technical Representative.
- Remove jewelry and other items that can reflect a laser beam into your eyes or those of others
- Wear proper eye protection and post a laser warning sign at the entrance to the laboratory if the laser protection is defeated for servicing
- DO NOT operate the laser when it cannot be cooled by its cooling fan; an overheated laser can cause severe burns on contact.

The following table lists laser safety symbols and alerts that may be present on the instrument.

Alert
<p> CAUTION! LASER HAZARD, Bar Code Scanner. The bar code scanner included with the instrument is a Class 2 laser. To avoid damage to eyes, do not stare directly into the beam or point into another person's eyes.</p>
<p> MISE EN GARDE ! RISQUE LIÉ AU RAYONNEMENT LASER, Lecteur de code-barres. Le lecteur de code-barres inclus dans l'instrument est un appareil laser de classe 2. Pour éviter toute lésion oculaire, ne regardez pas directement le faisceau et ne le dirigez pas vers les yeux d'une autre personne.</p>

Cleaning and decontamination



CAUTION! Cleaning and Decontamination. Use only the cleaning and decontamination methods that are specified in the manufacturer user documentation. It is the responsibility of the operator (or other responsible person) to ensure that the following requirements are met:

- No decontamination or cleaning agents are used that can react with parts of the equipment or with material that is contained in the equipment. Use of such agents could cause a HAZARD condition.
- The instrument is properly decontaminated a) if hazardous material is spilled onto or into the equipment, and/or b) before the instrument is serviced at your facility or is sent for repair, maintenance, trade-in, disposal, or termination of a loan. Request decontamination forms from customer service.
- Before using any cleaning or decontamination methods (except methods that are recommended by the manufacturer), confirm with the manufacturer that the proposed method will not damage the equipment.



MISE EN GARDE ! Nettoyage et décontamination. Utiliser uniquement les méthodes de nettoyage et de décontamination indiquées dans la documentation du fabricant destinée aux utilisateurs. L'opérateur (ou toute autre personne responsable) est tenu d'assurer le respect des exigences suivantes:

- Ne pas utiliser d'agents de nettoyage ou de décontamination susceptibles de réagir avec certaines parties de l'appareil ou avec les matières qu'il contient et de constituer, de ce fait, un DANGER.
- L'instrument doit être correctement décontaminé a) si des substances dangereuses sont renversées sur ou à l'intérieur de l'équipement, et/ou b) avant de le faire réviser sur site ou de l'envoyer à des fins de réparation, de maintenance, de revente, d'élimination ou à l'expiration d'une période de prêt (des informations sur les formes de décontamination peuvent être demandées auprès du Service clientèle).
- Avant d'utiliser une méthode de nettoyage ou de décontamination (autre que celles recommandées par le fabricant), les utilisateurs doivent vérifier auprès de celui-ci qu'elle ne risque pas d'endommager l'appareil.

Instrument component and accessory disposal

To minimize negative environmental impact from disposal of electronic waste, do not dispose of electronic waste in unsorted municipal waste. Follow local municipal waste ordinances for proper disposal provision and contact customer service for information about responsible disposal options.

Safety and electromagnetic compatibility (EMC) standards

The instrument design and manufacture complies with the following standards and requirements for safety and electromagnetic compatibility.

Safety standards

Reference	Description
EU Directive 2014/35/EU	European Union “Low Voltage Directive”
IEC 61010-1 EN 61010-1 UL 61010-1 CAN/CSA C22.2 No. 61010-1	<i>Safety requirements for electrical equipment for measurement, control, and laboratory use – Part 1: General requirements</i>
IEC 61010-2-010 EN 61010-2-010	<i>Safety requirements for electrical equipment for measurement, control and laboratory use – Part 2-010: Particular requirements for laboratory equipment for the heating of materials</i>
IEC 61010-2-020 EN 61010-2-020	<i>Safety requirements for electrical equipment for measurement, control and laboratory use – Part 2-020: Particular requirements for laboratory centrifuges</i>
IEC 61010-2-081 EN 61010-2-081	<i>Safety requirements for electrical equipment for measurement, control and laboratory use – Part 2-081: Particular requirements for automatic and semi-automatic laboratory equipment for analysis and other purposes</i>

EMC standards

Reference	Description
EMC Directive 2014/30/EU	European Union “EMC Directive”
IEC 61326-1	<i>Electrical Equipment for Measurement, Control and Laboratory Use – EMC Requirements – Part 1: General Requirements</i>
AS/NZS CISPR 11	<i>Limits and Methods of Measurement of Electromagnetic Disturbance Characteristics of Industrial, Scientific, and Medical (ISM) Radiofrequency Equipment</i>
ICES-003, Issue 6	<i>Industrial, Scientific and Medical (ISM) Radio Frequency Generators</i>
FCC Part 15 Subpart B (47 CFR)	<i>U.S. Standard Radio Frequency Devices</i>



Environmental design standards

Reference	Description
Directive 2012/19/EU	European Union “WEEE Directive”—Waste electrical and electronic equipment
Directive 2011/65/EU	European Union “RoHS Directive”—Restriction of hazardous substances in electrical and electronic equipment
SJ/T 11364-2014	“China RoHS” Standard—Marking for the Restricted Use of Hazardous Substances in Electronic and Electrical Products For instrument specific certificates, visit our customer resource page at www.thermofisher.com/us/en/home/technical-resources/rohs-certificates.html .

Radio compliance standards

Reference	Description
Directive 2014/53/EU	European Union “RE Directive”—Radio equipment Hereby, Thermo Fisher Scientific declares that the radio equipment type: QuantStudio™ 6 Pro Real-Time PCR System and QuantStudio™ 7 Pro Real-Time PCR System are in compliance with Radio Equipment Directive 2014/53/EU. The full text of the EU declaration of conformity is available at thermofisher.com .
FCC Part 15	Contains FCC ID: 2ADEZMRM102A FCC Notice (for U.S. Customers): This device complies with Part 15 of the FCC Rules: Operation is subject to the following conditions: <ol style="list-style-type: none"> 1. This device may not cause harmful interference, and 2. This device must accept any interference received, including interference that may cause undesired operation. Changes and modifications not expressly approved by Thermo Fisher Scientific can void your authority to operate this equipment under Federal Communications Commissions rules. Note: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense. This equipment complies with FCC radiation exposure limits set forth for an uncontrolled environment and meets the 47 CFR 2.1091 of the FCC radio frequency (RF) Exposure rules.

(continued)

Reference	Description
ICES (Innovation, Science and Economic Development Canada)	<p>This device contains licence-exempt transmitter(s)/receiver(s) that comply with Innovation, Science and Economic Development Canada's licence-exempt RSS(s). Operation is subject to the following two conditions:</p> <ol style="list-style-type: none"> 1. This device may not cause interference. 2. This device must accept any interference, including interference that may cause undesired operation of the device. <p>L'émetteur/récepteur exempt de licence contenu dans le présent appareil est conforme aux CNR d'Innovation, Sciences et Développement économique Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes:</p> <ol style="list-style-type: none"> 1. L'appareil ne doit pas produire de brouillage; 2. L'appareil doit accepter tout brouillage radioélectrique subi, même si le brouillage est susceptible d'en compromettre le fonctionnement. <p>This product complies to ICES-3 (A)/NMB-3(A). This equipment complies with ISED (IC) radiation exposure limits set forth for an uncontrolled environment and meets the ESS-102 of the ISED (IC) radio frequency (RF) Exposure rules.</p>
IMDA (Infocomm Media Development Authority) (Singapore)	Complies with IMDA standards DA107256
ANATEL (Agencia Nacional de Telecomunicacoes) (Brazil)	<p>Módulo Modelo: ID ISC.MRM102-A</p>  <p>06094-19-08681</p> <p>Este equipamento não tem direito à proteção contra interferência prejudicial e não pode causar interferência em sistemas devidamente autorizados.</p>
KCC (Korea Communications Commission) (Korea)	이 기기는 업무용 환경에서 사용할 목적으로 적합성평가를 받은 기기로서 가정용 환경에서 사용하는 경우 전파간섭의 우려가 있습니다.
MIC (Ministry of Internal Affairs and Communications) (Japan)	This product includes approved RFID module with MIC approval number AC-19073.
NCC (National Communications Commission) (Taiwan)	<p>低功率電波輻射性電機管理辦法</p> <p>第十二條經型式認證合格之低功率射頻電機，非經許可，公司、商號或使用者均不得擅自變更頻率、加大功率或變更原設計之特性及功能。</p> <p>第十四條低功率射頻電機之使用不得影響飛航安全及干擾合法通信；經發現有干擾現象時，應立即停用，並改善至無干擾時方得繼續使用。</p> <p>前項合法通信，指依電信規定作業之無線電信。低功率射頻電機須忍受合法通信或工業、科學及醫療用電波輻射性電機設備之干擾。</p>



RF transceiver specifications

RFID readers	
Description	ID ISC.MRM102-A RFID module with asynchronous RS232 interface
Frequency	13.56 MHz
Modulation	ASK
Transmit power	54.5 dBuV/m at 10m
Antenna	3674.000.00 (ID ISC. ANT40/30 HF PCB Antenna) Dimensions: 40 mm x 30 mm (1.57 inches x 1.18 inches) Antenna: 50 Ohm
Supply voltage	12 Vdc

Chemical safety



WARNING! GENERAL CHEMICAL HANDLING. To minimize hazards, ensure laboratory personnel read and practice the general safety guidelines for chemical usage, storage, and waste provided below. Consult the relevant SDS for specific precautions and instructions:

- Read and understand the Safety Data Sheets (SDSs) provided by the chemical manufacturer before you store, handle, or work with any chemicals or hazardous materials. To obtain SDSs, see the "Documentation and Support" section in this document.
- Minimize contact with chemicals. Wear appropriate personal protective equipment when handling chemicals (for example, safety glasses, gloves, or protective clothing).
- Minimize the inhalation of chemicals. Do not leave chemical containers open. Use only with sufficient ventilation (for example, fume hood).
- Check regularly for chemical leaks or spills. If a leak or spill occurs, follow the manufacturer cleanup procedures as recommended in the SDS.
- Handle chemical wastes in a fume hood.
- Ensure use of primary and secondary waste containers. (A primary waste container holds the immediate waste. A secondary container contains spills or leaks from the primary container. Both containers must be compatible with the waste material and meet federal, state, and local requirements for container storage.)
- After emptying a waste container, seal it with the cap provided.
- Characterize (by analysis if needed) the waste generated by the particular applications, reagents, and substrates used in your laboratory.
- Ensure that the waste is stored, transferred, transported, and disposed of according to all local, state/provincial, and/or national regulations.
- **IMPORTANT!** Radioactive or biohazardous materials may require special handling, and disposal limitations may apply.



AVERTISSEMENT ! PRÉCAUTIONS GÉNÉRALES EN CAS DE MANIPULATION DE PRODUITS CHIMIQUES.

Pour minimiser les risques, veiller à ce que le personnel du laboratoire lise attentivement et mette en œuvre les consignes de sécurité générales relatives à l'utilisation et au stockage des produits chimiques et à la gestion des déchets qui en découlent, décrites ci-dessous. Consulter également la FDS appropriée pour connaître les précautions et instructions particulières à respecter :

- Lire et comprendre les fiches de données de sécurité (FDS) fournies par le fabricant avant de stocker, de manipuler ou d'utiliser les matériaux dangereux ou les produits chimiques. Pour obtenir les FDS, se reporter à la section « Documentation et support » du présent document.
- Limiter les contacts avec les produits chimiques. Porter des équipements de protection appropriés lors de la manipulation des produits chimiques (par exemple : lunettes de sûreté, gants ou vêtements de protection).
- Limiter l'inhalation des produits chimiques. Ne pas laisser les récipients de produits chimiques ouverts. Ils ne doivent être utilisés qu'avec une ventilation adéquate (par exemple, sorbonne).
- Vérifier régulièrement l'absence de fuite ou d'écoulement des produits chimiques. En cas de fuite ou d'écoulement d'un produit, respecter les directives de nettoyage du fabricant recommandées dans la FDS.
- Manipuler les déchets chimiques dans une sorbonne.

- Veiller à utiliser des récipients à déchets primaire et secondaire. (Le récipient primaire contient les déchets immédiats, le récipient secondaire contient les fuites et les écoulements du récipient primaire. Les deux récipients doivent être compatibles avec les matériaux mis au rebut et conformes aux exigences locales, nationales et communautaires en matière de confinement des récipients.)
- Une fois le récipient à déchets vidé, il doit être refermé hermétiquement avec le couvercle fourni.
- Caractériser (par une analyse si nécessaire) les déchets générés par les applications, les réactifs et les substrats particuliers utilisés dans le laboratoire.
- Vérifier que les déchets sont convenablement stockés, transférés, transportés et éliminés en respectant toutes les réglementations locales, nationales et/ou communautaires en vigueur.
- **IMPORTANT !** Les matériaux représentant un danger biologique ou radioactif exigent parfois une manipulation spéciale, et des limitations peuvent s'appliquer à leur élimination.



WARNING! HAZARDOUS WASTE (from instruments). Waste produced by the instrument is potentially hazardous. Follow the guidelines noted in the preceding General Chemical Handling warning.



WARNING! 4L Reagent and Waste Bottle Safety. Four-liter reagent and waste bottles can crack and leak. Each 4-liter bottle should be secured in a low-density polyethylene safety container with the cover fastened and the handles locked in the upright position.

Biological hazard safety



WARNING! Potential Biohazard. Depending on the samples used on this instrument, the surface may be considered a biohazard. Use appropriate decontamination methods when working with biohazards.



WARNING! BIOHAZARD. Biological samples such as tissues, body fluids, infectious agents, and blood of humans and other animals have the potential to transmit infectious diseases. Conduct all work in properly equipped facilities with the appropriate safety equipment (for example, physical containment devices). Safety equipment can also include items for personal protection, such as gloves, coats, gowns, shoe covers, boots, respirators, face shields, safety glasses, or goggles. Individuals should be trained according to applicable regulatory and company/ institution requirements before working with potentially biohazardous materials. Follow all applicable local, state/provincial, and/or national regulations. The following references provide general guidelines when handling biological samples in laboratory environment.

- U.S. Department of Health and Human Services, *Biosafety in Microbiological and Biomedical Laboratories (BMBL)*, 6th Edition, HHS Publication No. (CDC) 300859, Revised June 2020; found at:
<https://www.cdc.gov/labs/pdf/CDC-BiosafetyMicrobiologicalBiomedicalLaboratories-2020-P.pdf>
- Laboratory biosafety manual, fourth edition. Geneva: World Health Organization; 2020 (Laboratory biosafety manual, fourth edition and associated monographs); found at:
www.who.int/publications/i/item/9789240011311



Documentation and support

Related documentation

Document	Pub. No.
<i>QuantStudio™ 6 Pro Real-Time PCR System and QuantStudio™ 7 Pro Real-Time PCR System Hands-Free Features Quick Reference</i>	MAN0018430
<i>TaqMan™ Array Plates with RFID for use with QuantStudio™ 6 Pro and 7 Pro Real-Time PCR Systems Quick Reference Guide</i>	MAN0018436
<i>QuantStudio™ 6 Pro and 7 Pro Real-Time PCR Systems Connectivity Quick Reference</i>	MAN0018431
<i>QuantStudio™ 6 Pro Real-Time PCR System and QuantStudio™ 7 Pro Real-Time PCR System Site Preparation Guide</i>	MAN0017992
Desktop software	
<i>QuantStudio™ Design and Analysis Software v2 User Guide</i>	MAN0018200
<i>QuantStudio™ Design and Analysis Software v2 Standard Curve Analysis Module User Guide</i>	MAN0018746
<i>QuantStudio™ Design and Analysis Software v2 Relative Quantification Analysis Module User Guide</i>	MAN0018747
<i>QuantStudio™ Design and Analysis Software v2 Presence Absence Analysis Module User Guide</i>	MAN0018748
<i>QuantStudio™ Design and Analysis Software v2 Genotyping Analysis Module User Guide</i>	MAN0018749
<i>QuantStudio™ Design and Analysis Software v2 High Resolution Melt Analysis Module User Guide</i>	MAN0018981
Connect cloud-based platform	
<i>QuantStudio™ Design and Analysis Software v2 User Guide</i>	MAN0018202
Security, Auditing, and E-signature (SAE) Administrator Console	
<i>SAE Administrator Console v2.0 User Guide</i>	MAN0017468
About the Assay Information File	
<i>Understanding Your Shipment</i>	MAN0017153



Customer and technical support

Visit thermofisher.com/support for the latest service and support information.

- Worldwide contact telephone numbers
- Product support information
 - Product FAQs
 - Software, patches, and updates
 - Training for many applications and instruments
- Order and web support
- Product documentation
 - User guides, manuals, and protocols
 - Certificates of Analysis
 - Safety Data Sheets (SDSs; also known as MSDSs)

Note: For SDSs for reagents and chemicals from other manufacturers, contact the manufacturer.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

Index

A

accessories 213
Add assay 66
Add data collection location 58
Add facial authentication 132
Add pause 62
Add PCR stage 58
Add PCR step 58
Add profile 134
Add reagent details 66, 68
Add target 66
Adjust speaker volume 169
Administrator privilege
 Disable 135
 Enable 135
Administrator profile, Create 131
Apply reagents 67, 68
array card
 centrifuge 205
 diagram 204
 guidelines 203
 load reaction mix 205
 seal 207
Assay
 Add 66
 Edit 65, 66
audit reason, specify 151
Auto Delta
 Remove 61
 Set up 61
Auto export setting
 Delete 55
 Import 55
AutoDelta, Overview 19
Automatic disk storage management 171
Automatic sign-out
 Edit 171
 Enable 171

B

Back up instrument 181
background calibration
 create plate 96

 plate, prepare 91
Background calibration
 Description 86
 Frequency 88
 Reminders 88
 Run 91, 100
 Status 87
 When to perform 85
Barcode scan 51
biohazard safety 231
Block
 Change 157
 Clean 159
 Remove 158
Block calibration 84

C

calibration, plate, prepare 91
Calibration
 Background 91, 100
 Custom dye
 Dilution guidelines 114
 Optimal concentration 116
 Perform 118
 Prepare plate 117
 Run dilution plate 116
 Workflow 114
 Custom dye dilution, Prepare dilution plate 115
 Custom melt curve 120
 Descriptions 86
 Dye 91, 100
 Export report 89
 Frequency 88
 Reminders, Connect 88
 Required materials 90, 98
 ROI/Uniformity 91, 100
 Schedule 85
 Status 87
 Transfer data files 94
 Transfer results 94
 View images 93, 101
 Workflow 90, 97
calibration failure, troubleshoot 95, 102
Calibration plates 90, 98

- Cancel scheduled run 75
- centrifuge the card 205
- Certificate, Upload 182
- Change block 157
- Change password
 - SAE, Change password 149
- Change PIN 133
- Clean block 159
- Clear notifications 188
- Cloud 26
- Comparative Ct run 32
- Comparison 16
- computer
 - connection to instrument or LAN 199
 - connections 197
 - requirements for desktop software 198
 - set up 194
- Configure hands-free features 185
- Configure network 169
- Configure ramp rates 62
- conformity symbols 220
- Connect
 - Calibration reminders 88
 - Link 130
 - Link to instrument 126
 - Roles
 - Administrator 124
 - Standard 124
 - Unlink from instrument 129
- Connect a proxy server 170
- Connect platform
 - Features 26
 - Privacy policy 167
- Connect platform connection 31
- Connect platform region 170
- Connect profile
 - Edit 133
 - Overview 123
- connect SAE server 144
- Connections
 - Options 31
 - Wired 31
 - Wireless 31
- consumables
 - 384-well 211, 212
 - 96-well 0.1-mL 211
 - 96-well 0.2-mL 210
- contamination, identify 95
- control and connection symbols 220
- Copy plate file 77
- Cover temperature 57
- Create and run a plate 47
- Custom dye, Add to instrument 118
- Custom dye calibration
- Dilution, Guidelines 114
- Dilution plate
 - Prepare plate 115
 - Run 116
- Optimal concentration 116
- Perform 118
- Prepare plate 117
- Workflow 114
- Custom dyes, Overview 113
- Custom melt curve, Calibrate 120
- Cycles 57

D

- Data
 - View during run 73
 - Zoom 74
- Data collection 18
- Data collection location
 - Add 58
 - Remove 58
- Data file
 - Delete 82
 - Transfer 81
- Data file name, Manage 185, 186
- Date
 - Format 168
 - Set 168
- Delete auto export setting 55
- Delete data file 82
- Delete plate file 78
- Delete profile 135
- Delete reagent details 67
- desktop software
 - download 198
 - download and install 198
 - install 199
- Desktop software, Features 26
- Direct network connection 31
- Disable administrator privilege 135
- Disable demonstration mode 168
- Disable email notifications 187
- Disable home screen notifications 187
- Disable remote monitoring 180
- Disable SAE 145
- Disk storage management 171
- disposal of instrument components and accessories 225
- documentation, related 232
- download desktop software 198
- Dye
 - Custom, Add 118
 - Edit 68
 - Passive reference 68

dye calibration, plate, prepare 91

Dye calibration

 Description 86

 Frequency 88

 Reminders 88

 Run 91, 100

 Status 87

 When to perform 85

Dyes

 Custom 17, 113

 System 16

E

Edit automatic sign-out 171

Edit guest access 122

Edit method 57–59, 61, 62

Edit passive reference dye 68

Edit plate file 71

Edit plate file during run 71

Edit plate layout 65, 66

Edit profile 133

Edit reagent details 66–68

Edit reagents 67, 68

Edit sleep mode 163

Email notifications 187

Enable administrator privilege 135

Enable automatic sign-out 171

Enable demonstration mode 168

Enable email notifications 187

Enable home screen notifications 187

Enable remote monitoring 180

Enable sleep mode 163

Ethernet port 202

EULA, View 167

Export calibration report 89

Export instrument log 179

F

Facial authentication

 Edit 132

 Indicators 23

 Overview 22

 Set up 131

 Sign in 44

Features 16

File locations 27

File name convention, Data file 185, 186

File names, System templates 34, 35, 37, 38

Files

 Description 30

 Features 28

firewall ports 201

fluorescence high

Functions, SAE 138, 139

Fuses

 Replace, Materials required 161

G

Genotyping run 33

Guest access

 Edit 122

 Overview 42

H

Hands-free, Configure 185

Hands-free features

 Facial authentication 22

 Indicators 23

 Voice activation 22

Home screen, Parts of 24

Home screen notifications 187

I

Import auto export setting 55

Import sample setup 55

Index Term 225

install, desktop software 199

Install block 156

installation, desktop software 198

instrument, connections 197

Instrument

 About 166

 Back up 181

 Edit name 168

 Features 15

 Guest access 42

 Log 178, 179

 Move 165

 Notifications, View 187

 Overview 15

 Parts of 20

 Power off 164

 Power on 43, 164

 Reset, Overview 183

 Restore backup 182

 Restore factory default, Overview 183

 Schedule 75

 Select language 186

 Set date and time 168

 Ship 165

 Software features 26

 Statistics 179

 Store 165

View details 166
 Workflow 40
 Instrument calibration 84
 Instrument comparison 16
 Instrument features, SAE enabled 137
 Instrument log
 Export 179
 View 178
 Instrument name 168
 Instrument profile
 Create 127
 Edit 133
 Facial authentication 132
 Instrument profiles, View all 133
 Instrument service 173
 instrument verification. See RNase P verification
 Interchangeable block calibration 84

L

Language 186
 Light guidelines 192
 limited product warranty 233
 List view 63
 Load plate 51, 52
 load the card 205
 Local area network (LAN) connection 31
 Local instrument profile
 Create 127
 Overview 123
 Lock touchscreen, During run 74

M

maintenance 156
 Manage reagents 66–68
 Manual disk storage management 171
 Melt curve run 33
 Method
 Auto Delta
 Remove 61
 Set up 61
 Data collection location
 Add 58
 Remove 58
 Definitions 57
 Edit 57
 Files 30
 Pause
 Add 62
 Remove 62
 Ramp rates 62
 Stage
 Add 58

Remove 58
 Step
 Add 58
 Remove 58
 VeriFlex Zone
 Remove 59
 Set up 59
 View during run 73
 Method elements 57
 Method tab
 Actions, Edit method 57
 Move instrument 165
 My instrument 27

N

network
 configurations supported 199
 guidelines and best practices 201
 network configuration and security 14
 Network connections 31
 Network drive 27
 networking 199
 No template control (NTC) 65, 66
 Notifications, Clear 188

O

Order to set up profiles 123

P

Pass criteria, RNase P verification 104, 106
 Passive reference dye, Edit 68
 Password, Change 133
 password security 14
 Pause
 Add 62
 Remove 62
 Pause run 72
 PCR cycles 57, 58
 PCR stage
 Add 58
 Remove 58
 PCR stages 58
 PCR step
 Add 58
 Remove 58
 performance verification. See RNase P verification
 PIN
 Change 133
 Sign in 43
 PIN reset 134
 Plate

- Files 30
- Load 51, 52
- Unload 52
- Plate file
 - Copy 77
 - Create and run a plate 47
 - Delete 78
 - Layout 63
 - Method 56
 - Properties 55
 - Repeat last run 49
- Plate layout, Edit 65, 66
- Plate tab, Actions 65, 66
- Plate view 63
- Plots
 - View during run 73
 - Zoom 74
- Post-read 27
- Power off instrument 164
- Power on 43, 164
- Power-saving options 169
- prepare the card
 - centrifuge 205
 - guidelines 203
 - load reaction mix 205
 - seal 207
- Presence/absence run 33
- Profile
 - Add 134
 - Administrator, Create 131
 - Connect
 - Link 126
 - Unlink 129
 - Create 127
 - Delete 135
 - Delete PIN 134
 - Edit 133
 - Facial authentication 131, 132
 - Reset PIN 134
 - Sign out 46
- Profiles
 - Connect
 - Administrator 124
 - Standard 124
 - Order to set up 123
 - SAE 141
 - Sign in 43, 44
 - View all 133
 - Voice commands 45
- Properties, Files 30
- Properties of a plate file 55
- Properties tab
 - Actions 55
 - Import sample setup 55
- Proxy server 170
- Public 27

R

- Ramp rates, Configure 62
- Reaction volume 57
- Reagents
 - Add 66, 68
 - Apply 67, 68
 - Delete 67
 - Edit 66–68
- Region, Connect platform 170
- related documentation 232
- Relative standard curve run 32
- Remote monitoring 180
- Remove Auto Delta 61
- Remove block 158
- Remove data collection location 58
- Remove pause 62
- Remove PCR stage 58
- Remove PCR step 58
- Remove VeriFlex Zones 59
- Repair software 181
- Repeat last run 49
- Replace fuses, Materials required 161
- Required materials, Calibration 90, 98
- Reset instrument 183
- Reset PIN 134
- Restore backup 182
- Restore factory default 183
- RNase P plates 103, 109
- RNase P verification
 - Pass criteria 106
 - Plate description 104
 - Run 107, 112
 - When to perform 85
- RNase P verification, plate, prepare 106
- ROI/uniformity calibration
 - Description 86
 - Frequency 88
 - plate, prepare 91
 - Reminders 88
 - Status 87
- ROI/Uniformity calibration
 - Run 91, 100
 - When to perform 85
- Roles, SAE 139
- Run
 - Lock touchscreen 74
 - Pause 72
 - Post-PCR 18
 - Real-time 18
 - Stop 72

Unlock touchscreen 75
 Run data file name 47
 Run queue 27
 Run type
 Comparative Ct 18
 Genotyping 18
 Melt curve 18
 Presence/absence 18
 Relative standard curve 18
 Standard curve 18
 Run types
 Comparative Ct 32
 Genotyping 33
 Melt curve 33
 Presence/absence 33
 Relative standard curve 32
 Standard curve 31

S

SAE
 dialog boxes 140
 disable 145
 enable 145
 Enable, Workflow 143
 enable on instrument 143
 Functions 138, 139
 instrument features 137
 overview 137
 Profiles 141
 Roles 139
 server, specify 143
 Workflow, Enable SAE 143
 SAE server, connect 144
 safety, biohazard 231
 safety symbols 215, 216
 sample block, decontaminate 159
 Sample information 65, 66
 Sample setup, Import 55
 Saved plate file, Create and run a plate 47
 Schedule, Overview 75
 Schedule run 75
 seal the card 207
 Self-verification test, Results 174
 Service, Request 173
 Set up Auto Delta 61
 Set up VeriFlex Zones 59
 Ship instrument 165
 Sign in 43, 44
 Sign out 46
 Sign-in required 122
 Sleep mode
 Edit 163
 Enable 163

Smart Help
 Instrument service 173
 Technical support 172, 183
 Software, Repair 181
 Software features 26
 Speaker volume, Adjust 169
 Standard 65, 66
 Standard curve run 31
 Standard profile 124
 Statistics, View 179
 Stop run 72
 Storage space 171
 Store instrument 165
 support, customer and technical 233
 symbols
 conformity 220
 control and connection 220
 safety 215, 216
 system components, SAE enabled 141
 System dyes 16
 System template, Create and run a plate 47
 System templates 34, 35, 37, 38

T

TaqMan Array Plates with RFID certificate, Upload 182
 Target
 Add 66
 Edit 65, 66
 Target/SNP 65, 66
 Technical support 172, 183
 Temperature 57
 Template 27
 terms and conditions 233
 Thermal protocol 57
 Thermo Fisher Cloud 26
 Thermo Fisher Connect 26
 third-party software 31, 202
 Time
 Format 168
 Set 168
 Time remaining, View during run 74
 Transfer calibration data files 94
 Transfer calibration results 94
 Transfer data file
 Immediately after run 81
 Later time 81

U

uniformity calibration, plate, prepare 91
 Uniformity calibration
 Description 86

Frequency 88
Reminders 88
Status 87

Unknown 65, 66
Unload plate 52
Unlock touchscreen, During run 75
Upload certificate 182
USB drive 27

V

Verification
Description 104
Pass criteria 106
RNase P 103, 109
Schedule 85

verification failure, troubleshoot 108

VeriFlex Zones
Overview 18
Remove 59
Set up 59

View
List of data files 82
Plate layout 63

View all profiles 133

View audit records (SAE) 153
View calibration images 93, 101
View data during run, Zoom 74

View EULA 167
View instrument details 166
View instrument log 178
View license agreement 167
View method during run 73
View notifications 187
View plots during run, Zoom 74
View privacy policy 167
View time remaining during run 74
Voice commands
Indicators 23
Overview 22
Supported 45

W

warranty 233
Well
Add assay 66
Add target 66
Assign 65, 66
Edit 65, 66
Wells, View 63
Wired connection 31
Wireless connection 31
Workflow
Calibration 90, 97
Custom dye calibration 114

thermofisher.com/support | thermofisher.com/askaquestion
thermofisher.com

28 July 2021

ThermoFisher
SCIENTIFIC